VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate www.vevor.com/support

LAB SCALES AND BALANCES

USER MANUAL

We continue to be committed to provide you tools with competitive price.

"Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

MODEL: SF-400D

Photo for reference

NEED HELP? CONTACT US!

Have product questions? Need technical support? Please feel free to contact us:

Technical Support and E-Warranty Certificate
www.vevor.com/support

This is the original instruction, please read all manual instructions carefully before operating. VEVOR reserves a clear interpretation of our user manual. The appearance of the product shall be subject to the product you received. Please forgive us that we won't inform you again if there are any technology or software updates on our product.

SAFETY INSTRUCTION AND PRECAUTIONS

SPECIFICATIONS

Capacity:600g

Accuracy: 0.01g

Units: g/ct/dwt/gn/T/oz

Tare range: Tare full capacity

Auto off (default): 60 seconds off

Operation temperature:5-35° C

Display: LCD 5 digits

OPERATION(ON/OFF)

1. Place Scale on a horizontal flat surface, press ON/OFF Switch
2. Wait until "0.0" is displayed
3. Put the object(s) on the weighing platform

COUNTING FUNCTION(PCS):

1. Put the relevant amount of goods that you are going to count sample (25,50,75,100).
2. Press and hold “pcs” key for two seconds, then you will see “100” .
3. Press “MODE” key and you will see “25,50,75,100” .Please select the number that you are willing to put as the sample number. Then press “PCS” key to confirm it. Now you can put the goods that you want to count on the center of the platform and the LCD will show the actual number of the goods on the platform.

MODE

According to various needs of the model unit, press the key of “mode” , and you can choose the unit of g/oz/lb/kg.

TARE WEIGHING

1. Turn on scale as described above
2. Place the "tare item" on the platform
3. Press "TARE" and wait until "0.0"(or "0")is displayed
4. Add the "net-weight-item"

CALIBRATION

Standard calibration :Turn on the scale by pressing "CAL" key and hold it until LCD displays then untie "CAL" key, press "CAL" again, after LCD flashes "CAL" and "100.0"(subject to the capacity of your scale)three times.Add 100g(subject to the scale of your scale) standard calweight, wait 3-4 seconds, LCD displays "PASS" , then "100.0" . Put away calibration

weight, press ON/OFF key to turn off scale, the calibration is finished*

Avoiding any exposure to extreme heat or cold, your scale works better when operated at normal room temperature. Always allow the scale to acclimate to normal room temperature for at least one hour before use.

* Keep your scale in a clean environment. Dust, dirt, moisture,vibration, air currents and a close distance to other electronic equipment can all cause an adverse effect on the reliability and accuracy of your scale.*Handle with care. Gently apply all items to be weighed onto tray top. Although this scale is designed to be quite durable, try to avoid rough treatment as this may permanently damage the internal sensor.

* Avoid shaking, dropping or otherwise shocking the scale. This is a precision instrument and must be handled with extreme care.

Special precautions: (If customers do not have their own standard weights for all of our products, do not press the "CAL" button to calibrate them. Our products are calibrated before leaving the factory)

Matching kit

1. Instruction manual * 1
2. Power adapter * 1;
3. Electronic scale body*1
4. Metal tray * 1
5. Plastic foot pads * 4,
6. USB to DC cable * 1(only for AUS model)

Choice of power source:

- 1、North America plug: AC100~240V - 50/60Hz
- 2、European plug: AC100~240V - 50/60Hz
- 3、Australian plug: AC100~240V - 50/60Hz

WARNING:

Install the battery or charge it. If the electronic scale requires battery power, install the battery correctly or charge it according to the instructions.

Placement position. Place the electronic scale on a flat, stable, and dry surface, avoiding damp and high temperature environments.

Avoid vibration. When using, avoid vibration sources around, such as direct sunlight or electromagnetic interference, to improve measurement accuracy.

Avoid overloading. Do not place objects on the electronic scale that exceed its load-bearing range. Pay attention to environmental factors. Avoid exposing electronic scales to extreme temperatures, humidity, or corrosive gas environments.

Keep clean. Regularly wipe the surface of the electronic scale with a dry cloth to avoid the accumulation of dust or dirt.

Proper use and shutdown. After use, turn off the electronic scale in the correct order or wait for it to automatically shut down to save 17 battery energy.

Regular calibration. Regularly calibrate the electronic scale to ensure measurement accuracy.

Properly stored. When not in use, store the electronic scale in a dry, well ventilated environment without direct sunlight. ten thousand

Check the power supply and battery. If the electronic scale cannot be turned on or charged, check if the power cord, battery, or external power supply is working properly.

FCC Information:

CAUTION: Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment!

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions:

- 1) This product may cause harmful interference.
- 2) This product must accept any interference received, including interference that may cause undesired operation.

WARNING: Changes or modifications to this product not expressly approved by the party responsible for compliance could void the user's authority to operate the product.

Note: This product has been tested and found to comply with the limits for a Class B digital device pursuant to Part 15 of the FCC Rules, These limits are designed to provide reasonable protection against harmful interference in a residential installation.

This product generates, uses and can radiate radio frequency energy, and if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this product does cause

harmful interference to radio or television reception, which can be determined by turning the product off and on, the user is encouraged to try to correct the interference by one or more of the following measures.

- Reorient or relocate the receiving antenna.
- Increase the distance between the product and receiver.
- Connect the product to an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for assistance.

CORRECT DISPOSAL

This product is subject to the provision of European Directive 2012/19/EU. The symbol showing a wheeled bin crossed through indicates that the product requires separate refuse collection in the European Union. This applies to the product and all accessories marked with this symbol. Products marked as such may not be discarded with normal domestic waste, but must be taken to a collection point for recycling electrical and electronic devices.

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YH CONSULTING LIMITED.
C/O YH Consulting Limited Office 147, Centurion House,
London Road, Staines-upon-Thames, Surrey, TW18 4AX

VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate

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We continue to be committed to provide you tools with competitive price. "Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

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Geïmporteerd naar AUS: SIHAO PTY LTD. 1 ROKEVA
Geïmporteerd naar de VS: Sanven Technology Ltd. Suite 250, 9166 Anaheim
E-CrossStu GmbH