VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate www.vevor.com/support

LAB DISTILLATION APPARATUS USER MANUAL

MODEL:YHEO-3

We continue to be committed to provide you tools with competitive price.

"Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

MODEL:YHEO-3

NEED HELP? CONTACT US!

Have product questions? Need technical support? Please feel free to contact us:

Technical Support and E-Warranty Certificate
www.vevor.com/support

This is the original instruction, please read all manual instructions carefully before operating. VEVOR reserves a clear interpretation of our user manual. The appearance of the product shall be subject to the product you received. Please forgive us that we won't inform you again if there are any technology or software updates on our product.

SAFETY INSTRUCTIONS

WARNING:

Read this material before using this product. Failure to do so can result in serious injury.

Assembly precautions

1. Assemble only according to these instructions. Improper assembly can create hazards.
2. Wear safety goggles and duty work gloves during assembly.
3. Keep the assembly area clean and well-lit.
4. Keep bystanders out of the area during assembly.
5. Do not assemble when tired or when under the influence of alcohol, drugs or medication.
6. The product capabilities apply to properly and completely assembled products only.
7. Assemble on a flat, level, hard and smooth surface capable of safely supporting the Lab Distillation Apparatus.
8. For additional information regarding the parts listed in the following pages, please refer to the Assembly Diagram of this manual. Unwrap and separate all parts in a clean work area.

Use precautions

1. DO NOT SIT OR STAND ON THIS ITEM.
2. This product is not a toy. Do not allow children to play with or near this item.
3. Do not exceed weight capacities.
4. Use only on a flat, level, hard and smooth surface capable of safely supporting a fully loaded Lab Distillation Apparatus.
5. Use as intended only.
6. Inspect before every use; do not use if parts are loose or damaged.

SAVE THIS MANUAL

INSTRUCTIONS

Thank you very much for choosing this model. Glass fittings should always be handled with care. Please read all of the instructions before using it. The information will help you achieve the best possible results.

Parts name	Number of parts	Description
Flask	1	500ml
Coil condenser	1	/
Pear-shaped Separatory funnel	1	125ml
Conical flask	1	100ml
Glass stopper	2	19#
beaker	1	100ml
Glue head dropper	1	Red glue head
Alcohol burner	1	150ml
thermometer	1	100°C
Rubber plug	4	/
Latex tubing	1	2m
Vacuum grease	1	Vaseline
PH test paper	1	/
Flask brush	1	Small size
Iron support stand	2	With 40cm supporting pole
Flask clamp	2	/
Cross clamp	2	/
Horn tube	1	/
Circular ring	2	1 of 5.1cm, 1 of 7.3cm
Terracotta net	1	/

OPERATION

1. Assemble the iron support stand;


2. Place alcohol burner;

3. Install the circular ring and place the terracotta net on the circular ring ;

4. Install the cross clamp and flask clamp at the appropriate height;

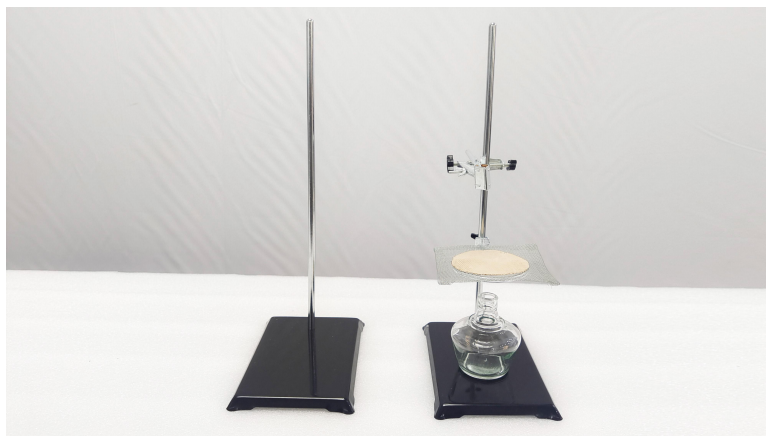5. Install the flask;

6. Install the coil condenser and fix it with the flask clamp;


7. Install the horn tube ;

8. Install the circular ring at the appropriate height of the iron support stand, and place the pear-shaped separatory funnel on it, then place the beaker below the outlet of the pear-shaped separatory funnel;

9. Screw the thermometer into the rubber plug and install the rubber plug in the mouth of the flask.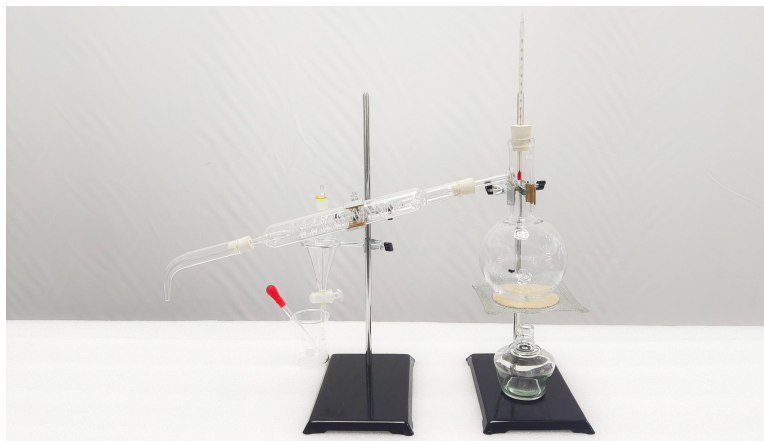

PRODUCT PARAMETER

Model	YHEO-3
Heating Method	Alcohol Burner
Flask Type	Distilling Flask

Manufacturer: Shanghaimuxinmuyeyouxiangongsi
Address: Shuangchenglu 803nong11hao1602A-1609shi,
baoshanqu, shanghai 200000 CN.
Imported to AUS: SIHAO PTY LTD. 1 ROKEVA STREET-
EASTWOOD NSW 2122 Australia
Imported to USA: Sanven Technology Ltd. Suite 250,
9166 Anaheim Place, Rancho Cucamonga, CA 91730

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YH CONSULTING LIMITED.
C/O YH Consulting Limited Office 147,
Centurion House, London Road,
Staines-upon-Thames, Surrey, TW18 4AX

VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate

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Importerad till AUS: SIHAO PTY LTD. 1 ROKEVA STREET-EASTWOOD
Importerad till USA: Sanven Technology Ltd. Suite 250, 9166
E-CrossStu GmbH