VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate www.vevor.com/support

LAB SCALES AND BALANCES USER MANUAL

We continue to be committed to provide you tools with competitive price.

"Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

VEVOR®

TOUGH TOOLS, HALF PRICE

LAB SCALES AND BAL
ANCES

MODEL: HZ5003/HZ-B20002/HZ-B30002

Photo for reference

NEED HELP? CONTACT US!

Have product questions? Need technical support? Please feel free to contact us:

Technical Support and E-Warranty Certificate
www.vevor.com/support

This is the original instruction, please read all manual instructions carefully before operating. VEVOR reserves a clear interpretation of our user manual. The appearance of the product shall be subject to the product you received. Please forgive us that we won't inform you again if there are any technology or software updates on our product.

Key Description:

ON/OFF - "On/Off" key PCS - "Count" key

TARE - "Peel off" key UNIT - "Unit conversion" key

CAL - "Calibration" key ZERO - "Zero" key

Specifications and performance indicators

The HZ-B series electronic balance is divided into the following specifications according to different ranges ()

Model HZ-B	5003	20002	30002
Maximum weighing	500g	2000g	3000g
resolving power	0.001g	0.01g	0.01g
Calibration scale	10d	10d	10d
tare range	500g	2000g	3000g
Calibration weight	200g	500g	2000g
Pan size	90mm	square 157*176cm	
Power	(AC110-220V±10% 50HZ/601HZ)AC/DC dual-purpose		
Operating temperature	-5-45℃		
Using humidity	50-85%		

Operation preparation ZERO - "Zero" key

Place the balance on a stable and flat workbench, adjust the adjustment feet so that the horizontal bubble is in the center position, and avoid balance vibration, sunlight exposure, airflow, and strong electromagnetic wave interference.

Power on

1. Connect the external power supply and turn on the power switch.
2. Press the "ON/OFF" button to display "8.8.8.8.", "maximum weighing value", and "--" in sequence. The display time of "-- --" depends on the stability of the sensor, so do not place the balance on the air vent or unstable workbench. Finally, display the weighing mode of "0, 0.0, or 0.00". If the "0" symbol in the upper left

corner flashes, it indicates that the workplace is unstable and the following 5 to 9 operations cannot be performed.

Calibration level

1. Place the balance on a stable and flat workbench, while avoiding vibration, sunlight exposure, airflow, and strong electromagnetic interference. Turn on the power of the balance and preheat for half an hour before proceeding to ensure accurate calibration results.

2. Calibration operation CAL - "Calibration" key

Calibration: Without adding any objects to the scale, hold down the "CAL Calibration" button and hold it down. After about 3 seconds, release it immediately when "- CAL -" is displayed. Wait for the flashing "Standard Weight Value" to be displayed. Place the shining "Standard Weight Value" weight on the scale and display the "-----" waiting state. Wait for the "Standard Weight Value" to be displayed. Remove the weight and the display will change to the "-----" waiting state. Wait for the display to be 0, 0.0, or 0.00, and the calibration will end. If the weighing is still inaccurate after calibration, repeat the above process several times for calibration.

3. Linear correction using 3000g/0.01g as an example

Press and hold the "CAL Calibration" button without releasing your hand. After about 3 seconds, when the "-- CAL --" display is displayed, release your hand immediately. Wait for the "Standard Weight Value" to flash, then press and hold the "PCS" button without releasing your hand. After about 3 seconds, when the "-- CAL --" display is displayed, release your hand immediately, and the 1000g flash will be displayed. At this time, you put in the 1000g weight, remove it when it stabilizes, and then the 3000g flash will jump out. At this time, you put in the 3000g weight, remove it when it stabilizes, and linear calibration is completed

Weighing

1. After the startup preheating is stable or calibrated, the weighing mode is displayed as "0 or 0.0 or 0.00".

2. Place the weighing object on the scale, which displays its weight.

Tare removal key

1. Place the container on the weighing pan, and the balance will display the weight of the container.
2. After pressing the "TARE Tare" button, it will display "0 or 0.0 or 0.00", indicating that the tare weight has been removed.
3. Place the object in the container, indicating the weight of the object being weighed.

Balance Counting Operation PCS - "Counting" Key

Enter the average state of setting a single counting object, Place the container needed for counting on the scale, and do not place any unnecessary containers. After pressing the "TARE Tare" button, the display will be "0, 0.0, or 0.00". Then, after pressing the "PCS Count" button, the display will be "10" flashing, and the display in the bottom right corner will change from "g" to "PCS". At this time, place 10 objects to be counted on the scale, then press the "PCS Count" button, and the display will change to "-- --" waiting state. After a while, the display will change to "10". After completing the average setting operation, the scale can proceed with counting. When performing this operation, the zero point must be "0 or 0.0 or 0.00", and the weight of a single count must be greater than the minimum reader of the balance, i.e. resolution.

SAFETY INSTRUCTION AND PRECAUTIONS

WARNING:

1. Before use, power should be turned on and preheated according to regulations.
2. The sum of the tare weight and the weight of the weighing object shall not exceed the weighing range.
3. If the weighing is not accurate, it is necessary to calibrate the balance with standard weights.
4. If you need to remove the circular scale plate from the balance, please turn the scale plate clockwise before removing it. Do not forcefully pull the scale plate upwards to avoid damaging the sensor.
5. Error message appears, please refer to the table below

Error message	Explain	Processing method
upper horizontal line	The weight on the balance is biased negative	Recalibrate with weights
Lower horizontal line	Overweight alarm, Article 10	Recalibrate with weights
ERR-1	Continuous power on/off caused	Turn off for 3 seconds before turning on again
ERR-2	Weighing unstable	Just wait for a moment

Matching kit

HZ-B20002 and HZ-B30002

1. Instructions * 1
2. Power adapter * 1;
3. 500g weight * 1;
4. Weighing plate * 1;
5. Weighing tray bracket * 1
6. Internal hexagonal wrench * 1
7. 4 bolts

HZ5003

- 1 Instructions*1
2. Power adapter*1
3. 200g weight * 1;
4. Weighing plate * 1;
5. Windproof hood base * 1
6. Windshield (composed of 5 acrylic panels)

Option: 1 standard weight

The relevant content of this manual may vary due to product updates or differences from the actual product. Please refer to the actual product without further notice.

Special precautions: (If customers do not have their own standard weights for all of our products, do not press the "CAL" button to calibrate them. Our products are calibrated before leaving the factory)

A scale with a specification of 3000g/0.01g and a weighing range of 0-3KG should be installed before powering on. When installing the scale, a bracket must be installed, with a distinction between the front and back sides. We have labels on it, and the scale can be placed freely in the front and back (scale size: 176*157mm stainless steel material). After installing the scale, the scale should be straightened. If it is uneven, the four adjustment feet at the bottom of the scale can be adjusted (usually by twisting the four feet to the innermost part) to prevent the scale from swinging. After completing these steps, turn on the machine and wait for "0.00" to appear before it can be weighed normally.

This product has a wide range of applications, mainly in laboratories, gold and jewelry stores, chemical, metallurgical, mining, textile, teaching, etc. It is important to avoid working environments (adverse environments include "wind, vibration, magnetic field", etc.)

The ZERO key of the product is the peeling and zeroing key, with the highest frequency of use. The other UNIT keys are some unit conversions, mainly including g, ct, oz, ozt, lb, %, etc. We will calibrate this product before shipping, and the accuracy of the product meets the requirements. Customers can use it with confidence after receiving it. If there are any problems during use, please feel free to contact us, and we will patiently guide and maintain it. In addition, this product cannot be weighed excessively ("-----" overweight warning), and should not be soaked in water or dropped. It is best to return it to its original packaging after use to ensure the long-term use of the product.

FCC Information:

CAUTION: Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment!

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions:

- 1) This product may cause harmful interference.
- 2) This product must accept any interference received, including interference that may cause undesired operation.

WARNING: Changes or modifications to this product not expressly approved by the party responsible for compliance could void the user's authority to operate the product.

Note: This product has been tested and found to comply with the limits for a Class B digital device pursuant to Part 15 of the FCC Rules, These limits are designed to provide reasonable protection against harmful interference in a residential installation.

This product generates, uses and can radiate radio frequency energy, and if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this product does cause harmful interference to radio or television reception, which can be determined by turning the product off and on, the user is encouraged to try to correct the interference by one or more of the following measures.

- Reorient or relocate the receiving antenna.
- Increase the distance between the product and receiver.
- Connect the product to an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for assistance.

CORRECT DISPOSAL

This product is subject to the provision of European Directive 2012/19/EU. The symbol showing a wheeled bin crossed through indicates that the product requires separate refuse collection in the European Union. This applies to the product and all accessories marked with this symbol. Products marked as such may not be discarded with normal domestic waste, but must be taken to a collection point for recycling electrical and electronic devices.

Manufacturer:Shanghaimuxinmuyeyouxiangongsi
Address: Shuangchenglu 803nong11hao1602A-1609shi, baoshanqu, shanghai 200000 CN.
Imported to AUS: SIHAO PTY LTD. 1 ROKEVA STREETEASTWOOD NSW 2122 Australia
Imported to USA: Sanven Technology Ltd. Suite 250, 9166 Anaheim Place, Rancho Cucamonga, CA 91730

EC	REP
----	-----

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YH CONSULTING LIMITED.
C/O YH Consulting Limited Office 147, Centurion House,
London Road, Staines-upon-Thames, Surrey, TW18 4AX

VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate

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We continue to be committed to provide you tools with competitive price. "Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

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ON/OFF - "On/Off"-tangent PCS - "Count"-tangent

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Importado a AUS: SIHAO PTY LTD. 1 ROKEVA STREET ASTWOOD NSW 2122 Australia
Importado a EE. UU.: Sanven Technology Ltd. Suite 250, 9166 Anaheim Place, Rancho Cucamonga, CA 91730
E-CrossStu GmbH