VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate www.vevor.com/support

HUMAN BODY TEACHING MODELS

USER MANUAL

MODEL:LD-203

We continue to be committed to provide you tools with competitive price.

"Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

VEVOR®

TOUGH TOOLS, HALF PRICE

HUMAN BODY TEACHING
MODELS

MODEL:LD-203

NEED HELP? CONTACT US!

Have product questions? Need technical support? Please feel free to contact us:

Technical Support and E-Warranty Certificate
www.vevor.com/support

This is the original instruction, please read all manual instructions carefully before operating. VEVOR reserves a clear interpretation of our user manual. The appearance of the product shall be subject to the product you received. Please forgive us that we won't inform you again if there are any technology or software updates on our product.

SAFETY INSTRUCTIONS

WARNING:

1. Read these instructions before using this product. Failure to do so can result in serious injury.
2. After use, please assemble all the parts and keep them vertically.
3. Avoid high temperature and direct sunlight.
4. Remove minor stains with neutral detergent. The use of alcohol, benzene and other substances may damage the product.

SAVE THESE INSTRUCTIONS

PRODUCT PARAMETER

Model	LD-203
Material	PVC
Removable Parts	15

INSTRUCTIONS

1. Lay the Torso Anatomy Model Flat on its Back on a Flat Surface

2. Add the Posterior(Back)Half of the Right Lung to the Right Chest

3. Add the Posterior(Back)Half of the Left Lung to the Left Chest

4. Add the Trachea / Esophagus / Aorta Piece Between the Lungs

5. Next Add The Heart on Top of the Trachea Piece

6. Add the Anterior(Front) Right Lung to the Right Cheae

7. Add the Anterior (Front) Left Lung to the Left Chest

8. Add the Diaphragm Piece Right Underneath the Lungs

9. Add the Liver Right Underneath the Diaphragm

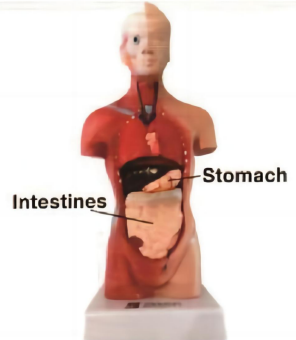
10. Add the Pancreas / Duodenum / Spleen to Upper Right Abdomen

10. CONT. Location of Pancreas / Duodenum/Spleen in Abdomen

11. Add the Intestines and Stomach

12. Face the Torso Model Downward on a Flat Surface

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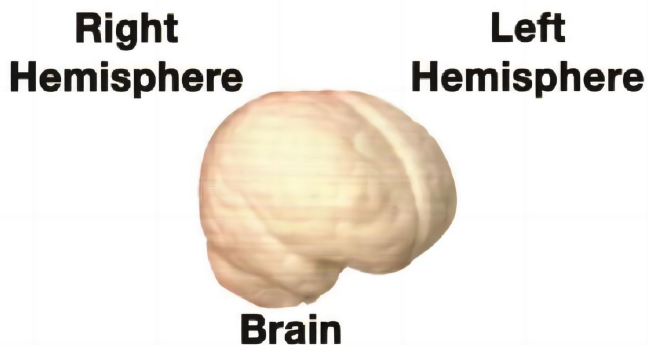
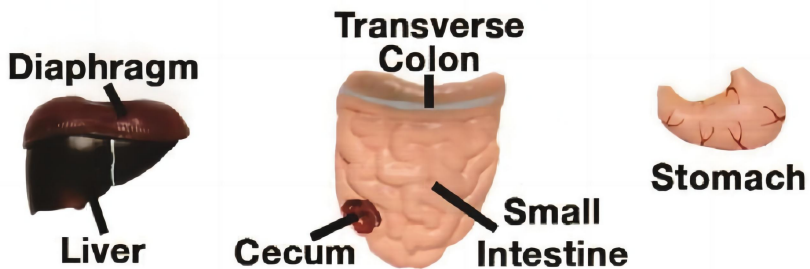
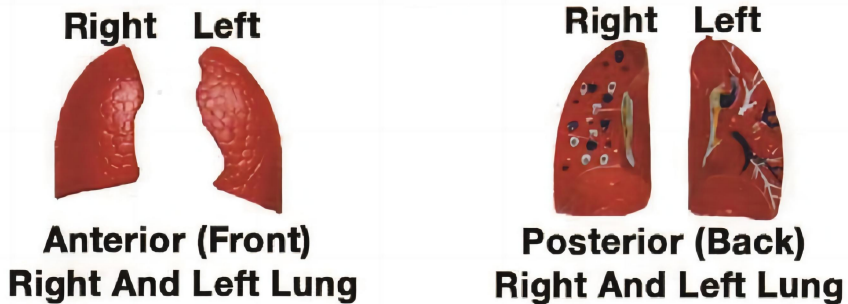
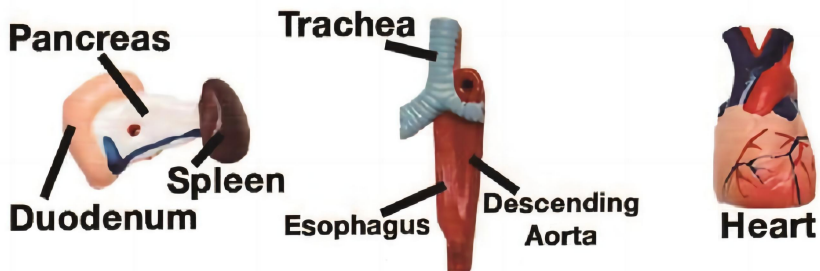
YH CONSULTING LIMITED.
C/O YH Consulting Limited Office 147, Centurion House,
London Road, Staines-upon-Thames, Surrey, TW18 4AX

VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate

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Imported to AUS: SIHAO PTY LTD. 1 ROKEVA STREET-
Imported to USA: Sanven Technology Ltd. Suite 250, 9166
E-CrossStu GmbH