## USER MANUAL

Bedienungsanleitung  
Instrukcja obsługi  
Návod k použití  
Manuel d'utilisation  
Istruzioni per l'uso  
Manual de instrucciones  
Használati útmutató  
Bruksanvisning  
Käyttöohje  
Gebruiksaanwijzing  
Bruksanvisning  
Instruções de utilização  
Používateľská príručka  
Ръководство за употреба  
Οδηγίες χρήσης  
Urpute za uporabu  
Naudojimo instrukcija  
Manual de utilizare  
Navodila za uporabo

---

---

## OVERHEAD STIRRER

<b>DE</b>	Produktname	Laborrührer
<b>EN</b>	Product name	overhead stirrer
<b>PL</b>	Nazwa produktu	Mieszadło laboratoryjne
<b>CZ</b>	Název výrobku	Horní míchačka
<b>FR</b>	Nom du produit	Agitateur en tête
<b>IT</b>	Nome del prodotto	Agitatore sopraelevato
<b>ES</b>	Nombre del producto	Agitador de cabeza
<b>HU</b>	Termék neve	Fej feletti keverő
<b>DA</b>	Produktnavn	Overhead omrører
<b>FI</b>	Tuotteen nimi	Yläsekoitin
<b>NL</b>	Productnaam	Bovenroerder
<b>NO</b>	Produktnavn	Overhead omrører
<b>SE</b>	Produktnamn	Överliggande omrörare
<b>PT</b>	Nome do produto	Agitador de sobrecabeça
<b>SK</b>	Názov produktu	Horná miešačka
<b>BG</b>	Име на продукта	Горна бъркалка
<b>EL</b>	Όνομα προϊόντος	Κορυφαίος αναδευτήρας
<b>HR</b>	Naziv proizvoda	Gornje miješalo
<b>LT</b>	Produktu pavadinimas	Viršutinis maišytuvas
<b>RO</b>	Numele produsului	Agitator superior
<b>SL</b>	Ime izdelka	Zgornje mešalo
<b>DE</b> Modell   <b>EN</b> Product model   <b>PL</b> Model produktu   <b>CZ</b> Model výrobku   <b>FR</b> Modèle   <b>IT</b> Modello   <b>ES</b> Modelo   <b>HU</b> Modell   <b>DA</b> Model   <b>FI</b> Tuotteen malli   <b>NL</b> Productmodel   <b>NO</b> Produktmodell   <b>SE</b> Produktmodell   <b>PT</b> Modelo do produto   <b>SK</b> Model   <b>BG</b> Модел на продукт   <b>EL</b> Μοντέλο προϊόντος   <b>HR</b> Model proizvoda   <b>LT</b> : Gaminio modelis   <b>RO</b> : Model de produs   <b>SL</b> : Model izdelka		SBS-MR-150
<b>DE</b> Hersteller   <b>EN</b> Manufacturer   <b>PL</b> Producent   <b>CZ</b> Výrobce   <b>FR</b> Fabricant   <b>IT</b> Produttore   <b>ES</b> Fabricante   <b>HU</b> Termelő   <b>DA</b> Producent   <b>FI</b> Valmistaja   <b>NL</b> Producent   <b>NO</b> Produsent   <b>SE</b> Tillverkare   <b>PT</b> Fabricante   <b>SK</b> Výrobca   <b>BG</b> Производител   <b>EL</b> Κατασκευαστής   <b>HR</b> Proizvođač   <b>LT</b> Gamintojas   <b>RO</b> Producător   <b>SL</b> Proizvajalec		expondo Polska sp. z o.o. sp. k.
<b>DE</b> Anschrift des Herstellers   <b>EN</b> Manufacturer Address   <b>PL</b> Adres producenta   <b>CZ</b> Adresa výrobce   <b>FR</b> Adresse du fabricant   <b>IT</b> Indirizzo del produttore   <b>ES</b> Dirección del fabricante   <b>HU</b> A gyártó címe   <b>DA</b> Producentens adresse   <b>FI</b> Valmistajan osoite   <b>NL</b> Adres producent   <b>NO</b> Produsentens adresse   <b>SE</b> Tillverkarens adress   <b>PT</b> Endereço do fabricante   <b>SK</b> Adresa výrobcu   <b>BG</b> Адрес на производителя   <b>EL</b> : Διεύθυνση κατασκευαστή   <b>HR</b> Adresa proizvođača   <b>LT</b> Gamintojo adresas   <b>RO</b> Adresa producătorului   <b>SL</b> Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra   Poland, EU

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

## Technical data

Parameter description	Parameter value
Product name	Overhead Stirrer
Model	SBS-MR-150
Rated voltage [V~] / frequency [Hz]	230 / 50
Rated power [W]	250
Dimensions [width x depth x height; cm]	30 x 20 x 70
Weight [kg]	10
Volume of agitated substance [L]	150
Max. rotation speed [rev/min]	1500
Disc holder dia. [Ø]	1.5 - 13
Distance from disc to upright central tube [mm]	160-270
Max distance of disc [mm]	560
Display	LCD
Stirring rod material	Stainless steel (304)

### Please read and save these instructions.

Read through this owner's manual carefully before using the product. Protect yourself and others by observing all safety information, warnings, and cautions. Failure to comply with instructions could result in personal injury and/or damage to product or property. Please retain instructions for future reference.

Operated with brush-less DC motor and micro-controller at constant speed by closed-loop control, the stirrer is equipped with speed LCD panel, rotary encoder and core paddle.

### **Application method:**

1. Speed setting: turn the knob to set the required speed (up by turning clockwise, and down counterclockwise).
2. Press down the knob when the speed is set to start the stirrer. The set value will keep unchanged until re-setting. Press the knob again to stop during operation.
3. During stirring, turn the knob to change the set speed. LCD panel shows the changing speed by turning. When the knob stops turning, the real speed will be indicated after 1 second. The speed changed during stirring will not be saved.

### **Notes:**

1. During application, the speed should be accelerated slowly. If the linearity of the stir bar is not good, the speed should not be fast or use flexible coupling shaft and fasten it with the central supporting ring.
2. Keep yourself away from the stirrer to prevent clothes or hair being caught in the machine.
3. Pay attention to spill during stirring.

### **Packing list:**

- one set of stirrer machine
- one base plate
- one universal joint
- one upright
- one drill chuck key
- one flexible coupling shaft

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