VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate <https://www.vevor.com/support>

OVERHEAD MIXERS

USER MANUAL

MODEL NO.: FSH-2A, JJ-1

We continue to be committed to provide you tools with competitive price.

"Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

VEVOR®

TOUGH TOOLS, HALF PRICE

Magnetic Stirrers

MODEL NO.: FSH-2A, JJ-1

NEED HELP? CONTACT US!

Have product questions? Need technical support? Please feel free to contact us:

Technical Support and E-Warranty Certificate
www.vevor.com/support

This is the original instruction, please read all manual instructions carefully before operating. VEVOR reserves a clear interpretation of our user manual. The appearance of the product shall be subject to the product you received. Please forgive us that we won't inform you again if there are any technology or software updates on our product.

WARNING:

Please read this manual carefully before using the product. Failure to do so may result in serious injury.

	Warning! ● Read the operating instructions carefully before use. ● Ensure that only trained staff works with the instrument.
	Protective ground contact ! ● Make sure that socket is grounded (protective ground contact) before use.

- When operating wear personal safety protection to avoid the risk of:
 - Splashing and evaporation of liquids
 - Release of toxic or combustible gases
- Set up the instrument in a spacious area on a stable, clean, non-slip, dry and fireproof surface. Do not operate the instrument in explosive atmospheres, with hazardous substances or around water.
- Gradually increase the speed, reduce the speed if:
 - Stirring bar breaks away due to high speed
 - The instrument is not running smoothly, or the container is not centered on the plate
- Be aware of hazards due to:
 - Flammable materials or media with a low boiling temperature
 - Overfilling of media
 - Unsafe container
- Process pathogenic materials only in closed vessels.
 - If the case of the stirrer bar is PTFE, please note:
 - Elemental fluorine, three fluoride and alkali metals will corrode the PTFE and Halogen alkanes making it expand at room temperature.
 - Molten alkali, alkaline earth metals or their solution, as well as the powder in the second and third ethnic of the Periodic Table of Elements will have a chemical reaction with PTFE when temperature reaches 300 ~ 400°C.
- Check the instrument and accessories before each use. Do not use

damaged components. Safe operation is only guaranteed with the accessories listed in the “Accessories” section. Accessories must be securely attached to the device and must be removed when not in use. Always disconnect the power before fitting accessories.

- When the external temperature sensor is needed, the tip of the measuring sensor must be at least 5-10 mm from vessel bottom and wall.
- The instrument can only be fully disconnected from the main power supply by turning off the main or disconnecting the plug.
- The voltage stated on the label must correspond to the main power supply.
- Ensure that the main power supply cable does not touch the hotplate. Do not cover the device.
- The instrument must only be operated by experts.

Keep away from high magnetic fields.

SAVE THESE INSTRUCTIONS

PROPER USE

The instrument is designed for mixing and/or heating liquids in schools, laboratories or factories

- Observe the minimum distances between multiple units, and distances to the rear wall and above the assembly (min. 100 mm).

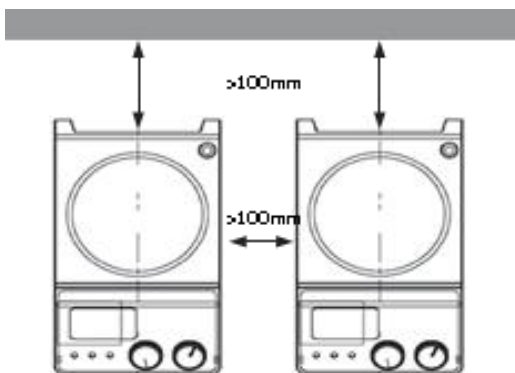
Figure 1

This device is not suitable for use in residential areas or other constraints mentioned in Safety Instructions.

TECHNICAL SPECIFICATIONS

Model	FSH-2A	JJ-1
Voltage [VAC]	220~240	220~240
Frequency [Hz]	50	50
Power [W]	185	100
Maximum stirring capacity [mL]	1000	1000
Product Dimensions (L*W*H)[mm]	265*165*460	200*125*200
Packing Size (L*W*H) [mm]	340*250*460	430*320*160
Net Weight [kg]	4.6	6.8
Gross Weight [kg]	5	7.3

*Products such as specifications, appearance, and design are subject to modification without prior notice.

PACKAGE CONTENTS

MODEL NO.: FSH-2A


Fixed Magnet*3

Glass*4

1 set of wrench


Mixing cutter head*2

MODEL NO.: JJ-1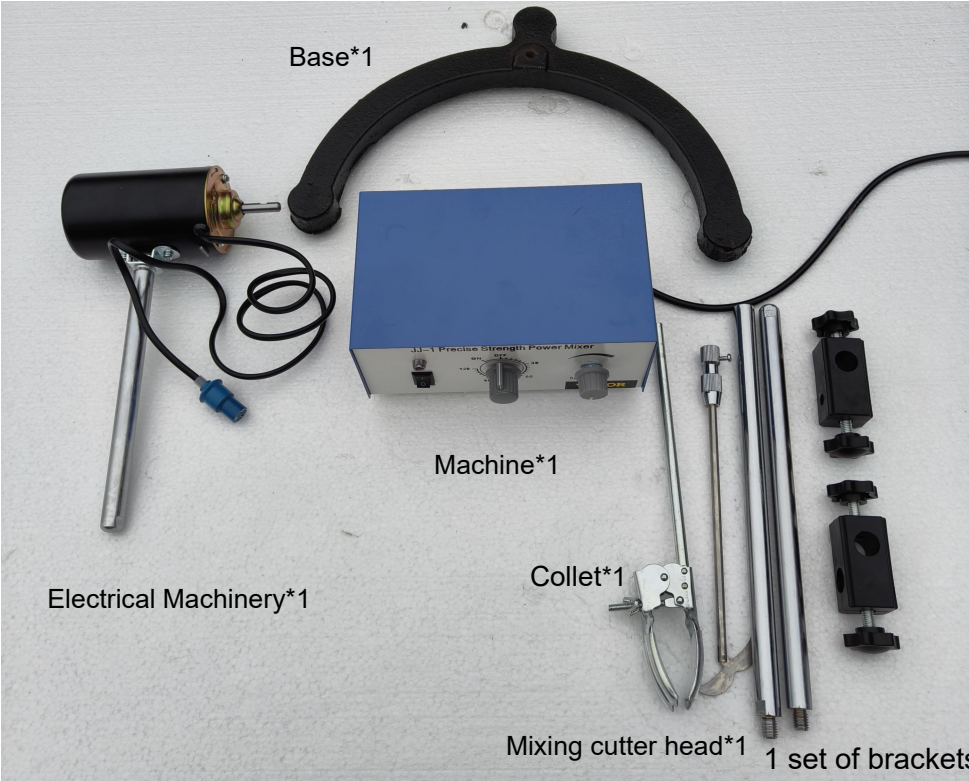

PRODUCT FUNCTIONS

MODEL NO.: FSH-2A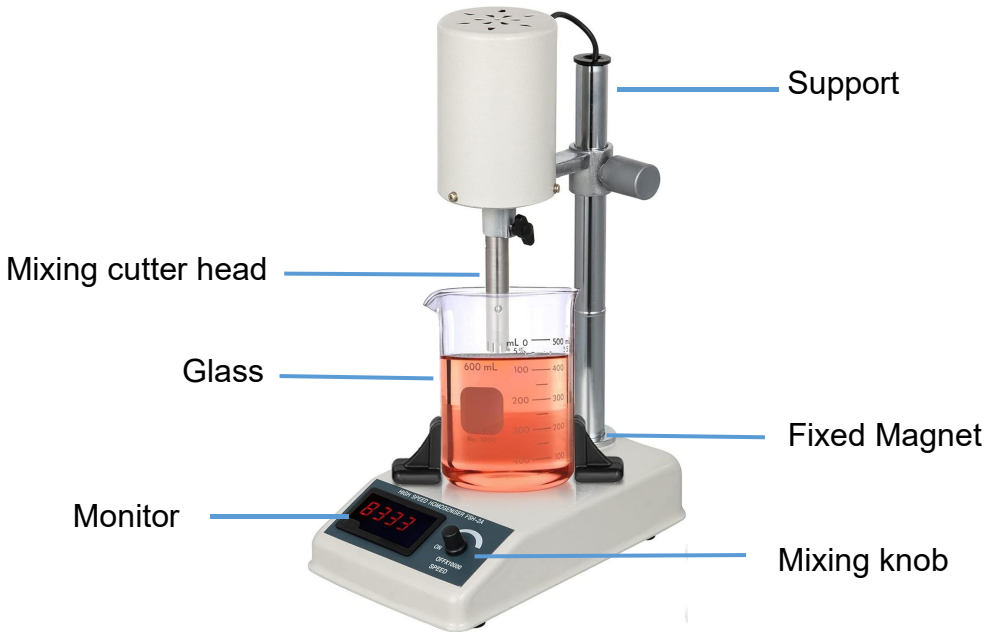

Operation mode:

1. Place the machine on a horizontal desktop, and then connect the power cord to the power outlet on the back of the base.
2. Plug the other end of the power cord into a standard socket.
3. Adjust the speed by turning the knob, and the current speed will be displayed on the display screen.

Precautions for use


The instrument is placed on a horizontal worktable, and the required homogenizing cutter must be in place. When the power supply is connected to the test machine, the plum knife head is submerged in the water with a cup or utensil filled with water. The power switch is turned on, the speed knob is adjusted slowly to fast, and there is no obvious metal friction (such as metal friction, switch off, loosen the cutter 6mm tightening bracelets to rotate the fixing jacket randomly). Turn up and install in place, then tighten up. You can't work until you've finished. Cut the plants or organisms that need to be tested into small pieces or pieces, place them in randomly equipped glassware, clamp them in the middle of the magnet bracket (heptagon), loosen the 8mm fastening bracelets of the fixed motor base, slowly move down the orifice of the utensil, make the head of the homogenate 2-3mm from the bottom of the utensil advantageous, and tighten the 8mm bracelets again. The working time of the experiment should not exceed 3minutes at a time.

Maintenance instructions:

After long-term use, the polyfluoroethylene in the tool of high-speed homogenizer is a wear part, which can be replaced by the user, and is the assembly procedure of the homogenizer.

Be careful: After each work, remove the cutter and rinse it with clean water. Firstly, insert the dispersing head into the polyfluoroethylene, when the step of the dispersing head exceeds the length of the sleeve by 0.2mm, if it is found that the sleeve is too long, one end of the small head can be lowered to size by a cylindrical knife, then the polyfluoroethylene sleeve can be pressed into the fixed head by hand, and then the dispersing head can be inserted into the sleeve. At this time, the dispersing head must rotate freely in the sleeve. If it is found that the sleeve is tight, it can be inserted and pulled in the sleeve with the drill bit of $\phi 5$. After pulling, insert the dispersing head into the sleeve and tighten the small rotating axle (when twisting, the force must be appropriate, otherwise it will cause slippage), and finally tighten the fixing sleeve and the fixing head.

MODEL NO.: JJ-1

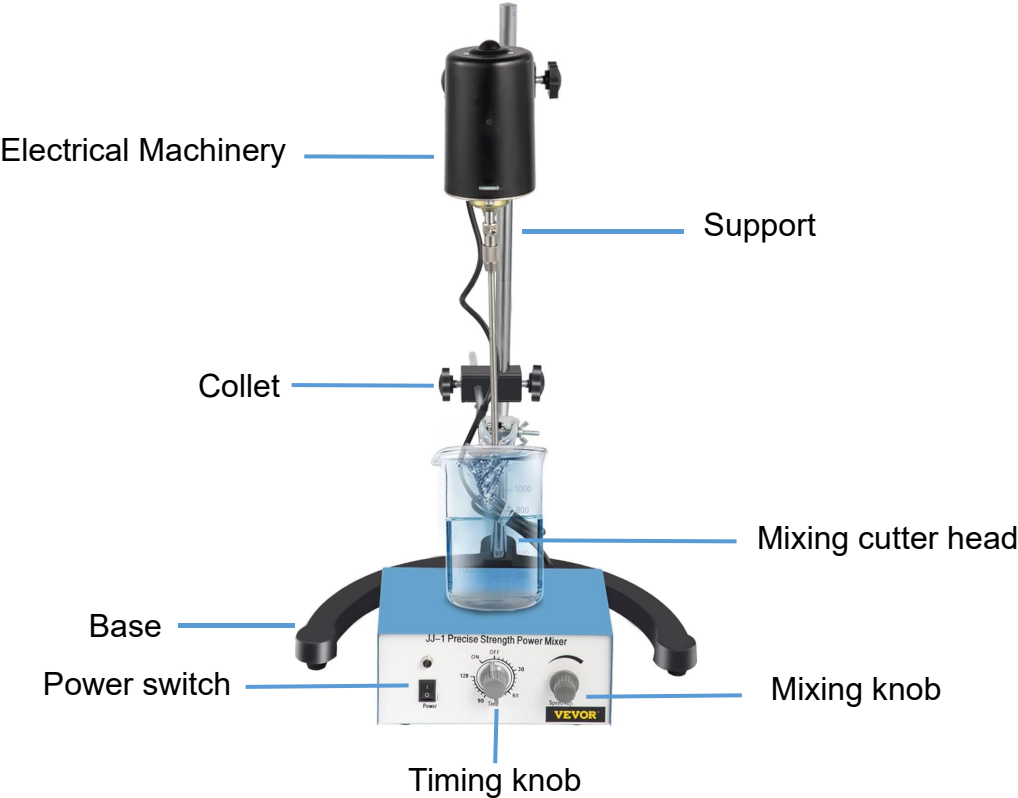Operation mode:

1. Place the machine on a horizontal desktop, and then connect the power cord to the power outlet on the back of the base.
2. Plug the other end of the power cord into a standard socket.
3. Pay attention to the power switch on the right side of the base. The line (-) indicates on, and (o) indicates OFF.
4. Place the machine stably, put the beaker with the liquid medicine in the center of the working plate, put an appropriate rotor in the beaker and make it in the center of the beaker, turn on the power supply, adjust it to the required speed by rotating the "Speed Adjust" knob.
5. If timing is required, turn the timing knob to the required timing time, and then start the countdown. Timing range: 0 - 120 min.

Precautions for use

The speed control knob is preset at the highest speed at the factory to protect the drive system from damage during transportation. Therefore, check the setting of the knob before use to ensure that it is suitable for the liquid being stirred: if the speed cannot be determined, turn the knob to a minimum. Mixer after a period of time is not used, or reused, the initial connection will hear friction noise, this is due to the friction wheel lining of the prestressed, no damage to the function of mixer, noise will disappear after a short run. The rotary chuck and mixing shaft allow the clamping mixing rod to be inserted with a diameter of up to 12 mm.

Maintenance instructions:

Electric mixers can only fail after long-term natural wear and aging of parts. When cleaning, use water only and contain surfactants. Oil or grease substances are prohibited from contacting the surfaces of the friction wheel and the tapered wheel. This greatly reduces the coefficient of friction between the two, thereby weakening the power transmission capability.

This product is subject to the provision of European Directive 2012/19/EC. The symbol showing a wheeled bin crossed through indicates that the product requires separate refuse collection in the European Union. This applies to the product and all accessories marked with this symbol. Products marked as such may not be discarded with normal domestic waste, but must be taken to a collection point for recycling electrical and electronic devices.

Manufacturer: Shanghaimuxinmuyeyouxiangongsi
Address: Baoshanqu Shuangchenglu 803long 11hao 1602A-1609shi
Shanghai
Imported to AUS: SIHAO PTY LTD. 1 ROKEVA STREETEASTWOOD
NSW 2122 Australia
Imported to USA: Sanven Technology Ltd.
Suite 250,9166 Anaheim Place, Rancho Cucamonga, CA 91730

EC	REP	SHUNSHUN GmbH Römeräcker 9 Z2021,76351 Linkenheim-Hochstetten,Germany
-----------	------------	---

UK	REP	Pooledas Group Ltd Unit 5 Albert Edward House, The Pavilions Preston, United Kingdom
-----------	------------	--

Made In China

VEVOR[®]

TOUGH TOOLS, HALF PRICE

Support and E-Warranty Certificate

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Importerad till Australien: SIHAO PTY LTD . 1 ROKEVA
Importerad till USA: Sanven Technology Ltd.