# Haier

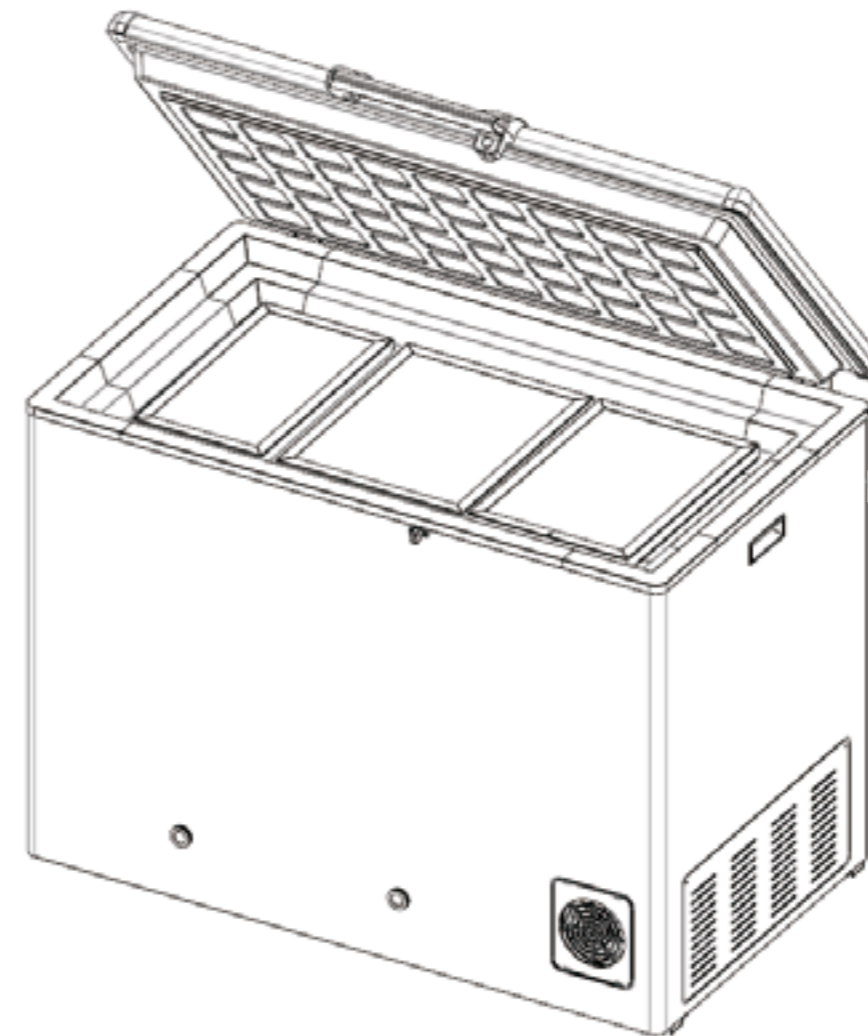
## Vaccine combined refrigerator&freezer Operation Manual

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
Français

El español

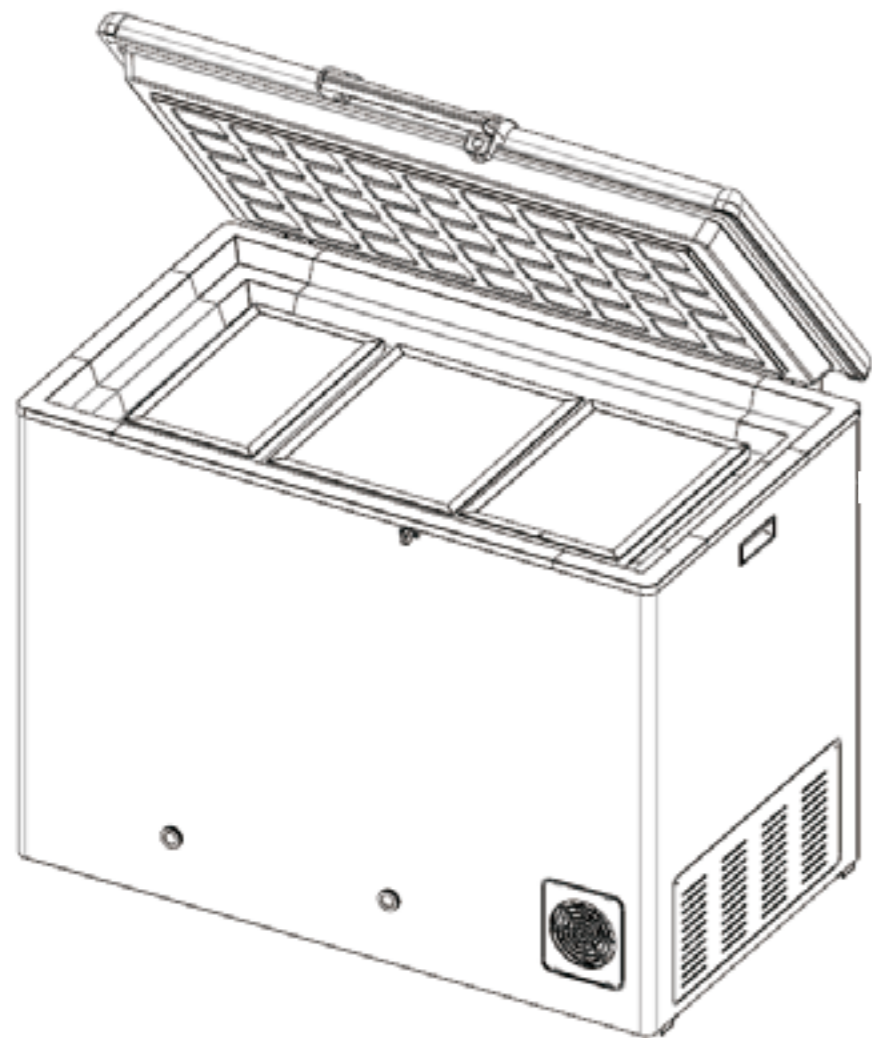
- Please read these instructions carefully before using this product.
- Our company reserves the right to interpret the instructions
- The appearance of the vaccine combined refrigerator&freezer shall be subject to the actual object.
- Please keep it properly after reading.
- In case of technology or software upgrade of the vaccine combined refrigerator&freezer, it is subject to change without further notice.

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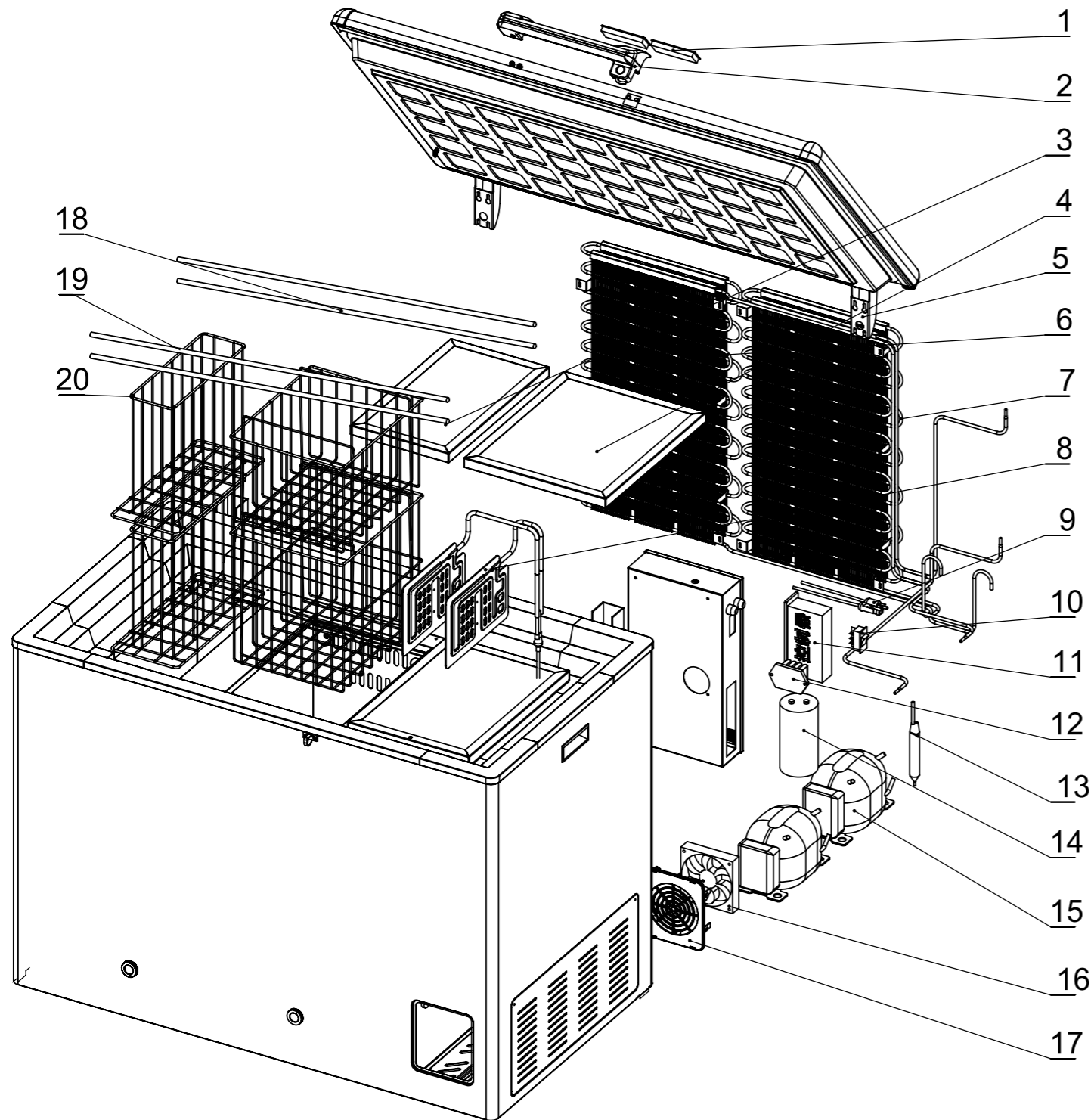# Haier quality, it merits your trust from beginning to end.

The vaccine combined refrigerator&freezer is applicable to pharmacy, pharmaceutical factory, epidemic prevention station, health center, hospital, etc. The refrigerator chamber is used to store biological products: vaccines that meet the storage conditions of + 2℃ ~ + 8℃ ; and the freezer is used for freezing and storing ice pack.

- Temperature Control**
  - Electronic thermostat control
  - Solar temperature displayer can use natural light source to drive display of inner temperature.
- Safety System**
  - Multiple fault alarms (high or low temperature alarm, power-off alarm, sensor fault alarm)
  - Two modes for alarms (sound buzzer alarm, light flashing alarm)
  - All independent components have safe grounding
  - Key components of high performance (compressor, fan, etc.) for security and reliability
- Excellent performance**
  - Authority certification of the International Health Organization (PQS certification)
  - It can meet the performance guarantee in the condition of only 8 hours of power supply per day.
  - Wide range ambient temperature design ensures that the performance is also outstanding at the ambient temperature of 5℃ ~ 43℃
  - Wide range voltage design ensures that the performance is also stable within the range of ±20% voltage.
- Ergonomic design**
  - Ergonomic design of computer control mode ensures intelligent control with no need to adjust.
  - High performance insulation layer contributes to a better thermal insulation effect
  - Safety door lock and other ergonomic designs

Haier vaccine combined refrigerator&freezer you got is probably not exactly the same as illustrated in the instructions due to upgrading of the product, we regret for this. The contents of the instructions are subject to change without prior notice.


Product Feature.....1  
 Contents .....2  
 Safety precautions.....3  
 Precautions for use.....6  
 Name of each component .....12  
 Technician’s manual .....13  
 Cleaning and Maintenance.....15  
 Wiring diagram .....17  
 Specification .....18  
 Package diagram.....19

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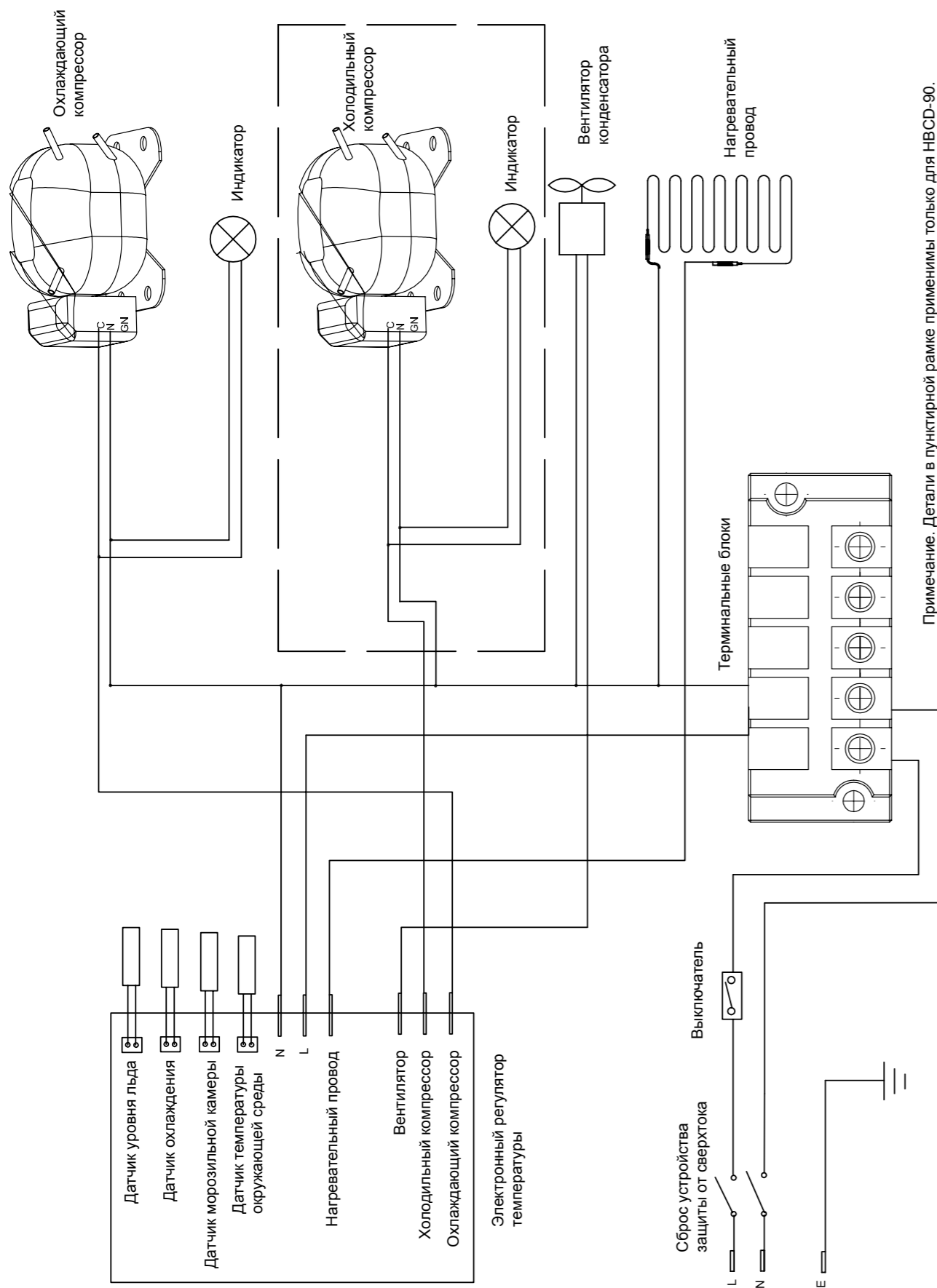Dear Haier users:

Thank you for using Haier vaccine combined refrigerator&freezer. In order to better read this manual and use the product to prevent personal injury and damage to the product, please read the instructions carefully and observe the contents marked with the following signs in the instructions.

 Actions or operations which are prohibited

 Actions or operations which must be followed

-  Close down the gas leaking valve and open the door and windows for ventilation in case of leakage of inflammable gas like coal gas etc., do not insert or pull out the power plug of the vaccine combined refrigerator&freezer, otherwise it may trigger an explosion and fire.
-  Never cut or remove the grounding pin of the power cord in any case. Make sure that the power plug and socket is inserted tightly and reliably, otherwise it may result in a fire.
-  Please use a power supply socket with a ground wire for fear of electric shock. If the power socket is not grounded, be sure to have a professional technician to install the grounding wire.
-  Do not use it in the open air. When the device is exposed to rain, it may cause electric leakage or electric shock.
-  Do not place the combined refrigerator&freezer in a damp site or a place where it may easily suffer from spraying water. Otherwise, it may cause leakage of electricity or electric shock due to degeneration of its insulation.
-  Do not pour water directly onto the vaccine combined refrigerator&freezer, otherwise it may cause an electric shock or a short circuit.
-  Do not place any vessel containing water or any heavy object on the combined refrigerator&freezer. Avoid power leakage or electric shock caused by the degeneration of insulation caused by water flowing out.
-  Do not make the vaccine combined refrigerator&freezer grounded via a gas-pipe, power supply conduit, telephone line or lightning rod. The above grounding will cause an electric shock or other dangers.
-  Do not contact any electric parts or any switches such as power plug with a wet hand, or otherwise it may cause an electric shock.
-  Neither set off hazardous articles that are inflammable, explosive or volatile in the vaccine combined refrigerator&freezer, nor use the flammable spraying agent near the vaccine combined refrigerator&freezer, or otherwise, it may trigger an explosion or fire.
-  Do not store corrosive articles such as acid, alkali etc. in the vaccine combined refrigerator&freezer. Otherwise, it may spoil the internal components or electric parts of the vaccine combined refrigerator&freezer.
-  Do not open holes on the vaccine combined refrigerator&freezer without permission, which would impact the refrigeration performance of the vaccine combined refrigerator&freezer.
-  It is not allowed to insert any article such as copper wire, iron wire or stick into any hole or gap of the vaccine combined refrigerator&freezer, otherwise, it may cause product damage or personal injury.

- ⊘ Do not climb up to the vaccine combined refrigerator&freezer or place articles on it, otherwise it may cause personal injury or product damage due to overturn of the vaccine combined refrigerator&freezer.
- ⊘ Do not place the plastic packing bags at the locations where children can easily take as plastic bags may cause occurrence of suffocation accident.
- ⊘ Children should be kept away from the combined refrigerator&freezer to prevent electric shock or other unknown injury.
- ⊘ When pulling out the plug from the power socket, hold the power plug firmly but do not pull the wire of the power plug. Pulling the wire with hand may cause an electric shock or a fire due to short circuit.
- ⊘ If the vaccine combined refrigerator&freezer runs out of order, please pull out the power plug. Continued running in an abnormal condition may cause an electric shock or fire.
- ⊘ Be sure to power off the vaccine combined refrigerator&freezer before you repair or maintain it for fear of electric shock or personal injury
- ⊘ Do not pull or carry the equipment by the door handle in case of damage to the combined refrigerator&freezer or personal injury.
- ⊘ Do not spoil the refrigerating circuit.
- ⊘ Do not use any electric appliance in the refrigerator of the vaccine combined refrigerator&freezer except for the types recommended by the manufacturer for use.
- ⊘ Do not place glass bottles or filling articles in the freezer, which may cause injury to personnel due to frost cracking.
- ⊘ Do not place the packing plastic bags at the locations where children can easily take as plastic bags may cause occurrence of suffocation accident.
- ⊘ Do not place the vaccine in the freezer.
- ⚠ If the vaccine combined refrigerator&freezer is left unused for a long time, disconnect the power supply to prevent electric shock, electric leakage or fire due to aging of power line.
- ⚠ If the vaccine combined refrigerator&freezer is left unused for a long time in an unattended area, make sure that children will not approach the combined refrigerator&freezer and the vaccine combined refrigerator&freezer door not be closed completely.
- ⚠ Only qualified personnel or after-sales service personnel can install this vaccine combined refrigerator&freezer, and the installation by non-professional personnel may lead to fire or damage.
- ⚠ The scrapping of the vaccine combined refrigerator&freezer shall be conducted by the relevant personnel. When scrapping, the door should be removed to prevent accidents such as suffocation.
- ⚠ Always place the vaccine combined refrigerator&freezer on a flat ground, as uneven ground may cause product damage or abnormal noise.
- ⚠ To store toxic, harmful or radioactive articles, please use this combined refrigerator&freezer in a safe place. Improper use may cause harm to human health or environment.
- ⚠ Make sure not to suck in medicine or suspended particles inside and surrounding the device at the time of repair and maintenance, otherwise, it may bring hurt to your health.
- ⚠ After moving the vaccine combined refrigerator&freezer, we should let it stand for 10 minutes, and do not power it on immediately, otherwise the refrigeration system may be damaged.

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- ! The door body handle should be held to close the door cover to avoid pinching fingers.
- ! Put on your gloves at the time of maintenance so as not to bump the sharp edges or corners, resulting in personal injury.
- ! The inclination angle shall be no greater than 45° when moving the vaccine combined refrigerator&freezer.
- ! Only the professional technicians or after-sale service men can install this refrigerator. Otherwise it may result in electric shock or fire.
- ! Make sure to place the vaccine combined refrigerator&freezer firmly on a solid and flat ground. If the ground is not stable or it is not properly placed, it will cause overturn of the combined refrigerator&freezer or personnel injury.
- ! Please use the dedicated power supply indicated on the nameplate of the vaccine combined refrigerator&freezer, or otherwise, it may cause a fire or electric shock.
- ! If the voltage is unstable, it is required to equip an automatic voltage regulator suitable for over 1000W motor load.
- ! If it is required to lengthen the power cord, the section area of the lengthen cord must not be less than 1mm<sup>2</sup> and the length shall not be longer than 3m. Otherwise, it may result in electric shock or fire.
- ! The power cord of this vaccine combined refrigerator&freezer is provided with a three-wire (grounding) plug that conforms to the standard 10A three-wire (grounding) socket.

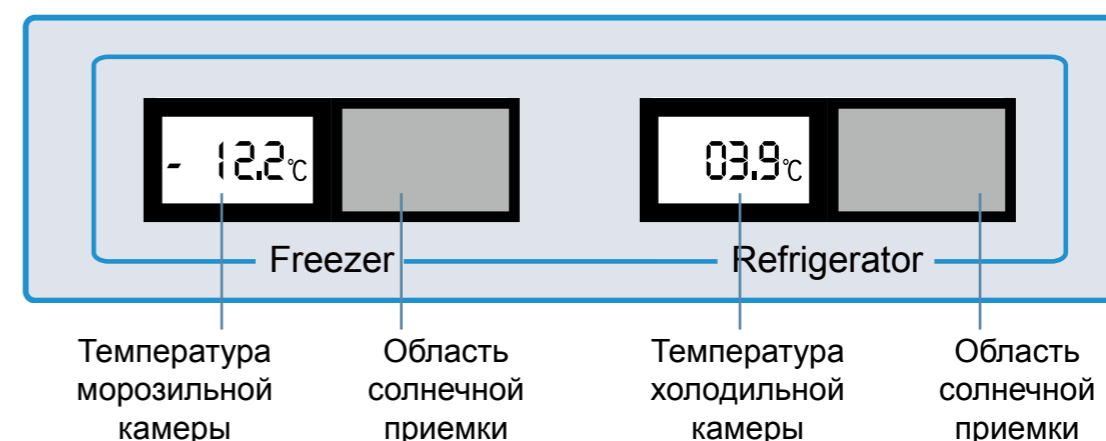
# Precautions for use

- Make sure firstly that the temperature inside the vaccine combined refrigerator&freezer has reached the setting temperature before putting articles into the refrigerator, and then put them in batches.
- The indicated temperature value of the vaccine combined refrigerator&freezer is the value indicated inside the box at the temperature sensing probe. Although the temperature displayed may sometimes be somewhat different from the actual temperature of the central portion of the vaccine combined refrigerator&freezer, it will gradually approach the real temperature.
- Use the diluted neutral detergent to clean the vaccine combined refrigerator&freezer. Do not use brushes, acid, gasoline, soap powder, polishing powder or hot water to clean the vaccine combined refrigerator&freezer as the above materials may damage the painting surface as well as the plastic and rubber parts and components. Take particular care not to use volatile solvent such as gasoline to wipe plastic and rubber parts and components.
- Cut off the power of the vaccine combined refrigerator&freezer when it is not in service for a long time.
- Every time when putting in and taking out the articles, please try to reduce the door opening time so as not to cause larger fluctuations of the temperature and humidity inside the refrigerator.
- After opening the door, the temperature inside the refrigerator will rise steeply in a short time, which is normal, and the temperature will gradually recover after the door is closed.
- The temperature range of this series of the vaccine combined refrigerator&freezer is +5°C ~ +43°C . If it is lower than or higher than this range, it may cause the temperature of the vaccine storage area in the box to exceed the range of +2 ~ +8°C .
- In the high-temperature and high-humidity environment, the product's cabinet may have condensation, which is a normal phenomenon without affecting the storage temperature in the refrigerator. If it happens, please improve the ventilation condition as soon as possible, and dry it with a clean rag.
- The vaccine combined refrigerator&freezer shall be equipped with a storage basket, and the articles shall be kept in the basket as far as possible.
- The indicated temperature of the vaccine combined refrigerator&freezer may have a certain difference with the actual temperature inside the box due to refrigeration inertia, which is normal.
- It is strictly prohibited to place the drugs such as vaccine which should have been in the refrigerator in the freezer.

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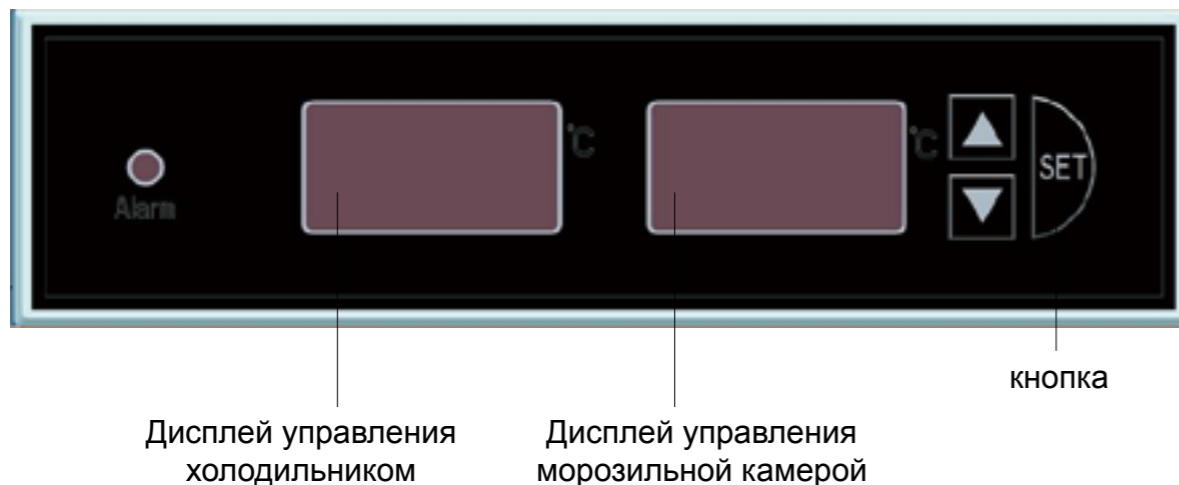
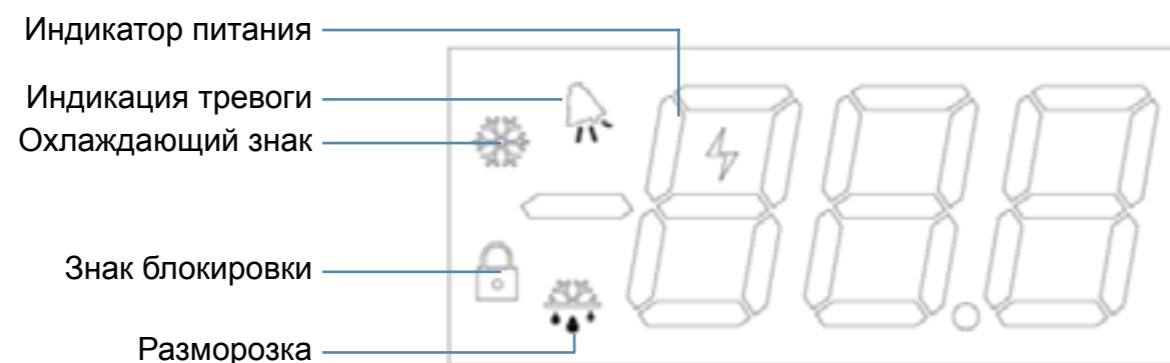
## Installation environment

- Ambient temperature: 5°C to 43°C , good ventilation, and air conditioning system shall be used if necessary.
- Ambient humidity: 15%~90%.
- Avoid large amounts of dust.
- Avoid mechanical sway or vibration.
- Rated power: 220V-240V~/50Hz , referring to each model specification list for detail.

Caution

- It is prohibited to install the vaccine combined refrigerator&freezer in the outdoor environment. When the vaccine combined refrigerator&freezer gets wet by rainwater, it may cause leakage of electricity or electric shock accident.
- The ambient temperature of the equipment shall not be lower than + 5°C , and shall not be higher than + 43°C .

## Installation site

For normal operation and optimum performance of the vaccine combined refrigerator&freezer, where it is installed shall conform to the following conditions:

- It can only be installed alone and shall not be installed in a recess or installed in an embedded manner, and not in a narrow, small and airtight space. The door frame of the room for placing it shall not be less than and lower than this product's, and shall at least guarantee its normal passing in and out so as to avoid any maintenance difficulty and failure of in-time repair of the machine, causing spoilage to the stored articles.
- It shall be installed on the solid and flat ground.
- It shall be well-ventilated without direct sunshine.
- It shall not share a socket with other any equipment, and its plug shall be well connected with the socket.
- Do not distort or press the power cord.
- If it is required to lengthen the power cord, the section area of the lengthened wire shall not be less than 1mm<sup>2</sup> and the length shall not be longer than 3m.
- Check the operating voltage before use, and the use of a voltage regulator suitable for the motor load shall be considered for voltage stabilization at the area where the voltage is not stable. So that the normal input voltage is stabilized at 220V±10%, and the power of the voltage regulator is more than 1000W.
- Be sure that the vaccine combined refrigerator&freezer is grounded reliably, and it is prohibited to cut or remove the third (grounding) plug in the power line in any case.
- If the power socket has a grounding wire, please check whether it is good before use.
- If the socket is not installed with a grounding wire, please be sure to have a professional engineering and technical personnel to install the grounding wire.
- Every 8g of R600a refrigerating fluid shall have a space of ≥ 1m<sup>3</sup> to avoid the formation of combustible mixed gas in case of accidental leakage of it.

Warning

- No ground connection for the vaccine combined refrigerator&freezer via gas piping, water supply pipe, telephone line or lightning rod that may easily cause an electric shock.
- After installed, the power bus or power plug must be accessible in order to disconnect power supply in case of emergency. Do not block the ventilation of the vaccine combined refrigerator&freezer.

## Preparation before use

### 1. Take down the packing materials and the packing straps

Take down all the packing materials and packing straps used for transportation.

- The vaccine combined refrigerator&freezer can be loaded and unloaded with a forklift or a package-clamping forklift before it is unpacked.
- If a forklift is used, please insert it from the bottom of the wooden bottom bracket on the front or back of the vaccine combined refrigerator&freezer and then move the vaccine combined refrigerator&freezer .
- If a package-clamping forklift is used, please clamp it from the bottom of the wooden bottom bracket, and only the side of the vaccine combined refrigerator&freezer can be clamped.
- After the package is removed, please strengthen the protection to avoid hitting your foot when carrying it or changing its position.
- Be sure that the tilt angle of the vaccine combined refrigerator&freezer no greater than 45° when it is carried.
- Packaging plastic bags shall be kept away from children.

### 2. Check the accessories

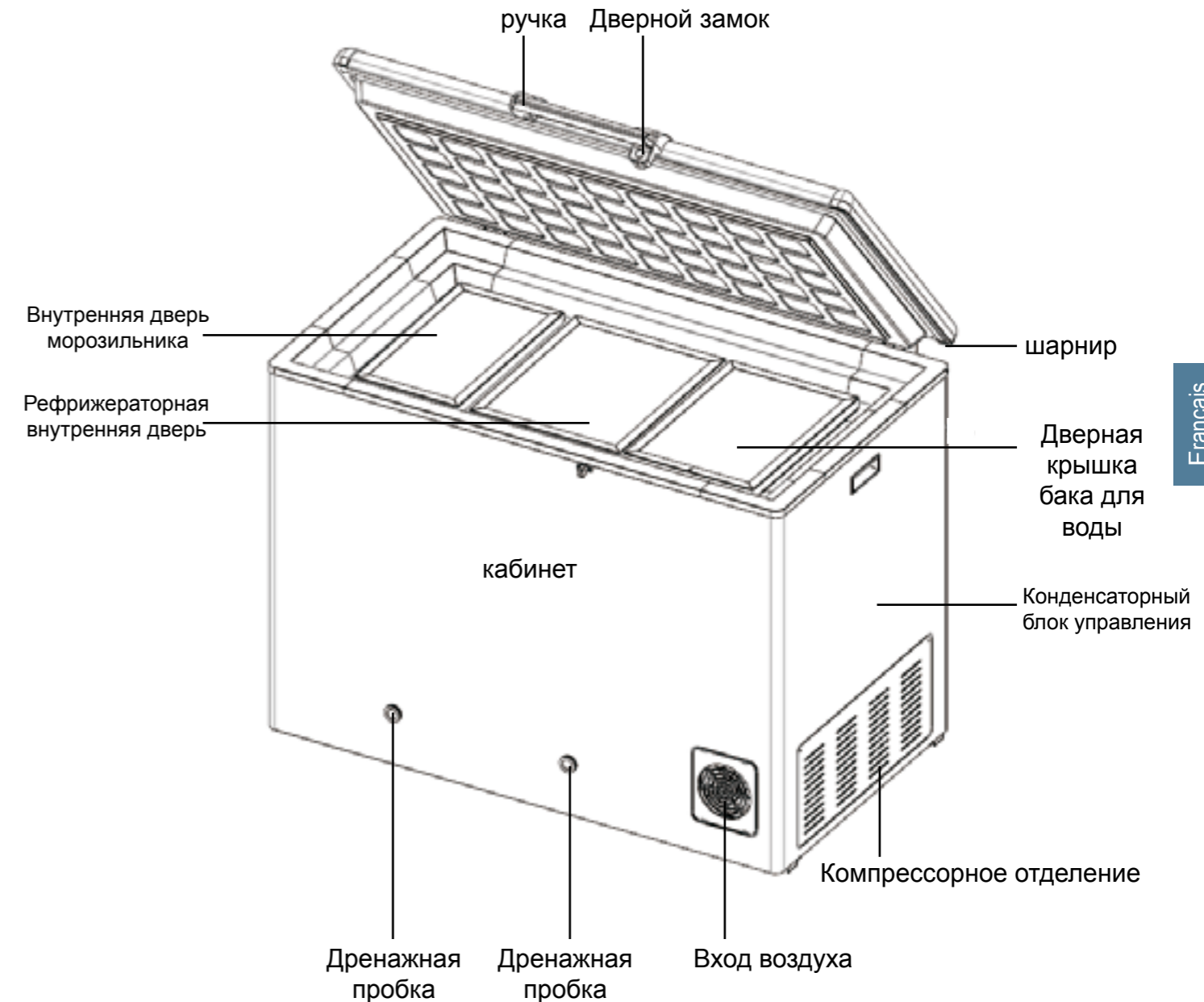
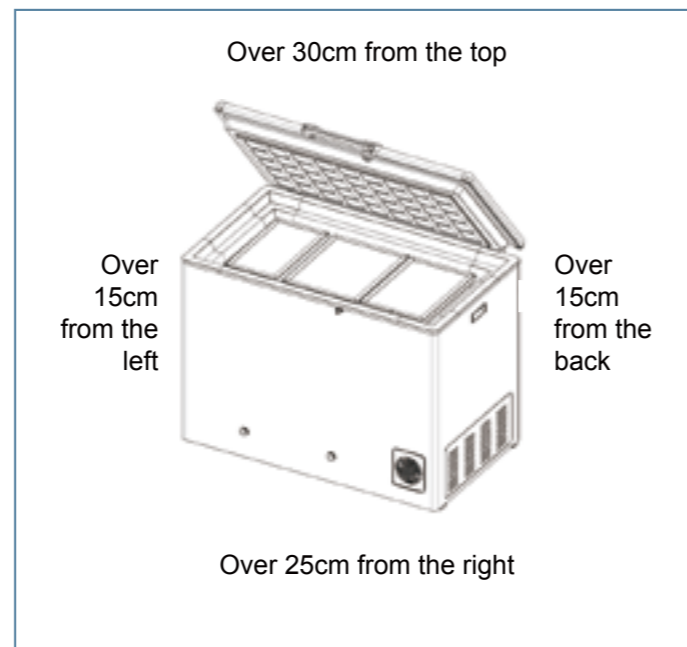
Please check the articles in the vaccine combined refrigerator&freezer according to the packing list. If there is any discrepancy, please contact the after-sales service in time.

### 3. Placement condition

- Leave a space of at least 15cm around the vaccine combined refrigerator&freezer to ensure good circulation of surrounding air and facilitate heat dispersion.
- In order to ensure the normal opening of the door, a space of at least 80cm shall be reserved above the vaccine combined refrigerator&freezer.

sufficient space (at least 25cm away from the wall surface) shall be reserved at the engine-room grating to facilitate maintenance and ventilation.

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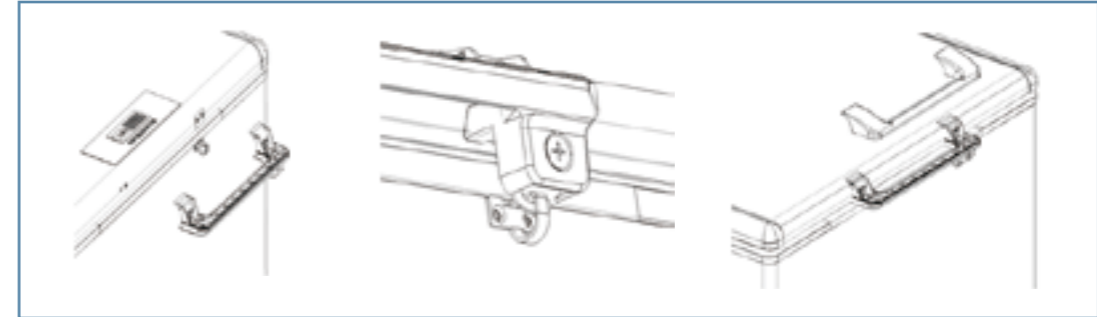
## 4. Handle installation

- Take out the handle and remove the four bolts from the door
- Align the lower part with the screw hole on the door cover
- Make sure that the locking hooks on the lock nose and machine body can be accurately matched
- Tighten screws in turn.
- Buckle the handle cover and tighten the bottom bolts.



## 5. Standing


Do not power on the device firstly before it stands stably and is cleaned. Let it stand for over 24 hours before power it on so as to ensure normal running of the vaccine combined refrigerator&freezer.

## 6. Add water to the water tank

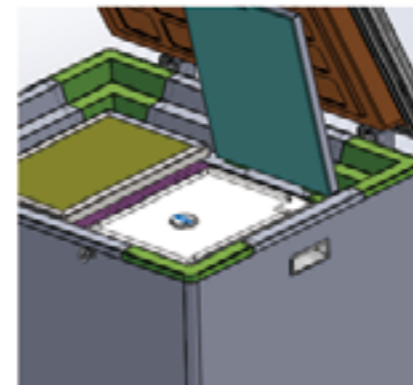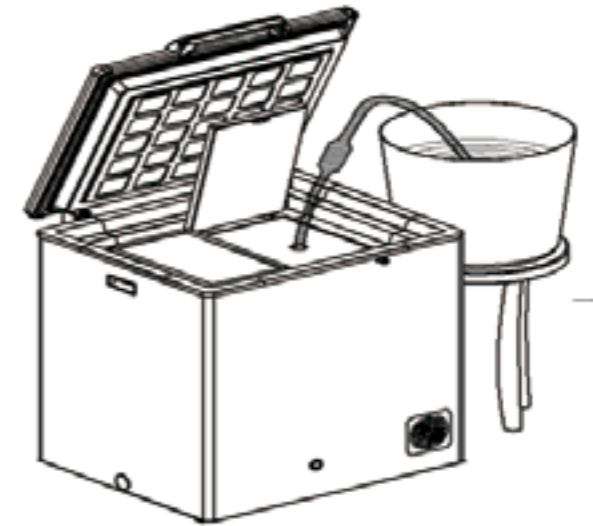- Open the water tank cover on the right side of the box and unscrew the cover bolts.
- Pour the whole bottle of corrosion inhibitor into the water tank first;
- Put a pure water container with a certain volume (about 25L) of water beside the combined refrigerator&freezer and higher than the water tank
- Insert the water inlet of the suction device into the water container, and insert the water outlet into the water tank. If the air bag of the suction device is extruded continuously by hand, the water flows into the water tank from the container. After the water begins to flow, it is unnecessary to squeeze the air bag, and the water will automatically flow into the water tank;

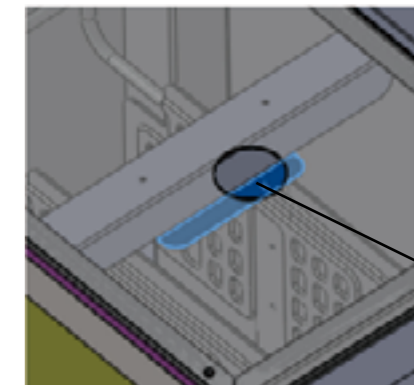a. Open the right cover of the water tank of the cabinet.

b. Stop water injection when the liquid level is even with this level.

- 5). After injecting a certain volume of water, observe the level of the liquid. If the level is even with the liquid level indicated in the diagram, stop water injection;
- 6). If the injection is completed, please screw in the water filling nozzle plug.
- 7). Cover the top cap and tighten the screws.

#### 7. Main power switch

Please confirm the followings before powering it on:

- a. Verify whether the voltage is 220 V-240V and 50Hz alternating current;
- b. Be sure to let the vaccine combined refrigerator&freezer stand for more than 24 hours at the first time of electrifying.
- c. Open the power switch on the side of the electric control cabinet after confirming no errors above.

#### 8. Device operation

Check whether the compressor and fan in the engine room are operating normally during power-on operation. The cooling time of each vaccine combined refrigerator&freezer is different due to different operating environment.

#### 9. Conditions for placing drugs

- a. Continuous electricity supplying for 48 hours; temperature in the refrigerator has been stable within the range of (+ 2°C -+ 8°C ) for 24 hours;
- b. In case of intermittent electricity supplying, ensure that the total electrifying time is over 60 hours, and the temperature in the refrigerator has been stable within the range of (+ 2°C -+ 8°C )for 24 hours; The vaccine can be loaded only after the above conditions are met, and the vaccine should be placed in the vaccine basket as far as possible.

Pay attention to observe the temperature change of the freezer's display screen. When the temperature is  $\leq -5^{\circ}\text{C}$  , the freezer can be loaded with ice pack.

- Do not connect it with the power source that does not conform to the provision.
- The display on the door cover is driven by solar energy without other electricity supplying. There may be no temperature displaying or screen flicker in case of weak light. If you use a flashlight or other light to light it for about 3 seconds, the screen will display the real-time temperature inside. And the screen will return to the original condition without light source.

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## Initial Power-on

Please observe the following rules when conducting initial start-up and continuous running:

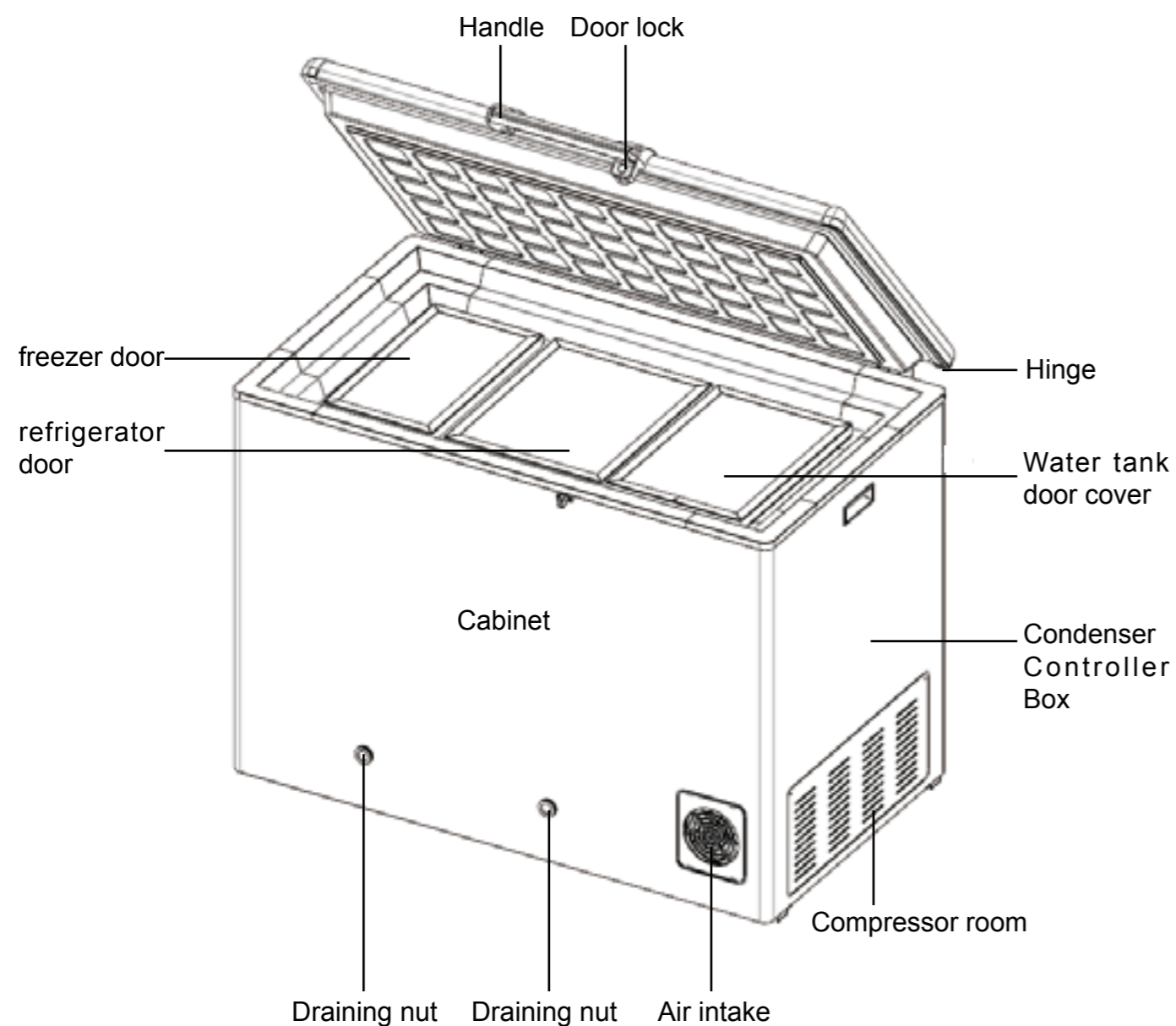
When the vaccine combined refrigerator&freezer is empty, connect it to the power source of the appropriate specification (220V-240V~, 50Hz).

After it is powered on, the temperature control display will be illuminated.

The corresponding parameters have been set before the vaccine combined refrigerator&freezer leaves the factory, and there's no need for self-setting.

After continuous electrifying for 48 hours or intermittent electrifying (total electrifying time is 60 hours), the temperature inside the refrigerator is constant between + 2°C and + 8°C .

After checking the running of the vaccine combined refrigerator&freezer, the temperature shall be reduced to a reasonable range, and the articles can be loaded in.

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
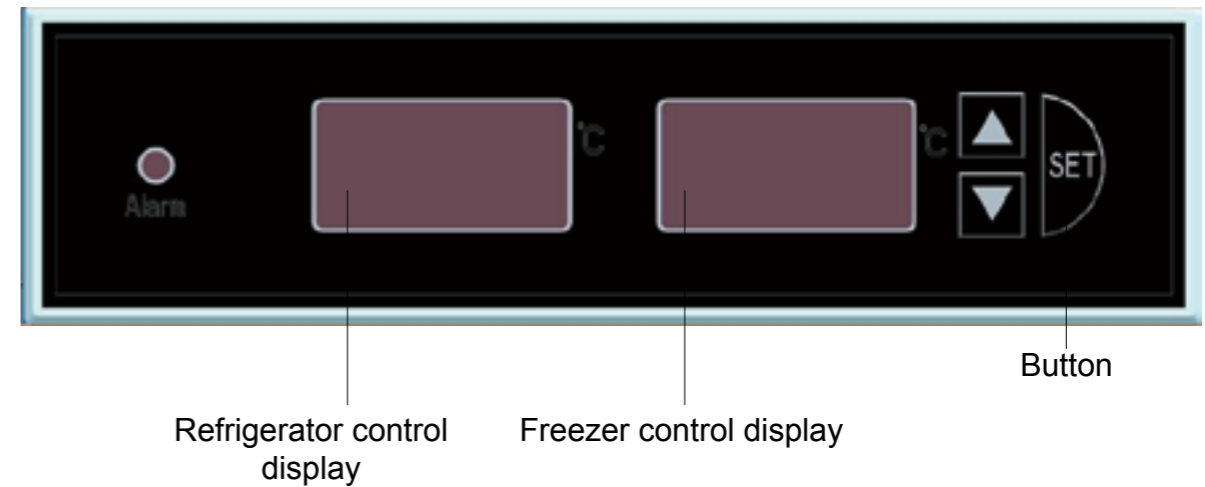
## Thermostat

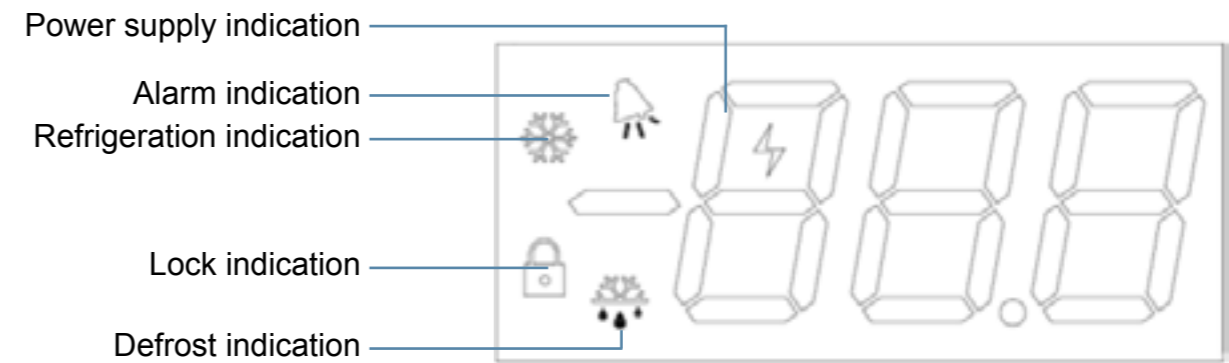The Controller installed in the electric cabinet is used for controlling compressor, fan and heater strip. The parameters have been set when the machine is delivered and does not support user's change.

1. Power source mark: This symbol is displayed when the external voltage reaches a certain value.
2. Warning sign: When the Controller is giving an alarm, this symbol flickers to remind the user.
3. Refrigeration mark: This symbol is displayed when refrigeration is output.
4. Lock sign: This symbol is displayed when the controller is locked.
5. Defrost sign: This symbol is displayed when the controller is defrosting. This product does not have this function.

## Caution

For the HBCD-90 vaccine combined refrigerator&freezer, the ice storage temperature of the water tank in the refrigerator is displayed on the right screen, and the temperature in the freezer shall be displayed on the left screen.

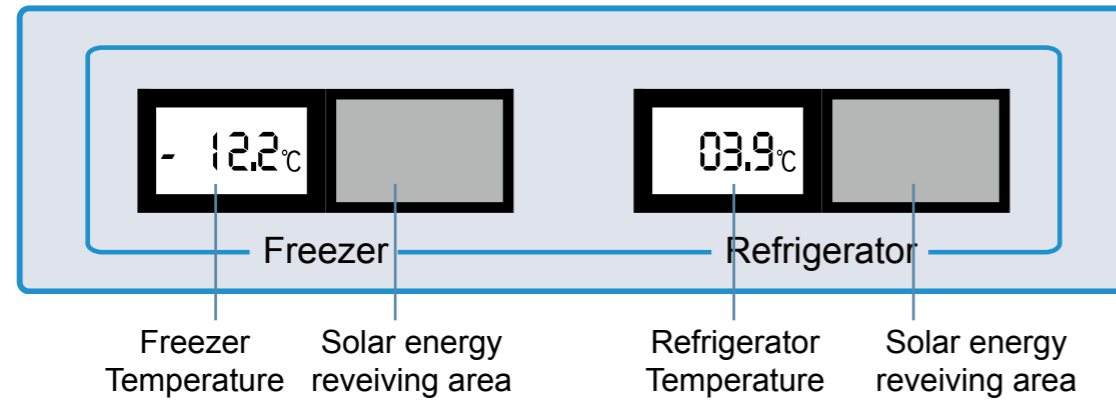
When the temperature controller is powered on, the temperature controller screen is illuminated, and the buzzer sounds twice; if there is no operation of any button within 2 minutes, the controller will automatically turn off the display, and only the power source mark flickers.

Press the \* SET \* button again, and the screen will turn on again.

Alarm function introduction (for after-sales service reference only)

Indication code	Code Descriptions	Indication code	Code Descriptions
Er1	Ice storage sensor failure	AL1	Ice storage low temperature alarm
ER2	Refrigerating sensor failure	AH2	Refrigerating high-temperature alarm
Er3	Freezing sensor failure	AL2	Refrigerating low temperature alarm
Er4	Ambient temperature sensor failure	AH3	Freezing high temperature alarm
AH1	Ice storage high temperature alarm	AL3	Freezing low temperature alarm

## Temperature indicator

REFERERATOR---Show the temperature of the vaccine storage area;

FREELESER---Display the temperature of the freezer;

The displayer is powered by solar energy and requires no extra connection to electricity. The displayer only displays the temperature inside the box, and does not control the refrigeration system of the product.

When it is used, the solar energy receiving area cannot be shaded, otherwise, the temperature display area will not work.

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## Component Cleaning

Warning

- Turn off the master switch and disconnect power before cleaning.
- The detergents for cleaning the product must be diluted, otherwise the surface of the plastic parts or the housing will be discolored.
- If the product will not be used for a long time, turn off the master switch and disconnect power, and the articles in the box shall be emptied.

### Daily maintenance

- Wipe the inner container all around to prevent leakage of drugs or corrosion by other liquids;
- When the container of the inner container or the ice pack freezing bracket is frosted by more than 5 mm, it shall be cleaned in time to avoid excessive frost, which impacts refrigerating performance.

### Monthly maintenance

- Remove dust for engine-room grating, condensate fan and compressor baseboard to prevent negative influence on cooling effect.
- Wipe the door seal to prevent dirt impact on the door seal.
- Wipe off the dust from the product's housing and internal space with a dry cloth. If the product is very dirty, use a piece of cloth dipped with neutral detergent to clean off the dirt, and use a piece of wet cloth to wipe off the residual detergent, and then wipe it with a piece of dry cloth.
- Clean the condenser to prevent the impact of dust accumulation on heat dissipation.

### Quarterly maintenance

Check the water level of the water tank. When it is lower than the water line, proper water supplement is necessary.

## Scrapping of vaccine combined refrigerator&freezer

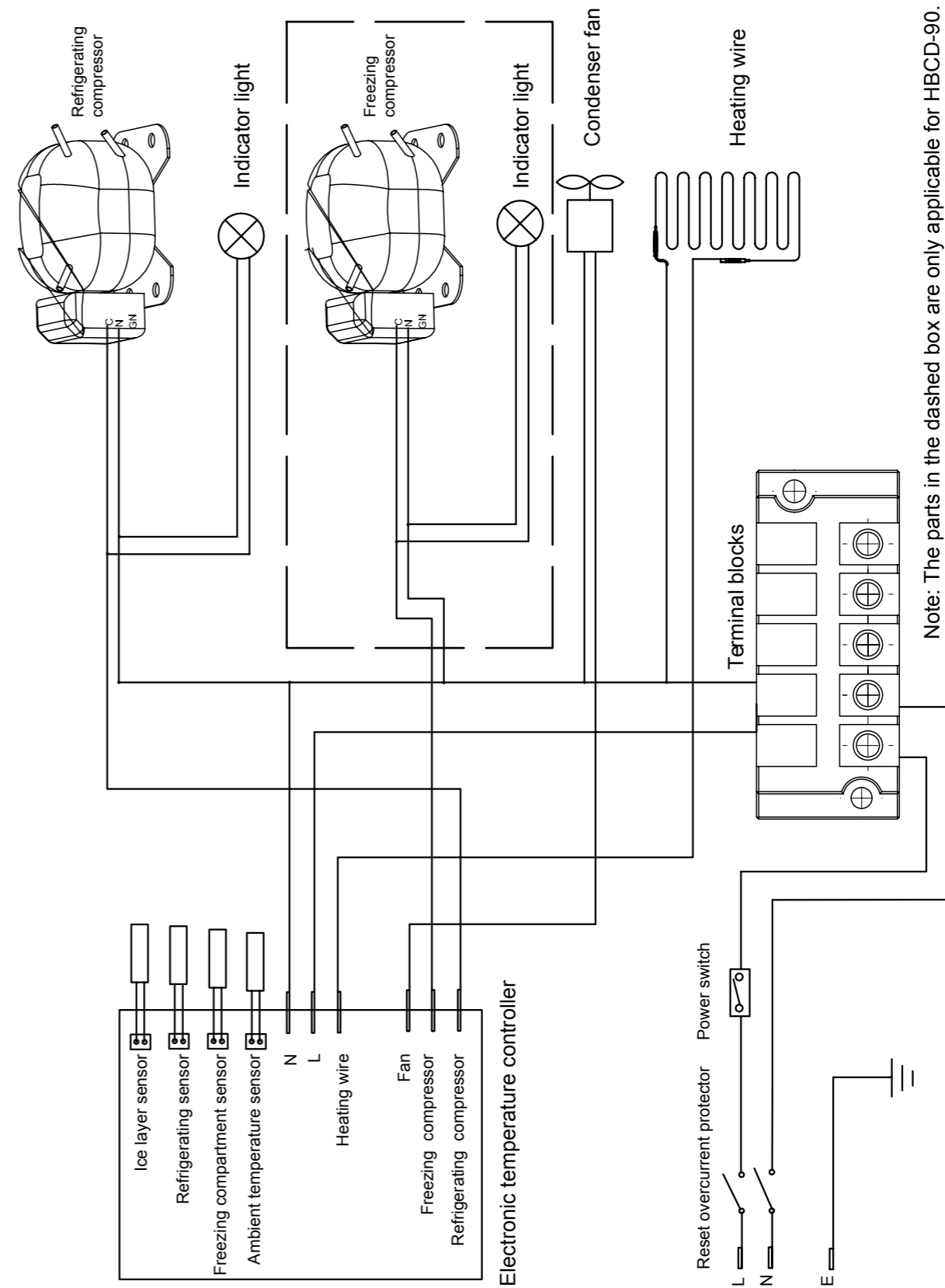
 There is a symbol on the vaccine combined refrigerator&freezer, its package and its accompanying documents, which indicates that the used electrical, electronic product and battery cannot be equated with the general household garbage. In order to properly handle, recover and reuse the old product, please follow the relevant regulations of the host country and the 2002/ 96/ EC and 2006/ 66/ EC to send it to the designated recovery place. Proper Treatment of these articles is to preserve useful resources to prevent them from causing damage to human health and environment, and to avoid evil consequences caused by improper treatment of the waste products. For more information about the recovery and re-use of articles, please contact the local government authority, the waste grounding center, or the sites for selling the product. According to the national regulations, inappropriate disposal of these wastes may be subject to disciplinary action.

For the EU business users, if it's needed to dismantle the electrical and electronic device, please contact the wholesaler or the supplier for further information. Information on disposal of these wastes in other countries other than EU is valid only in EU. If it's necessary to dismantle the vaccine combined refrigerator&freezer, please contact the local government authority or distributor for the correct method of disposal.

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Note: The parts in the dashed box are only applicable for HBCD-90.

## Specification

Model	HBCD-90
Surrounding temperature	+5 - +43
Temperature range of refrigerator (°C)	+2 - +8
Volume of refrigerator (L)	42
Volume of freezer (L)	32
Refrigerator compressor refrigerant R600a (g)	45
Freezer compressor refrigerant R600a (g)	40
Rated Voltage (V)	220V/240V
Rated frequency (Hz)	50
Rated power (W)	190
Dimensions W x D x H (mm)	1128*654*875

## Packing List

Model	The description	Keys	Handle attachment	Refiller	Corrosion inhibitors	Defroster spade
HBCD-90	1	2	1	1	1	1

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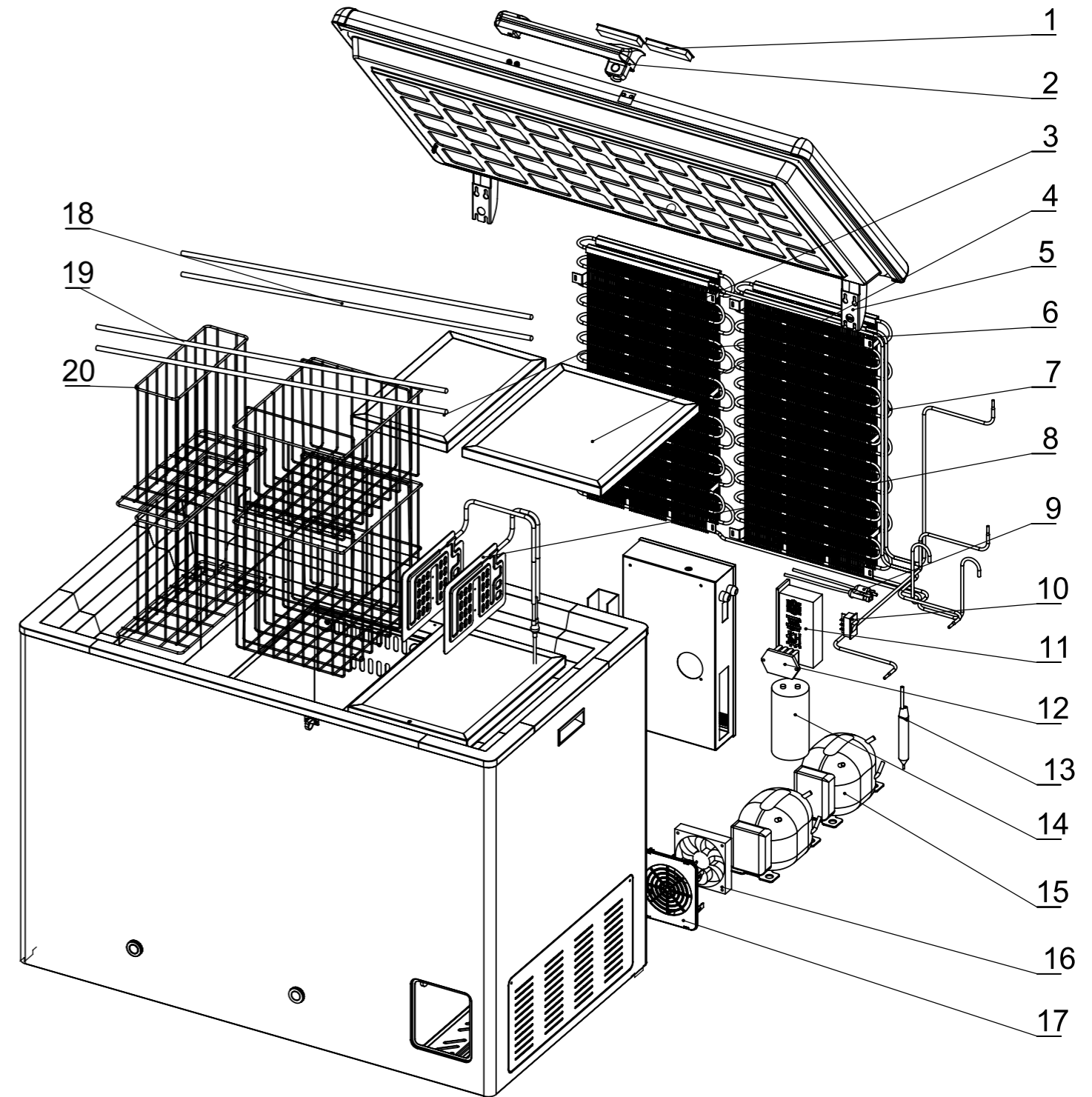
Haier quality,

it merits your trust from beginning to end.

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