

Cannabis Testing Solutions

Terpene Analysis by HS-GCMS

Dr. Manuel Braun
Shimadzu Europa GmbH



Terpenes

- Terpenes – organic, fragrant compounds produced by plants
 - Distinctive flavor and aroma
 - Primary constitute of essential oils of medical plants
 - Medicinal hydrocarbon building blocks
- Terpenes in cannabis
 - Terpenes are produced in trichomes (like THC)



- Synergistic effect with cannabinoids





GI Grow Biomedical Farm Video, available at Shimadzu's www.GrowYourLab.com

The Synergistic Effect

- Cannabis has over 140 terpene components, many of which are of medicinal interest



“Pine needles”

- Anti-inflammatory
- Enhanced concentration
- Bronchodilator for better cannabinoid absorption



“Lavender”

- Anti-anxiety
- Analgesic
- Anticarcinogenic





“Citrus”

- Anti-bacterial
- Anticarcinogenic
- Anti-depression


The Synergistic Effect

- “Entourage Effect”
- Terpenes have a synergistic relationship with cannabinoids, which further enhance the therapeutic effect
- The various compounds in cannabis work together to produce a **synergy** of effects

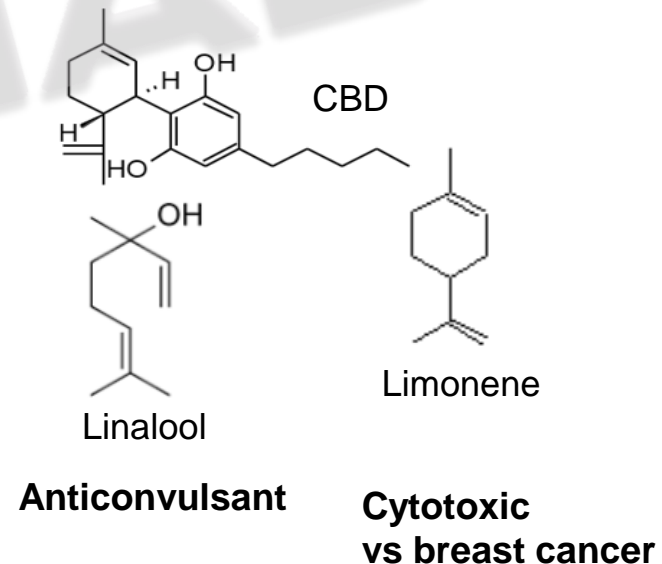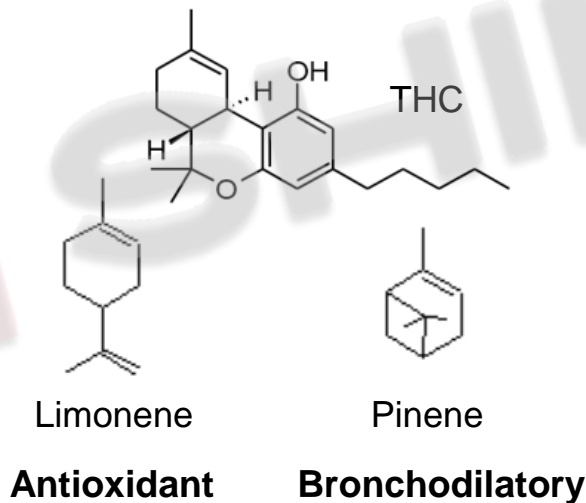

Terpene Profiling by GCMS

- Different strains of cannabis plants can be distinguished by different terpene contents
 - Due to the uniqueness of terpene profiles, they can be used by cultivators as a “fingerprint” to partially ID the specific strain in question
- Sativa vs. Indica
 - Typical Sativa Terpenes: α -Phellandrene and α -Terpinolene
 - Typical Indica Terpenes: myrcene, camphene and limonene

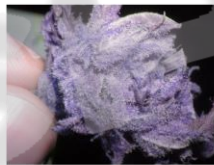

Purple DanX



Purple Erkle



Super Silver Haze



Trainwreck



Bubba Kush



OG Kush




Purple Kush



Sour Diesel

Terpene Profiling by GCMS

- Isoprene rule - General formula of terpenes is $(C_5H_8)_n$

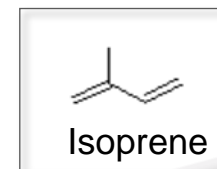

Monoterpenes (n = 2)

Sesquiterpenes (n = 3)


acyclic

monocyclic


bicyclic

acyclic

monocyclic

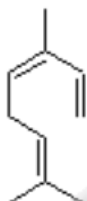


Ocimene

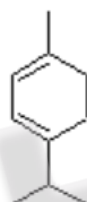


α -Terpinene



α -Pinene

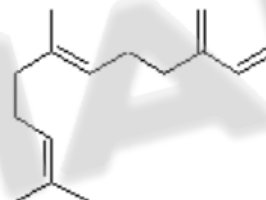


β -Farnesene

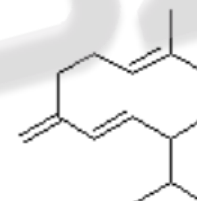


Germacrene D



Myrcene

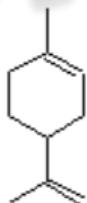


Limonene

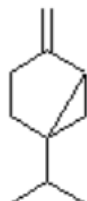


Sabinene

bicyclic

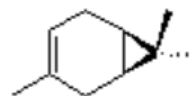


Carene

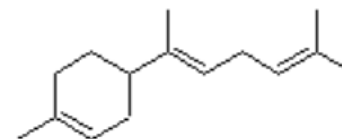

α -Bisabolene

Terpene Profiling by GCMS

- Gas chromatography is the preferred way to test for terpenes
 - Terpenes are volatile which makes them ideal candidates for GC
 - Boiling point between 119 – 200 °C and therefore easily converted to the gas phase
- Liquid chromatography
 - Limited chromophore
 - Low UV sensitivity
 - Coelution of terpenes and cannabinoids

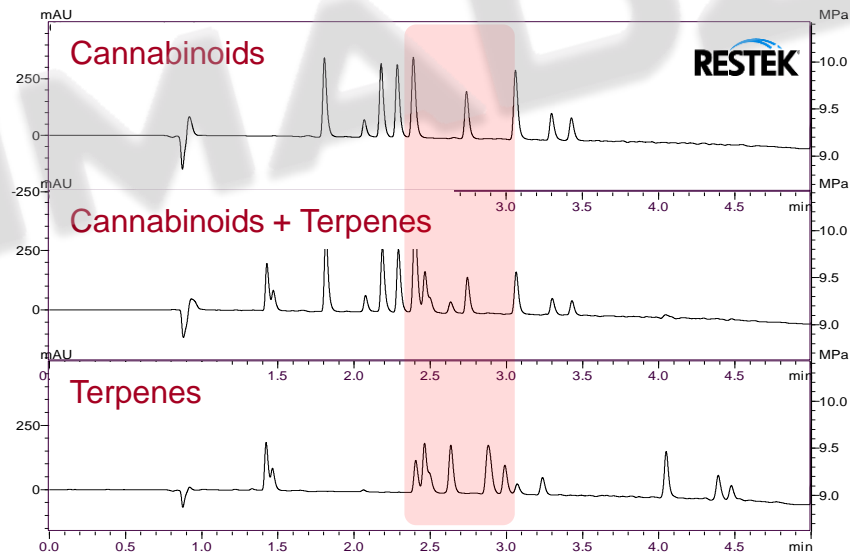

Headspace-Analysis with GCMS

- Mass Spectrometry (MS) or Flame Ionization Detector (FID)

MS

- Mass spectrum library for structural identification
- Can identify unknown peaks or contaminants


FID

- No library information
- Increased in linear range
- Good if you are only quantifying the same terpenes




Shimadzu GCMS-QP2020NX
with HS20NX (headspace) autosampler
for the analysis of terpenes

Headspace-Analysis with GCMS


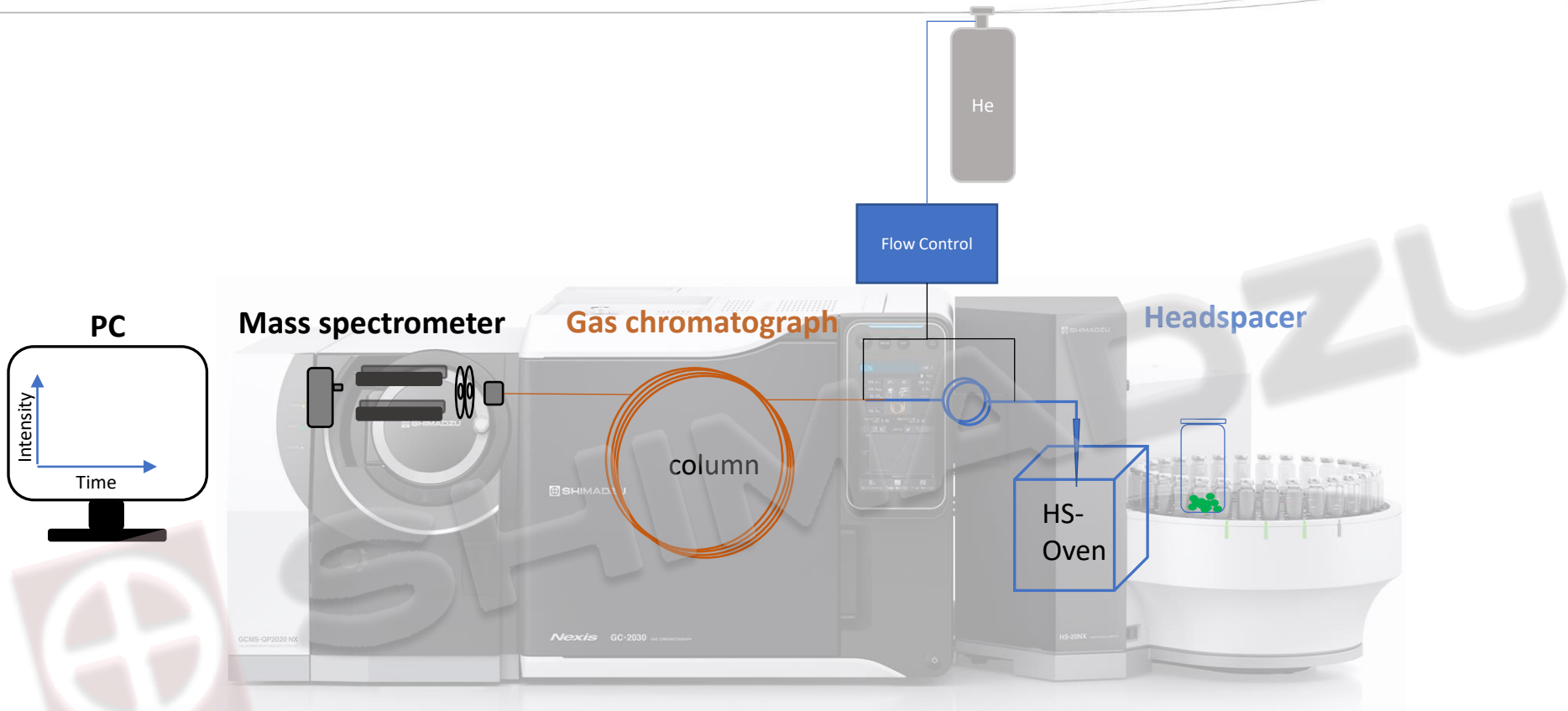


Headspace-Analysis with GCMS

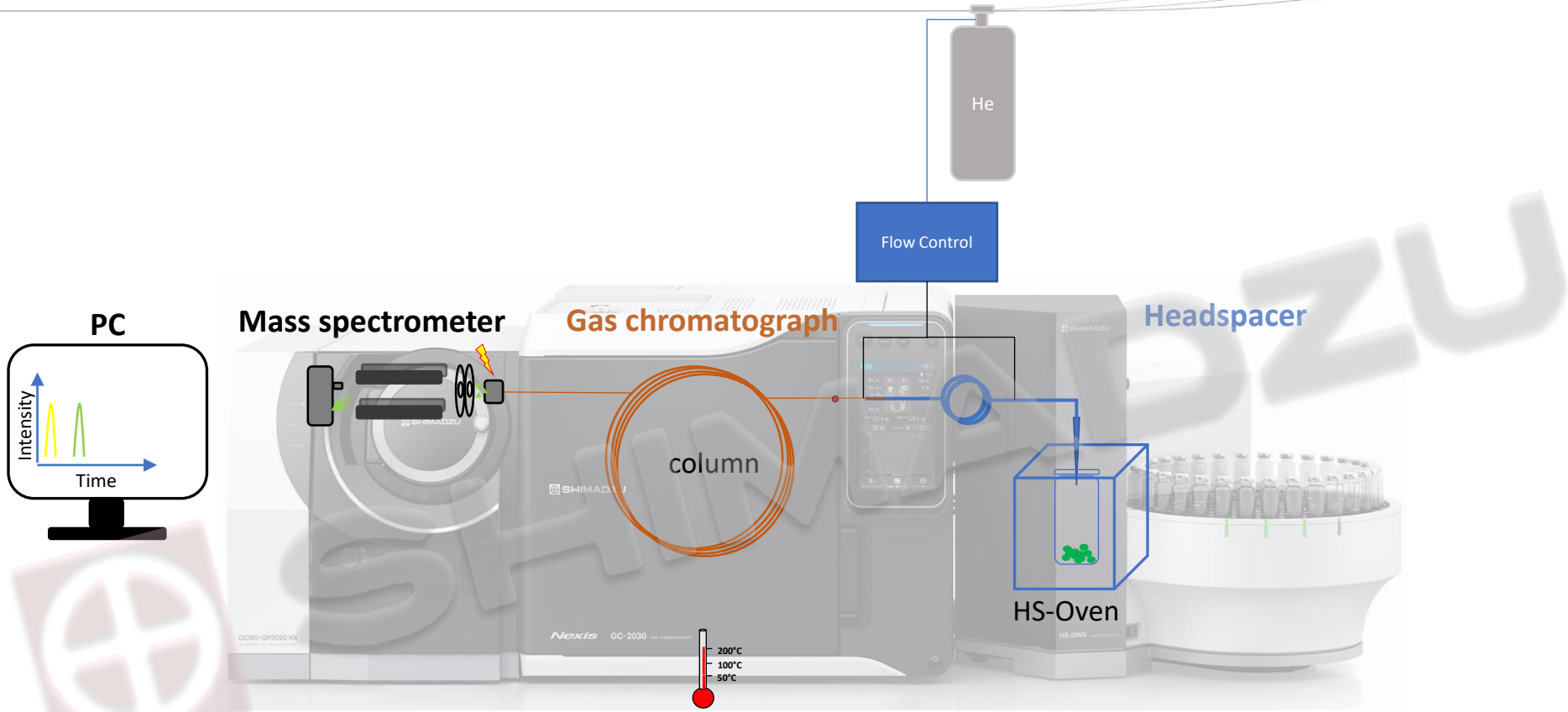

Terpene Profiling by GCMS

- HS is the sampling technique used to extract volatile compounds from a liquid or solid sample

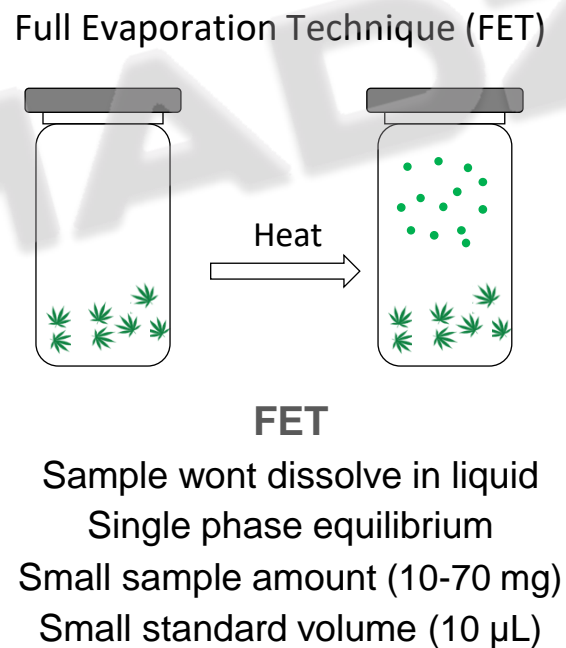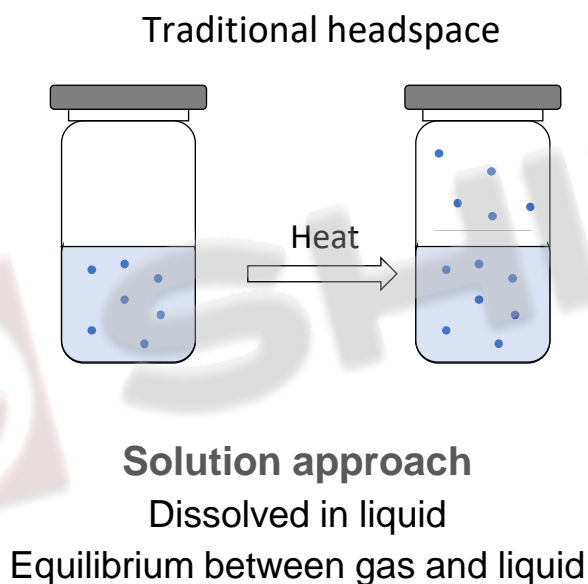


Terpene Profiling by GCMS

- Standard samples

Dilute standard solution



10 μ L volume

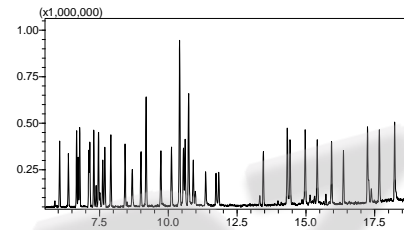

Analyze by GCMS

- Cannabis samples




Weigh out cannabis sample



Crushing



Weigh 10 - 70 mg in HS vial

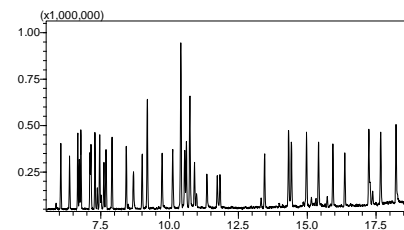

Analyze by GCMS

Hemp tea measured by HS-GCMS



Crushing


- At RT (ZM300)
- At -100°C (MM400+MM500)
- At -196°C (CryoMill)





Weighing



Analyze by GCMS

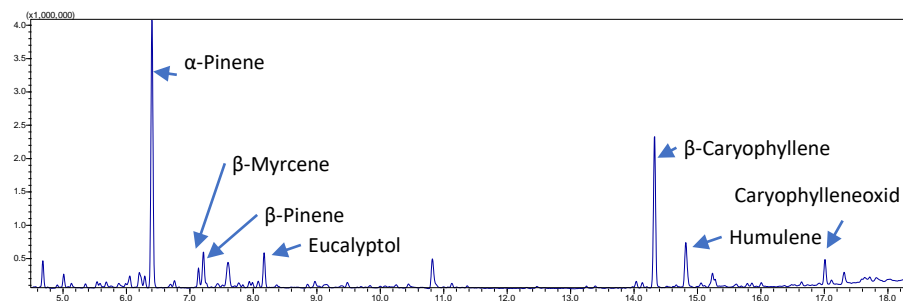

Hemp tea measured by HS-GCMS

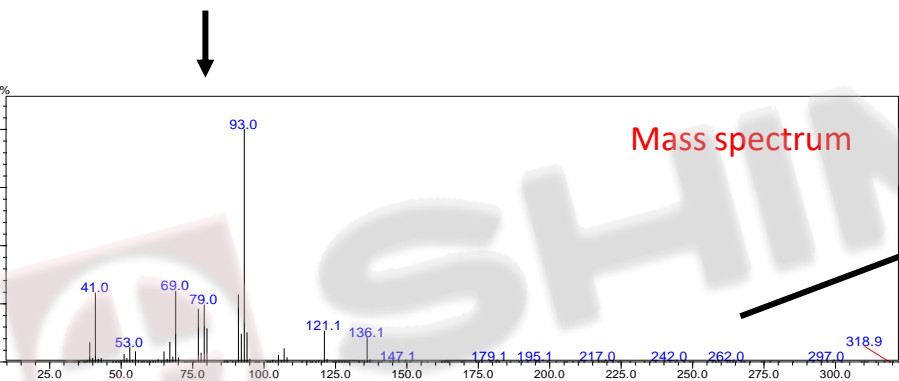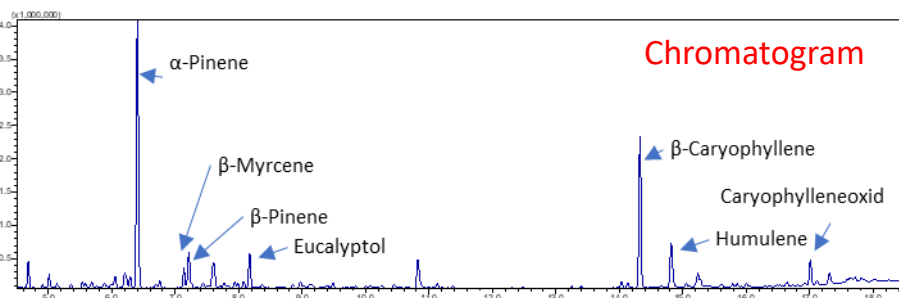


Similarity Search Results


Report View Compound Info Process Help

Hit#	Similarity	Register	Compound Name	Mol Wt	Formula	Library
1	94	<input checked="" type="checkbox"/>	Pinene <beta-> \$\$ Bicyclo[3.1.1]heptane, 6,6-	136	C10H16	FFNSC 4.0.lib
2	92	<input type="checkbox"/>	Sabinene \$\$ Bicyclo[3.1.0]hexane, 4-methylen	136	C10H16	FFNSC 4.0.lib
3	91	<input type="checkbox"/>	Mentha-1(7),8-diene <para-> \$\$ Cyclohexane,	136	C10H16	FFNSC 4.0.lib
4	90	<input type="checkbox"/>	Phellandrene <beta-> \$\$ Cyclohexene, 3-meth	136	C10H16	FFNSC 4.0.lib
5	89	<input type="checkbox"/>	Pinene <alpha-> \$\$ Bicyclo[3.1.1]hept-2-ene,	136	C10H16	FFNSC 4.0.lib
6	89	<input type="checkbox"/>	Apopinene <3-methyl-> \$\$ Bicyclo[3.1.1]hept-	136	C10H16	FFNSC 4.0.lib
7	88	<input type="checkbox"/>	Carene <delta-3-> \$\$ Bicyclo[4.1.0]hept-3-ene	136	C10H16	FFNSC 4.0.lib
8	86	<input type="checkbox"/>	Tricyclene \$\$ Tricyclo[2.2.1.02,6]heptane, 1,7	136	C10H16	FFNSC 4.0.lib

Target:

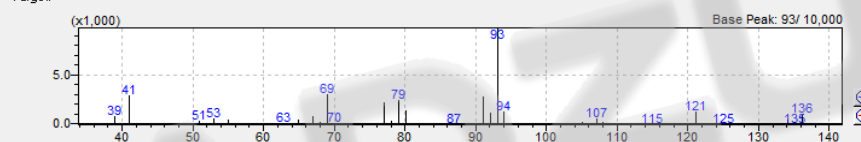

1 : 136 : Pinene <beta-> \$\$ Bicyclo[3.1.1]heptane, 6,6-dimethyl-2-methylene-

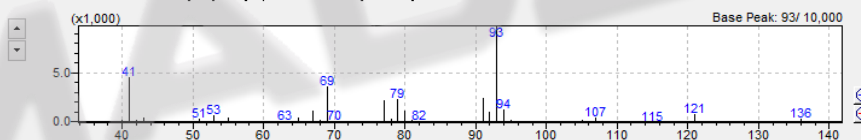

CAS#: 127-91-3 Mol Wt: 136 Serial#: 837
 Cmpd Name: Pinene <beta-> \$\$ Bicyclo[3.1.1]heptane, 6,6-dimethyl-2-methylene-
 Formula: C10H16 Class Flag: No Class Flags.
 Description: SAFC Cat. n.W290300 Column: SLB-5ms part#28471-U:Supelcowax-10 part#24079:Equity-1 part#28046-U; www.sigmaaldrich.com |
 Ret.Index: 978 / SLB-5MS(Hydro) T:0.25um L:30.0m D:0.25mm
 651 / SLB-5MS(FAMES) T:0.25um L:30.0m D:0.25mm
 505 / Supelcowax10(FAMES) T:0.25um L:30.0m D:0.25mm
 460 / Supelcowax10(FAEES) T:0.25um L:30.0m D:0.25mm
 966 / Equity-1(Hydro) T:0.25um L:30.0m D:0.25mm

Library search

• **Registered compounds**
4,030 flavor and fragrances compounds


• **Linear Retention Indices (LRI)**
Non-polar column (SLB®-5ms)
Highly-polar column (SUPELCOWAX® 10)
Non-polar column (Equity®-1)

• **Registered information**
Mass spectrum, retention index for each column,
CAS number, compound name, molecular weight,
compositional formula

FFNSC4 Library

Flavour & Fragrance Natural & Synthetic Compounds

Crushing affects Terpene analysis


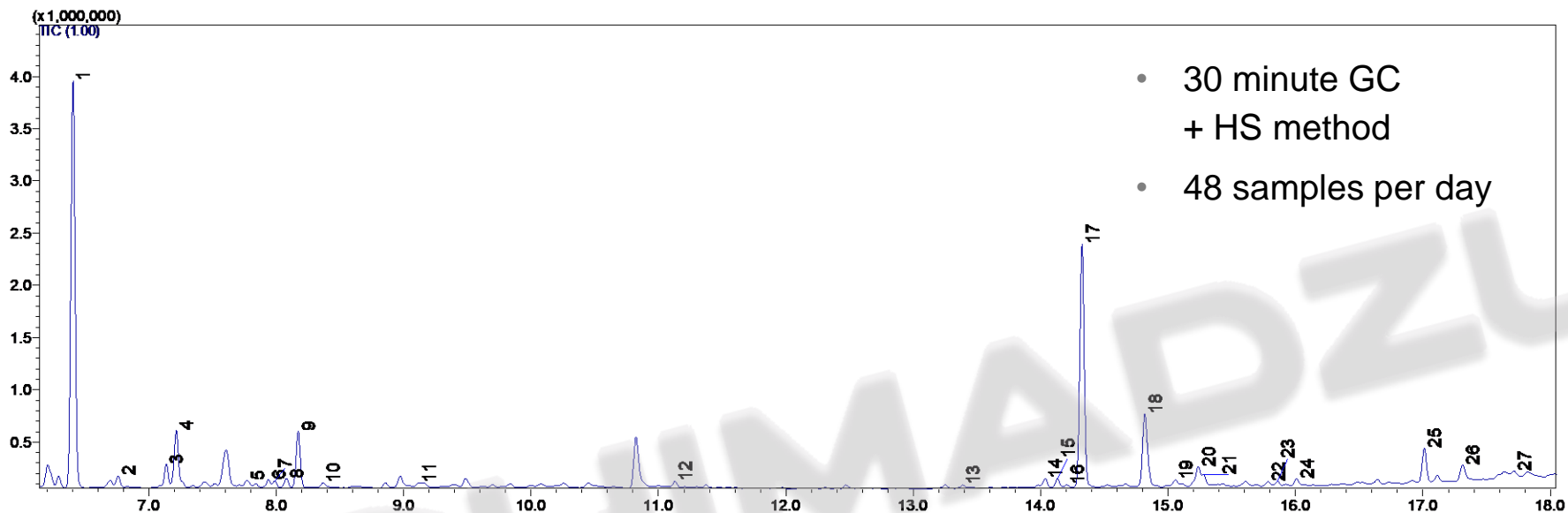


- 30 minute GC + HS method
- 48 samples per day

Peak #	Compound	Peak #	Compound
1	Pinene <alpha->	14	Bicyclo[7.2.0]undec-4-ene, 4,11,11-trimethyl-8-methylene-, [1R-(1R*,4Z,9S*)]-
2	Camphene	15	Bergamotene <alpha-, cis->
3	.beta.-Myrcene	16	Farnesene <(E)-, beta->
4	Pinene <beta->	17	Caryophyllene <(E)->
5	Carene <delta-2->	18	.alpha.-Humulene
6	Limonene	19	Selina-4,11-diene
7	Cymene <para->	20	Selinene <beta->
8	Carene <delta-3->	21	.alpha.-Selinene
9	Eucalyptol	22	Selina-4(15),7(11)-diene
10	Terpinene <gamma->	23	Selina-3,7(11)-diene
11	Sabinene hydrate <trans->	24	Nerolidol <(E)->
12	Terpineol <alpha->	25	Caryophyllene oxide
13	Ylangene <alpha->	26	Humulene epoxide II
		27	.alpha.-Bisabolol

Milling affects Terpene analysis

Terpenes in Hemp tea (differently crushed)

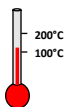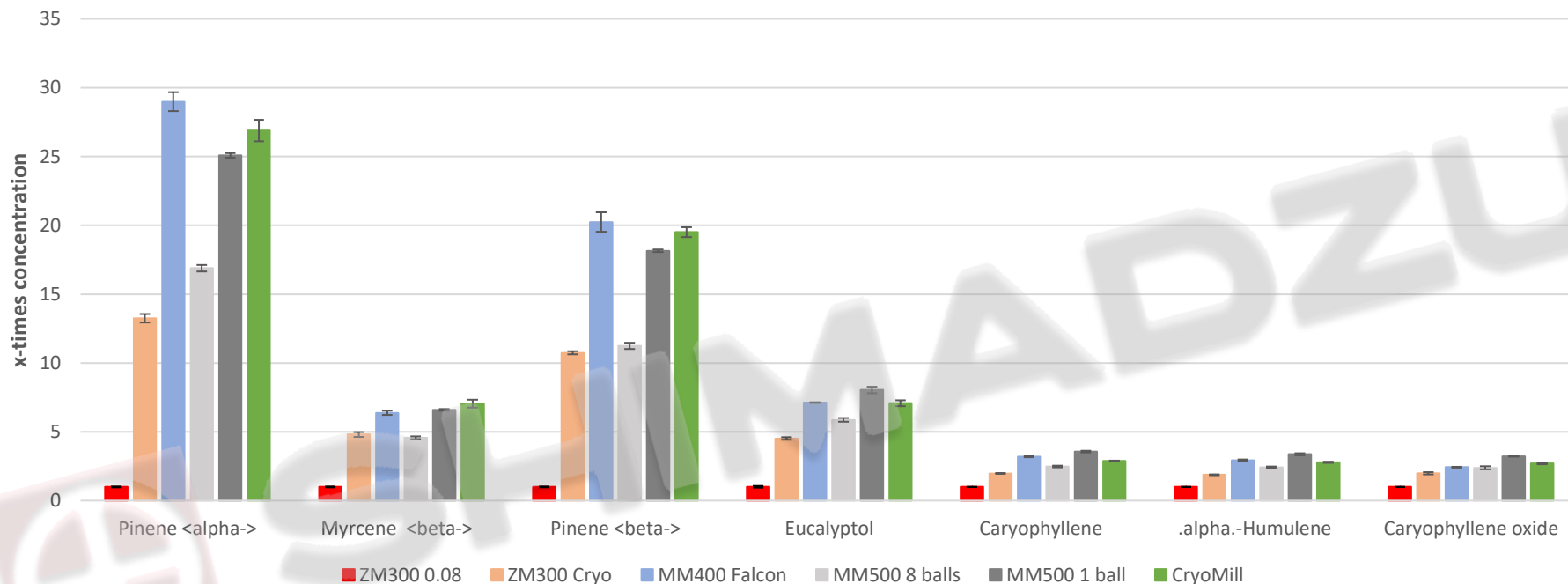

-190°C at the start



-20°C



-100°C



-100°C




-196°C

Smart Aroma Database

Database for GC-MS(MS) Aroma Analysis

Smart Aroma Database Database with 500 Aroma compounds

- A method (SIM or MRM) could be created without measuring standards first
- Only an n-Alkanstandard is necessary to calculate retention times

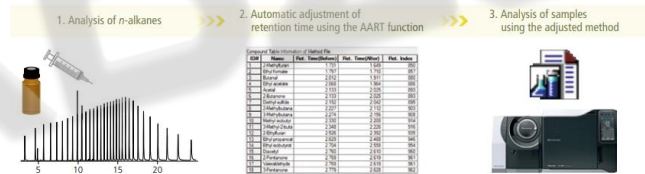

Create Method File

Instrument Type: TO Series Lang: []

Parameter

Ret. Index for AART: [Ret. Index 1]

n-alkane data file: []

Template Method File: []

Divide Method into: [1]


Advanced

Acq. Mode	Compound Name (E)	Ret. Index 1	Cas#	Comment (E)
Scan	alpha-Pinene	935	80-56-8	pine, turpentine
SIM	Myrcene	990	123-35-3	balsamic, must, spice
MRM	delta-3-Carene	1009	13466-78-9	lemon, resin
Scan	alpha-Terpinene	1017	99-86-5	lemon
Scan	p-Cymene	1025	99-87-6	solvent, gasoline, citrus
Scan	Limonene	1030	138-86-3	lemon, orange

MRM Transition				m/z for SIM or Scan									
Type	m/z	CE	Ratic	Type	m/z	CE	Ratic	Type	m/z	Ratic	Type	m/z	Ratic
T	93.00>77.00	12	100.00	Ref.1	93.00>51.00	27	39.91	T	93.0	100.00	Ref.1	121.0	10.43
T	93.00>77.10	12	100.00	Ref.1	69.00>41.10	6	87.54	T	93.1	100.00	Ref.1	91.1	30.08
T	136.10>93.10	9	100.00	Ref.1	136.10>121.10	9	76.90	T	136.1	100.00	Ref.1	121.1	106.83
T	121.10>93.10	6	100.00	Ref.1	136.10>121.10	12	71.07	T	136.1	100.00	Ref.1	121.1	156.15
T	134.10>119.10	9	100.00	Ref.1	119.10>91.10	15	80.53	T	134.1	100.00	Ref.1	119.1	321.93
T	136.00>93.00	15	100.00	Ref.1	107.00>91.00	12	104.64	T	93.0	100.00	Ref.1	121.0	29.51

Hemp tea vs. Cannabis bud

- Measured Hemp tea had a lower concentration of terpenes, and the terpene profile was less complex as in cannabis buds.
- Possible reasons: Different Hemp strain (Finola), Mixture from leaves and buds, different drying procedure, Age of the samples...

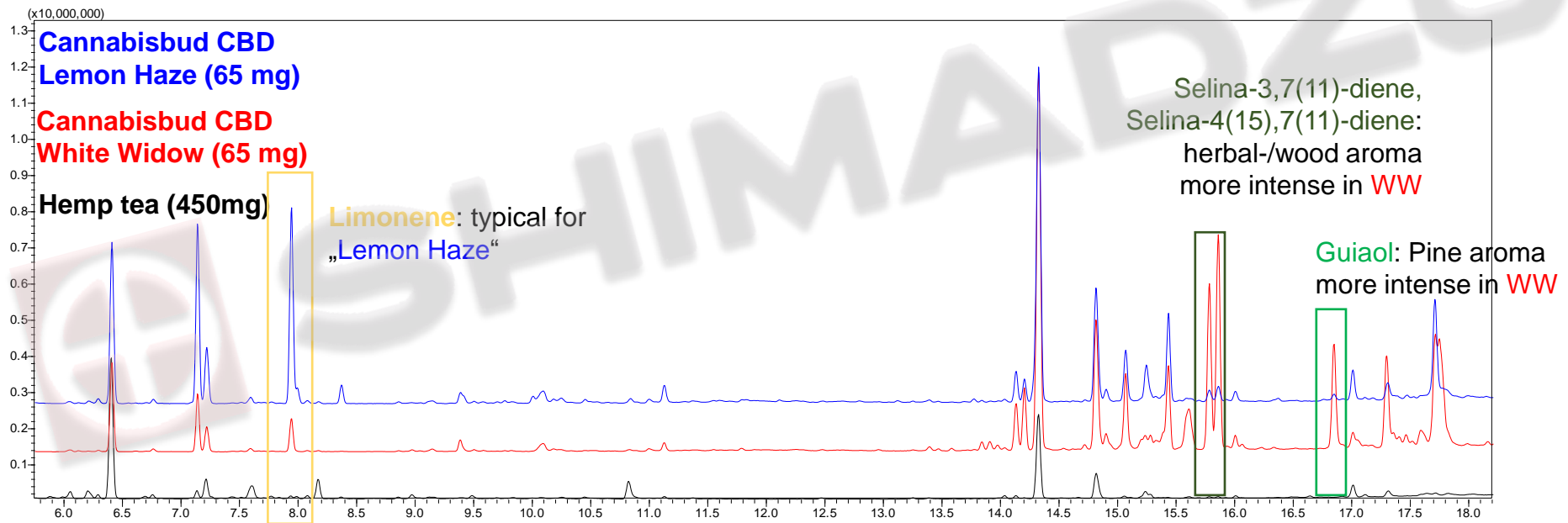

Summary

- Terpene analysis is easier over Headspace-GC than via LC
- MS or FID are both suitable detectors
- Cannabis samples
 - The temperature during the milling process of the cannabis samples has a strong effect on the terpene profile and their concentrations
 - Terpene content depends on harvesting, storage and drying conditions.
 - Terpenes can provide a reference to the variety (sativa vs. indica) and can provide a "variety-specific" fingerprint



Disclaimer



Shimadzu does not support or promote the use of its products or services in connection with illegal use, cultivation or trade of cannabis products. Shimadzu products are intended to be used for research use only purposes or state approved medical research. Shimadzu is not condoning the use of recreational nor medical marijuana, we are merely providing a market summary of the cannabis testing industry.