



C₁ – C₄ hydrocarbons

Application Note

Energy & Fuels

Authors

Agilent Technologies, Inc.

Introduction

GC analysis of C₁ - C₄ hydrocarbons is achieved in 15 minutes using the Agilent Select Al₂O₃ MAPD.



Agilent Technologies


Conditions

Technique : GC
Column : Agilent Select Al₂O₃ MAPD, 0.53 mm x 50 m fused silica (Part no. CP7432)
Temperature : 40 °C, 5min → 160 °C, 10 °C/min → 200 °C, 20 °C/min, hold 1 min
Carrier Gas : He, 4 psig, 4 min → 11 psig, 0.5 psig/min, 2 min
Injector : Split 60 mL/min
Detector : FID
Concentration Range : approx 100 ppm in nitrogen, synthetic standard

Courtesy : J. Luong, Dow Chemical Canada

Peak identification

1. methane
2. propane
3. propylene
4. iso-butane
5. trans-2-butene
6. 1-butene
7. iso-pentane
8. cis-2-butene
9. methyl acetylene
10. vinyl acetylene

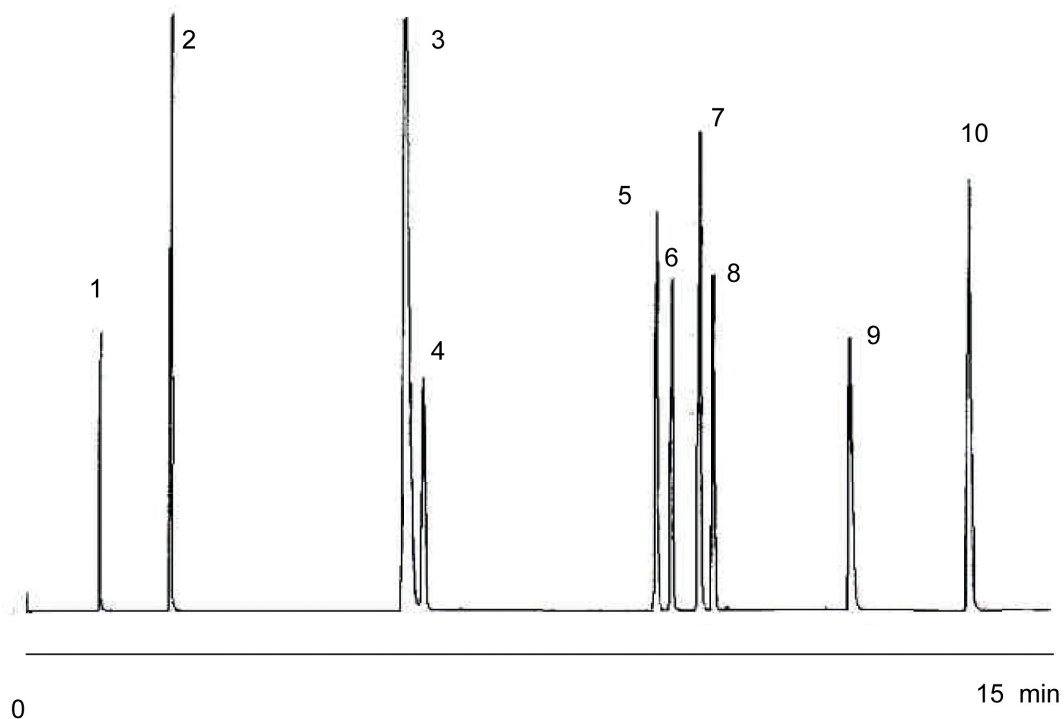

www.agilent.com/chem

This information is subject to change without notice.

© Agilent Technologies, Inc. 2011

Printed in the USA

31 October, 2011

First published prior to 11 May, 2010

A02042



Agilent Technologies