

# Glycosylation and Sialic Acid Analysis of Biotherapeutic Glycoproteins

Tom Rice  
Bioconsumables Product Manager

August 20, 2020

# Outline

Importance of N-glycan analysis for biotherapeutic development

Dye options for released N-glycan analysis

- InstantPC for HILIC (rapid, high FLD and MS signal)
- 2-AB for HILIC (well-established)

HILIC separations of labeled N-glycans

- FLD and MS detection

Sialic acid quantitation and analysis workflow

- Total sialic acid quantitation (plate based)
- Sialic acid species profiling (LC/FLD/MS)

# N-Glycans on Biotherapeutics

1/5<sup>th</sup> of all proteins in SwissProt are glycoproteins

More than 60% of biotherapeutics are glycosylated

N-Glycan structure can affect pharmacokinetics, pharmacodynamics, and immunogenicity

Glycosylation can be a critical quality attribute (CQA)

Even if not a CQA, glycosylation still monitored as a product quality attribute

Monoclonal antibody (mAb) and antibody-derived biotherapeutic proteins are glycosylated

Other molecules including blood factors (EPO, Factor VIII) also glycosylated

[Khoury, et al., \*Nat. Sci. Rep.\* 1 \(90\), doi:10.1038/srep00090 \(2011\)](#)

[Planinc et al., \*Anal. Chim. Acta\* 921, 13–27 \(2016\)](#)

[Walsh, \*Nat. Biotechnol.\* 32\(10\), 992–1000 \(2014\)](#)

[Liu, \*J. Pharm. Sci.\* 104 \(6\), 1866–1884 \(2015\)](#)

# Top Selling Pharmaceuticals (2010)

	Trade name	Type	Main Indications	Company	Sales (USD millions/year)
1	Lipitor	Small molecule	Hypercholesterolemia, Dyslipidemia	Pfizer Inc.	7,244
2	Nexium	Small molecule	Symptomatic Gastroesophageal Reflux Disease	AstraZeneca Pharmaceuticals LP	6,309
3	Plavix	Small molecule	Reduction of atherosclerotic events	Bristol-Myers Squibb Company	6,129
4	Advair Diskus	Small molecule	Chronic obstructive pulmonary disease	GlaxoSmithKline	4,711
5	Abilify	Small molecule	Bipolar Mania	Bristol-Myers Squibb Company	4,551
6	Seroquel	Small molecule	Bipolar Depression	AstraZeneca Pharmaceuticals LP	4,349
7	Singular	Small molecule	Asthma	Merck & Co., Inc.	4,072
8	Crestor	Small molecule	Hypercholesterolemia, dyslipidemia	AstraZeneca Pharmaceuticals LP	3,758
9	Actos	Small molecule	Type 2 diabetes	Takeda	3,534
10	Epogen	Biologic	Anemia	Amgen	3,323

 Glycosylated

- Only 1 glycosylated molecule in 2010 Top 10

\*IMS Institute for Healthcare Informatics. The use of medicines in the United States: review of 2010.

# Top Selling Pharmaceuticals (2018)

Rank	Drug	Trade names	Type	Main indications	Company	Sales (USD billions/year)
1	adalimumab	Humira	Biologic	Autoimmune	AbbVie	19.936
2	apixaban	Eliquis	Small molecule	Anticoagulation	Bristol-Myers Squibb/Pfizer	9.872
3	lenalidomide	Revlimid	Small molecule	MDS, multiple myeloma, MCL	Celgene	9.685
4	nivolumab	Opdivo	Biologic	Cancer (anti-PD-1 immunotherapy)	Bristol-Myers Squibb/Ono Pharmaceutical	7.570
5	pembrolizumab	Keytruda	Biologic	Cancer (anti-PD-1 immunotherapy)	Merck & Co.	7.171
6	etanercept	Enbrel	Biologic	Autoimmune including RA	Amgen/Pfizer	7.126
7	trastuzumab	Herceptin	Biologic	Breast cancer	Roche (Genentech)	6.981
8	bevacizumab	Avastin	Biologic	Cancer types, AMD	Roche (Genentech)	6.847
9	rituximab	Rituxan, MabThera	Biologic	Autoimmune, cancer types	Roche (Genentech)/Biogen	6.750
10	rivaroxaban	Xarelto	Small molecule	Anticoagulation	Bayer/J&J	6.589



 Glycosylated

- 7/10 glycosylated biotherapeutics

Top-selling drugs are ranked based on sales or revenue reported for 2018 by bio/pharma companies in press announcements, annual reports, investor materials, and/or conference calls.

<https://www.genengnews.com/a-lists/top-15-best-selling-drugs-of-2018/>

# Effect of N-Glycan Structure on Biotherapeutics

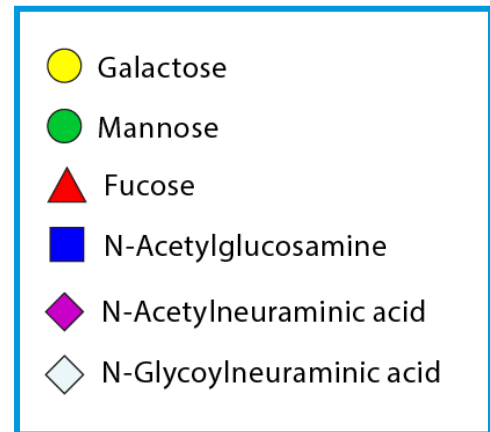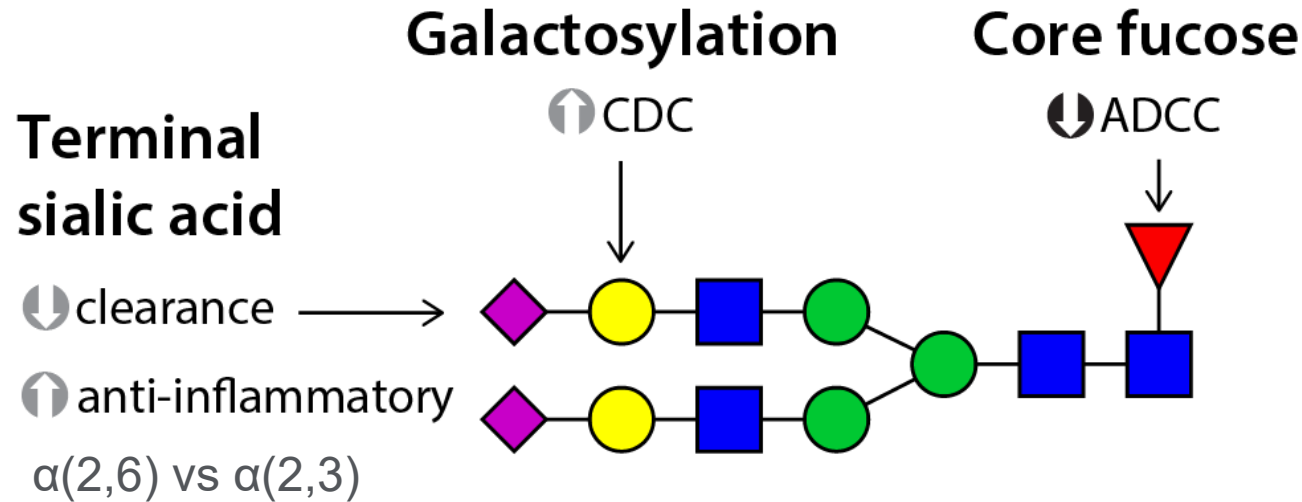

- ADCC: Antibody-Dependent Cell-mediated Cytotoxicity
- CDC: Complement-Dependent Cytotoxicity

Reviewed in:


Jones, *BioPharm Intl.* 2017; 30:20-25

Higel et al., *Eur J Pharm Biopharm.* 2016; 100: 94

Liu, *J Pharm Sci.* 2015; 104: 1866-1884

Abes & Teillaud, *Pharmaceuticals* 2010; 3:146-57

# Effect of N-Glycan Structure on Biotherapeutics

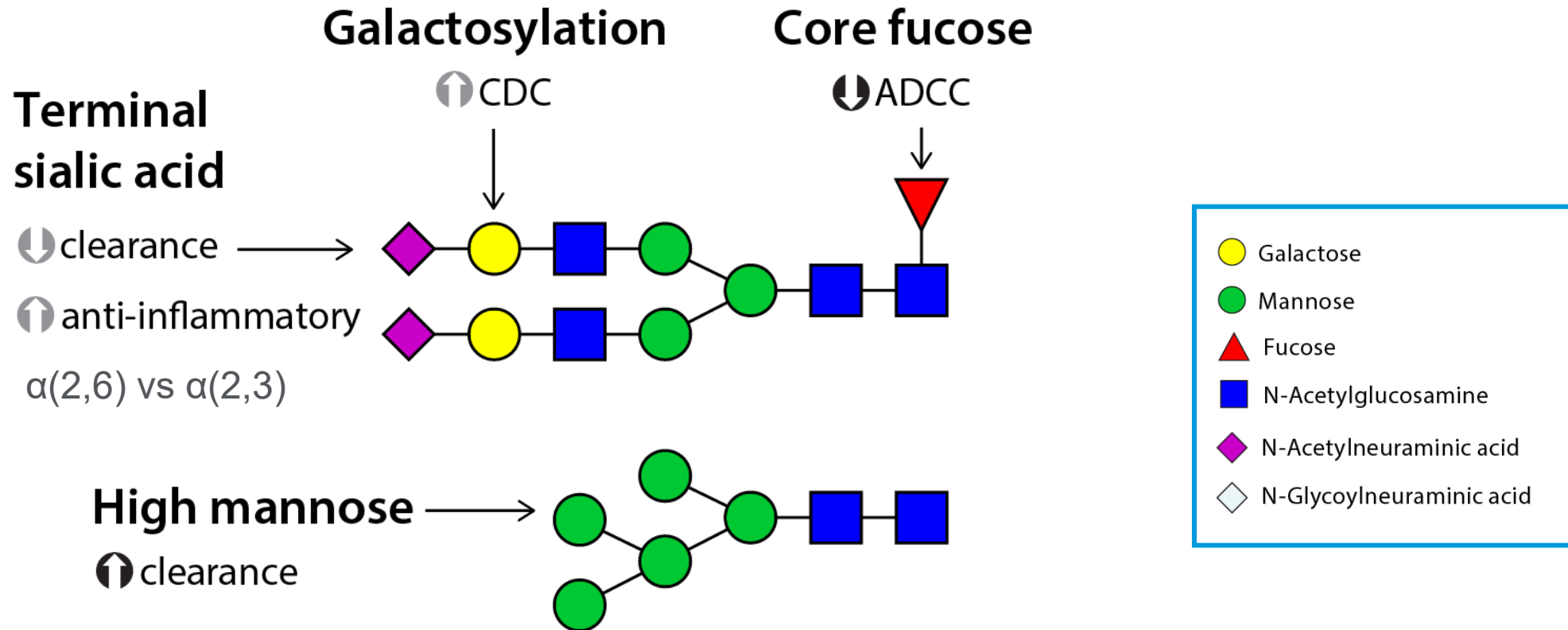

Reviewed in:


Jones, *BioPharm Intl.* 2017; 30:20-25

Higel et al., *Eur J Pharm Biopharm.* 2016; 100: 94

Liu, *J Pharm Sci.* 2015; 104: 1866-1884

Abes & Teillaud, *Pharmaceuticals* 2010; 3:146-57

# Non-Human/Immunoreactive N-Glycans

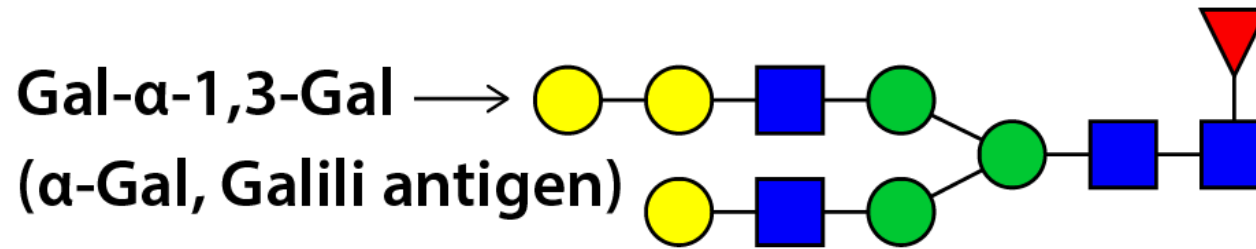


Cetuximab - N-glycans with  $\alpha$ -Gal on Fab region

- Associated with adverse immunological responses

Abatacept - N-glycans with  $\alpha$ -Gal on Fc region

- No reported adverse effects

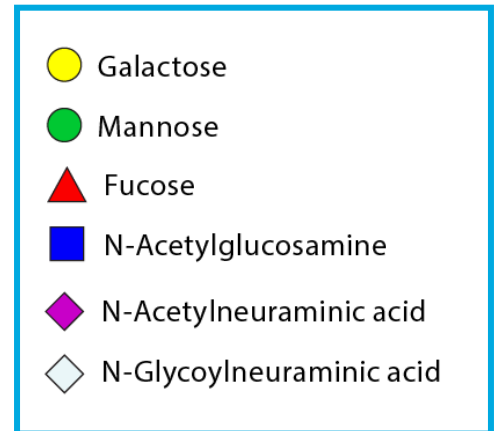

Macher & Galili, *Biochim Biophys Acta*. 2010; 1780 (2): 75-88




Van Bueren et al., *Nat Biotechnol*. 2011; 29: 574-576

Bosques et al., *Biotechnol*. 2010; 28: 1153-1156

Chung et al., *N Eng J Med*. 2008; 358(11): 1109-1117




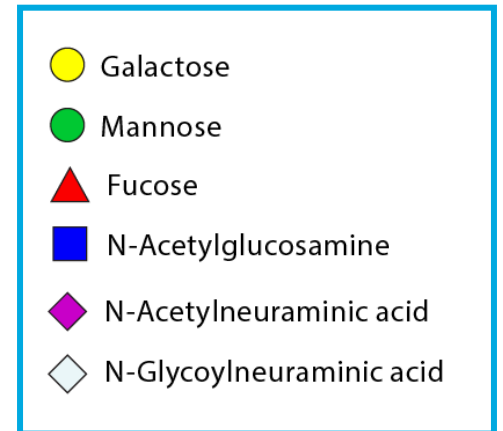
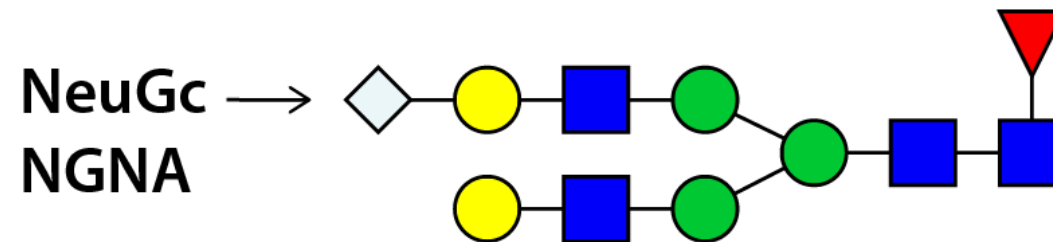
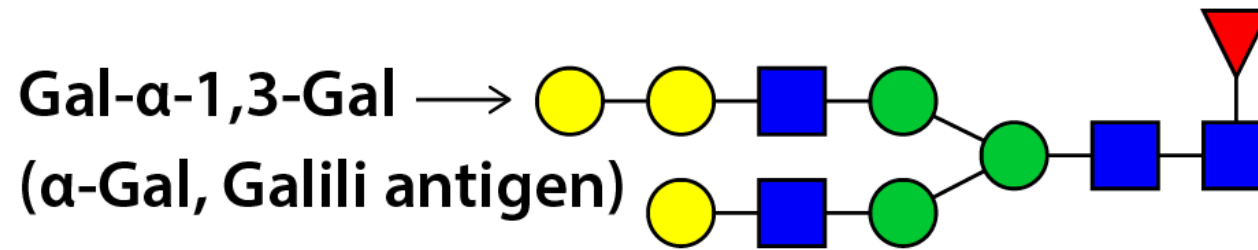


# Non-Human/Immunoreactive N-Glycans



Potentially immunogenic

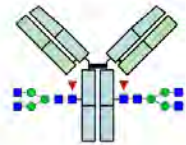
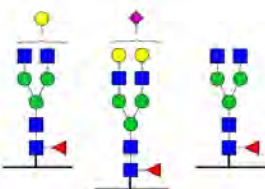
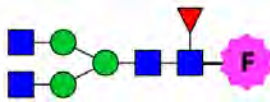
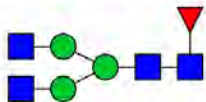


Ghaderi et al., *Biotechnol Genet Eng Rev.* 2012; 28:147-75

# N-Glycan Analysis Options


Material	Structure	Analytical Methods
Intact glycoproteins		Electrophoresis, lectin arrays, LC-ESI-MS: HRAM (+ reduction)
Glycopeptides		LC-MS, MALDI, CE
Released glycans		
▶ Derivatized		LC-FLR, LC-MS, CE-LIF Exoglycosidases (linkage assignment)
▶ Underivatized		MALDI-TOF, HPAE-PAD
Monosaccharides		
▶ Derivatized		RP LC-FLR
▶ Underivatized		HPAE-PAD

[Jones, BioPharm Intl. 2017; 30:20-25](#)

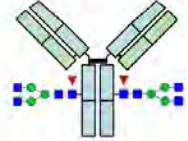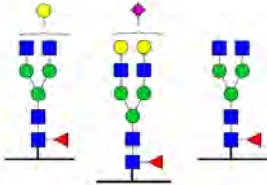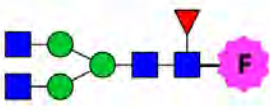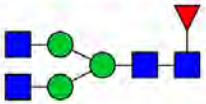
# N-Glycan Analysis Options

Material	Structure	Analytical Methods	
Intact glycoproteins		Electrophoresis, lectin arrays, LC-ESI-MS: HRAM (+ reduction)	
Glycopeptides		LC-MS, MALDI, CE	
Released glycans			
<ul style="list-style-type: none"> <li>▶ Derivatized</li> </ul>		LC-FLR, LC-MS, CE-LIF Exoglycosidases (linkage assignment)	→ Relative % area of glycan species
<ul style="list-style-type: none"> <li>▶ Underivatized</li> </ul>		MALDI-TOF, HPAE-PAD	
Monosaccharides			
<ul style="list-style-type: none"> <li>▶ Derivatized</li> </ul>		RP LC-FLR	→ Sialic acid quantitation
<ul style="list-style-type: none"> <li>▶ Underivatized</li> </ul>		HPAE-PAD	

# Agilent Released N-Glycan Workflow

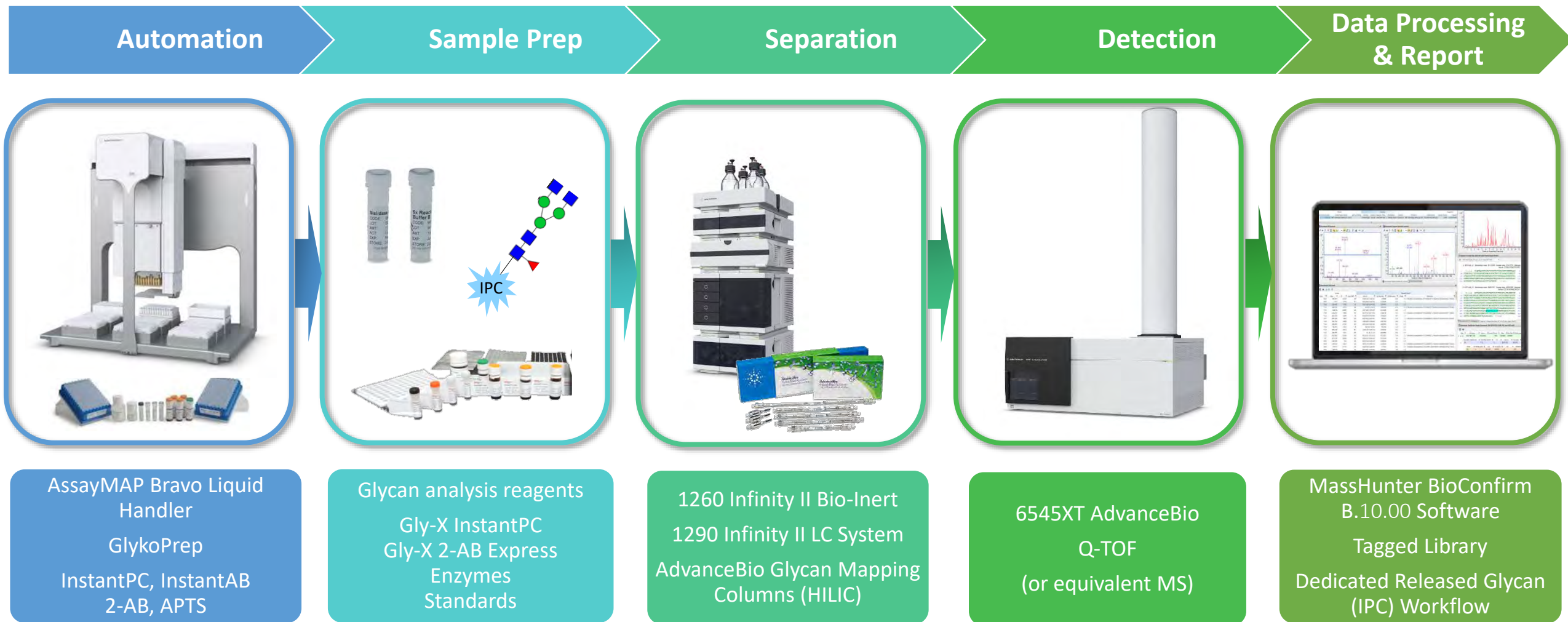


# N-Glycan Sample Prep Evolution

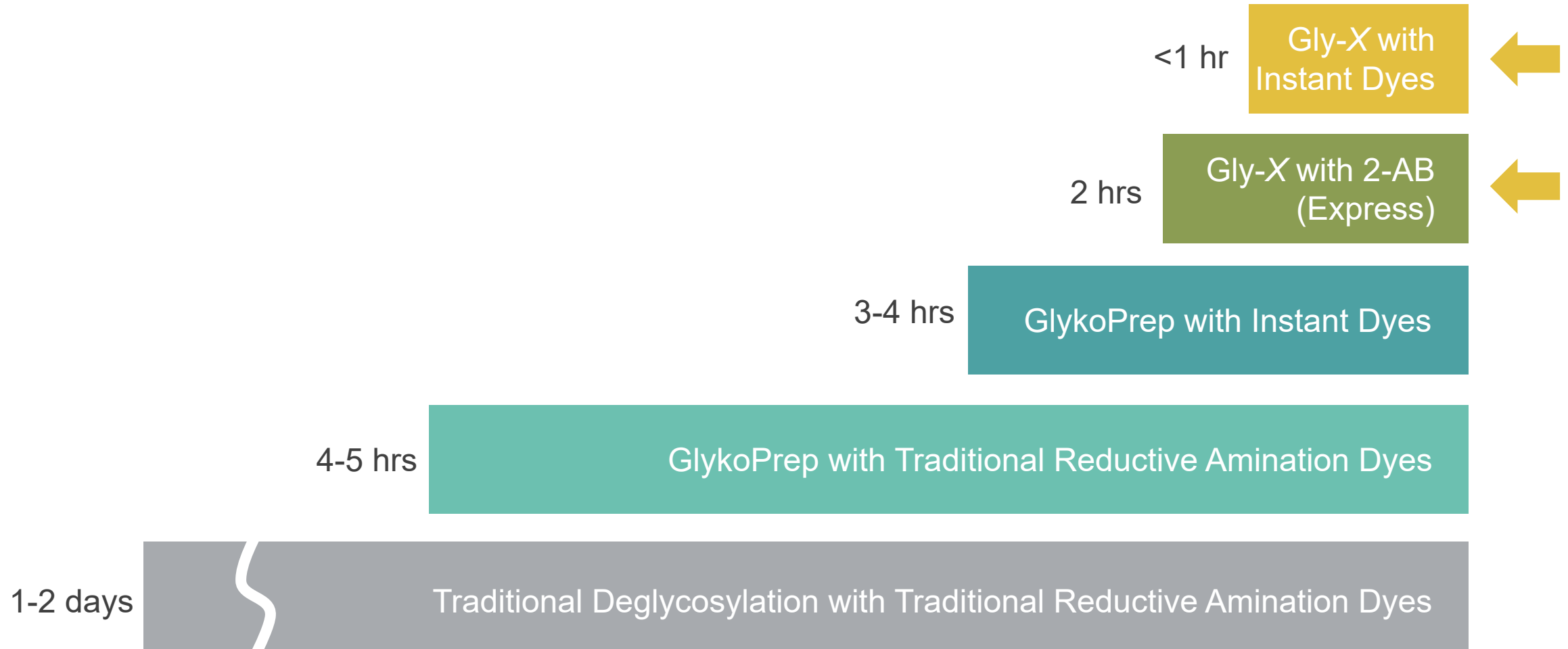


# Gly-X N-Glycan Sample Prep Technology


1-40 µg glycoprotein sample\*, 0.05-2 mg/ml

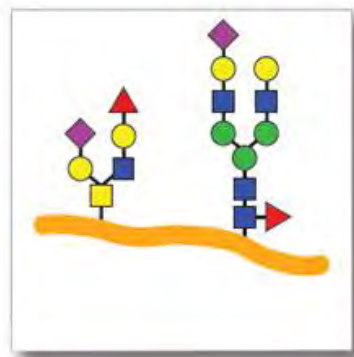


Denature  
3 min at 90 °C

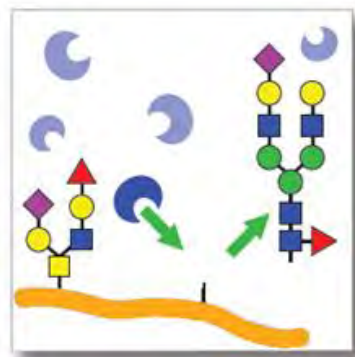


Deglycosylate  
PNGase F, 5 min  
at 50 °C

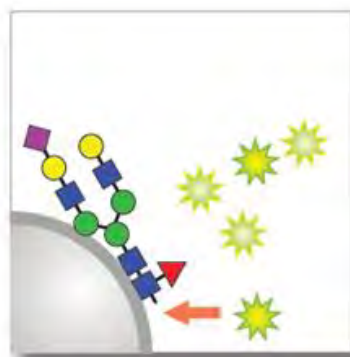

Label  
InstantDye, 1 min  
*InstantPC, InstantQ, InstantAB*  
Reductive amination dye, 60 min  
(no dry down)  
*2-AB, APTS*

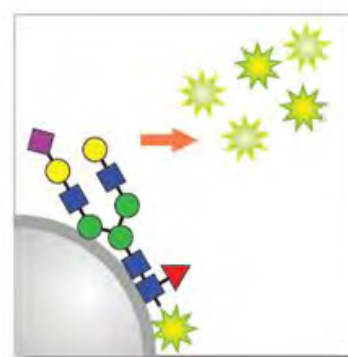

Clean up  
15-20 min at  
RT

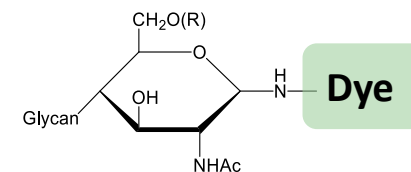

Labeled N-Glycans  
LC/FLD, LC/MS, CE




Workflow time: 45 min – 120min (16 samples)

\*Loading range depends on protein, can go up to 100 µg for mAbs e.g. Rituxan

# Novel In-Solution Enzymatic Protein Denaturation & Deglycosylation


Gly-X in-solution 5-minute deglycosylation with PNGase F achieves > 99% N-glycan release for most proteins tested

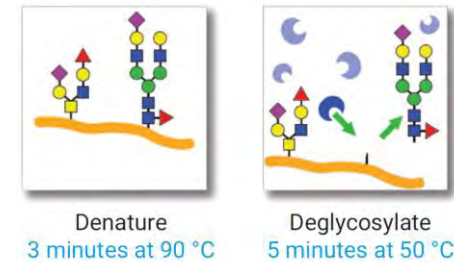

Protein	% Deglycosylation
Rituximab	>99%
Cetuximab	>99%
Enbrel	>99%
Zaltrap	>99%

## Analytical Method:



Agilent 2100 Bioanalyzer, Protein 230 Kit  
(electropherograms not normalized)

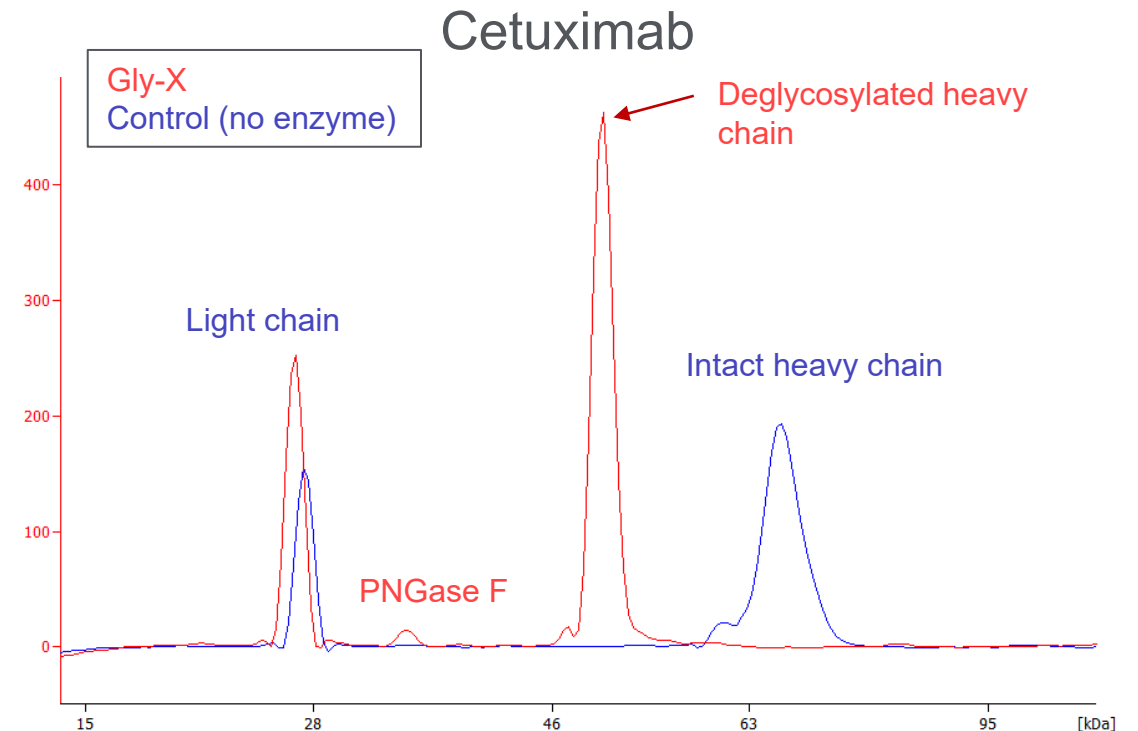


# N-Glycan Label Choices

## InstantDyes:


Glycoprotein  $\xrightarrow{\text{N-Glycanase}}$

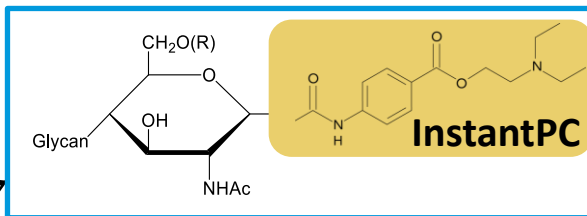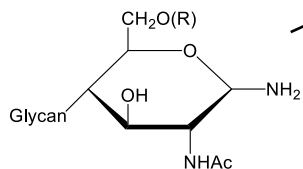

- InstantDye for LC/FLD/MS
- Strong FLD and MS signal

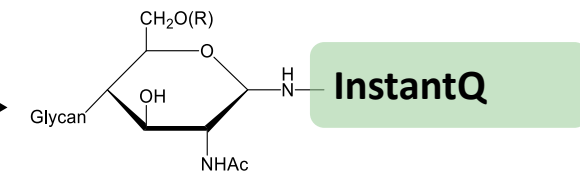



- InstantDye for CE on Gly-Q System
- Introduces negative charge for CE

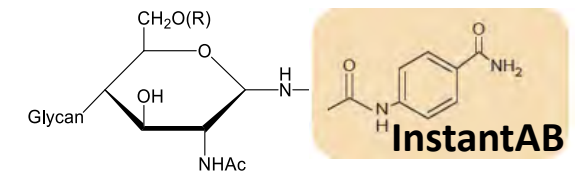


- 1<sup>st</sup> generation InstantDye for LC/FLD
- Lower MS signal

## Reductive Amination:


Glycoprotein  $\xrightarrow{\text{N-Glycanase}}$

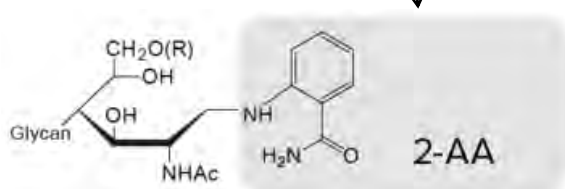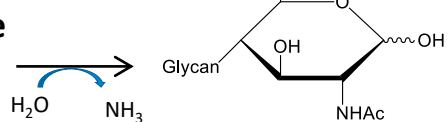

- Standards available
- Labeling not supported

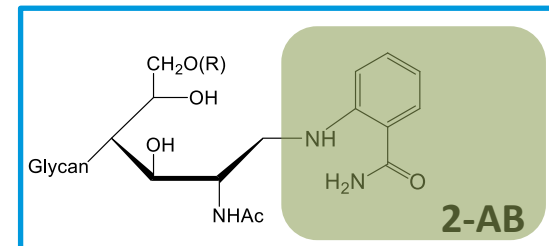



- Label makes N-Glycan *less* polar
- Improves HILIC separation for UHPLC
- Well-established glycan label

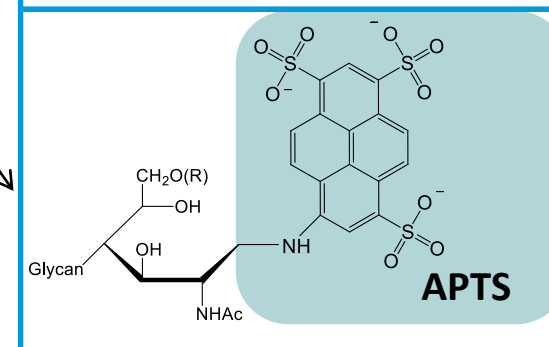

- Label makes N-Glycan *more* polar
- Introduces negative charge for CE




# InstantPC Dye (IPC)

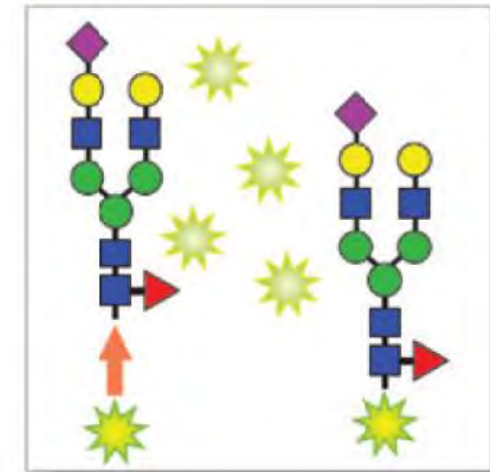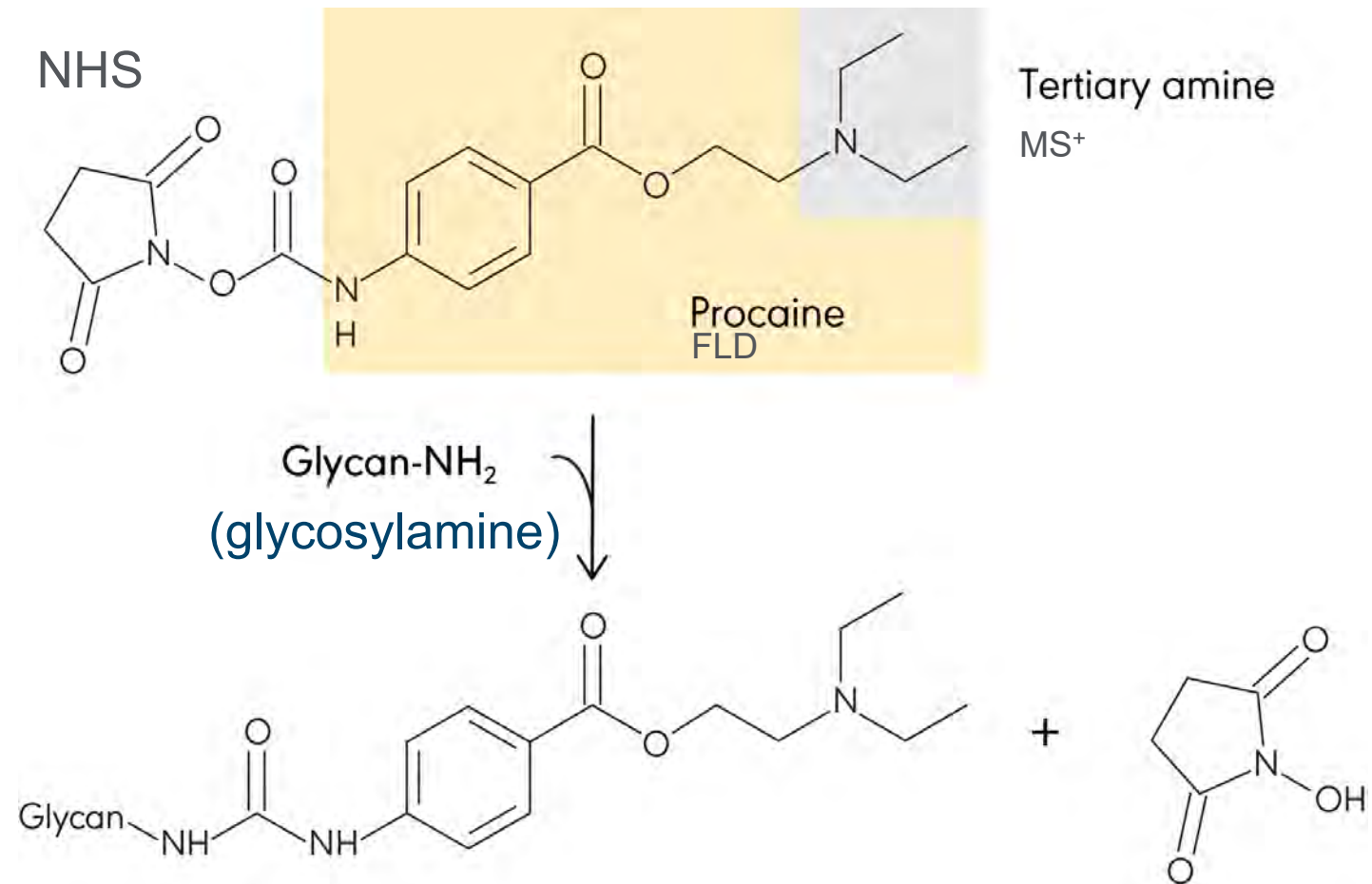

Label  
1 minute at 50 °C

Workflow time: 45 min (16 samples), 100 min (96 samples)

# FLD and MS Response Comparison

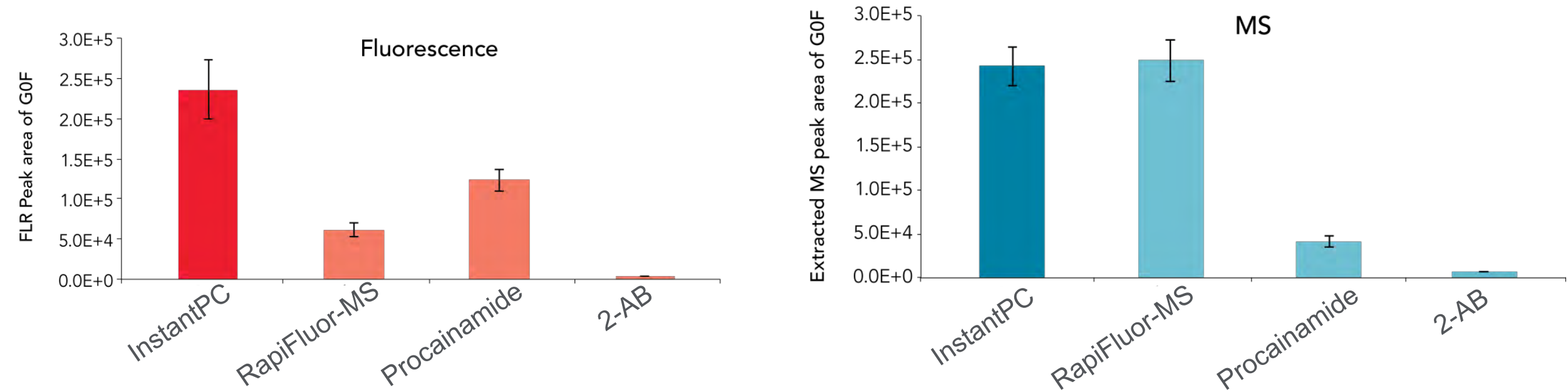


Labeling uses same amount of released glycans from GlykoPrep digestion


*Reference: ProZyme Poster Presentation at ASMS, St. Louis, MO, May 31 - June 4, 2015*

# UHPLC-HILIC Separation of Labeled N-Glycans

Separation

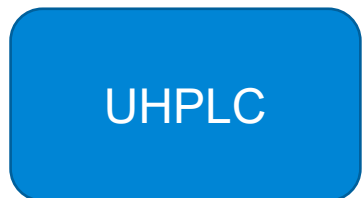

Detection

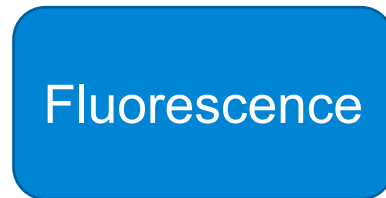

Amide HILIC column



- 1.8  $\mu\text{m}$  fully porous for speed and performance (1200 bar)
- 2.7  $\mu\text{m}$  superficially porous for resolution at lower pressures (600 bar)
- Unique hydrophilic bonding
- MS and FLD compatibility

- InstantPC: ex/em 285/345 nm
- 2-AB: ex/em 260/430 nm



- LC/MS mass confirmation for glycan ID
- Positive mode

HILIC: hydrophilic interaction liquid chromatography  
Example Method: 1 hour

Time (min)	%A	%B	Flow (ml/min)
0.00	20.0	80.0	0.50
2.00	25.0	75.0	0.50
48.00	38.0	62.0	0.50
49.00	60.0	40.0	0.50
51.50	60.0	40.0	0.50
52.00	18.0	82.0	0.50
60.00	18.0	82.0	0.50


AdvanceBio Glycan Mapping Column  
2.1 x 150 mm, 1.8  $\mu\text{m}$

A: 50 mM ammonium formate, pH 4.5  
B: Acetonitrile

Column temperature: 40  $^{\circ}\text{C}$

# Rituximab InstantPC Glycans

	Average Rel % Area	Standard Deviation	%CV
G0F-N	0.75	0.01	1.55
G0	1.47	0.02	1.18
G0F	46.82	0.07	0.15
Man5	1.21	0.01	0.83
G1[6]	0.75	0.02	2.67
G1F[6]	31.21	0.11	0.35
G1F[3]	9.27	0.05	0.54
G2F	7.04	0.04	0.51
G2FS1[6]	0.67	0.02	2.29
G2FS1[3]	0.37	0.06	15.98
G2FS2	0.45	0.03	6.67

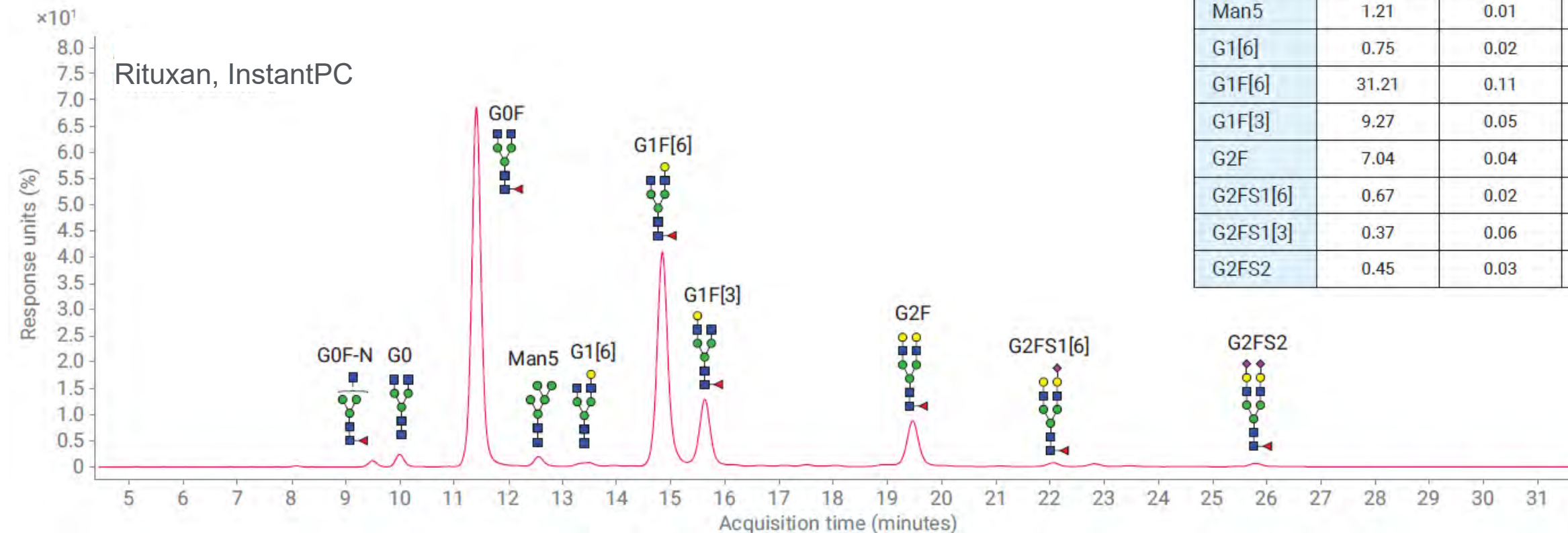

UHPLC-HILIC, 60 minute method

AdvanceBio Glycan Mapping column

2.1 x 150 mm, 1.8  $\mu$ m


FLD: 285/345 nm

1  $\mu$ l injection (equivalent to glycans from 0.4  $\mu$ g protein)

N = 4



# Etanercept InstantPC N-glycans

- Etanercept: Fc fusion protein
- Tumor necrosis factor receptor (TNFR) fused to Fc of IgG1
- 3 N-glycan sites: 1 in Fc, 2 in receptor
- N = 4

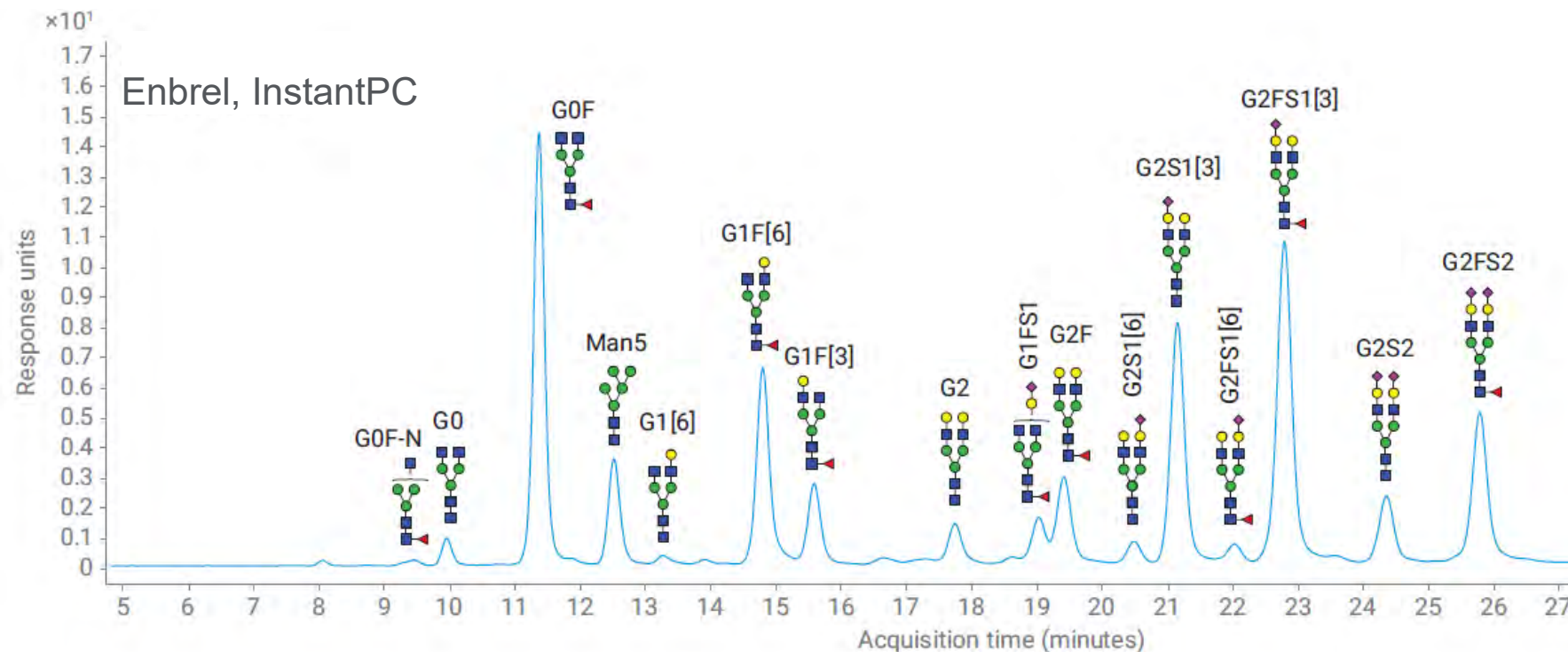

	Average Rel % Area	Standard Deviation	%CV
G0	1.10	0.02	2.09
G0F	19.36	0.16	0.84
Man5	5.08	0.03	0.52
G1[6]	0.48	0.00	0.00
G1F[6]	10.48	0.04	0.39
G1F[3]	3.97	0.01	0.25
G2	2.08	0.01	0.55
G1FS1	1.84	0.05	2.49
G2F	4.26	0.09	1.99
G2S1[6]	1.18	0.01	0.49
G2S1[3]	13.91	0.04	0.31
G2FS1[6]	0.89	0.00	0.00
G2FS1[3]	20.54	0.08	0.37
G2S2	4.26	0.01	0.14
G2FS2	10.54	0.08	0.78

# MS of InstantPC N-Glycans from Enbrel

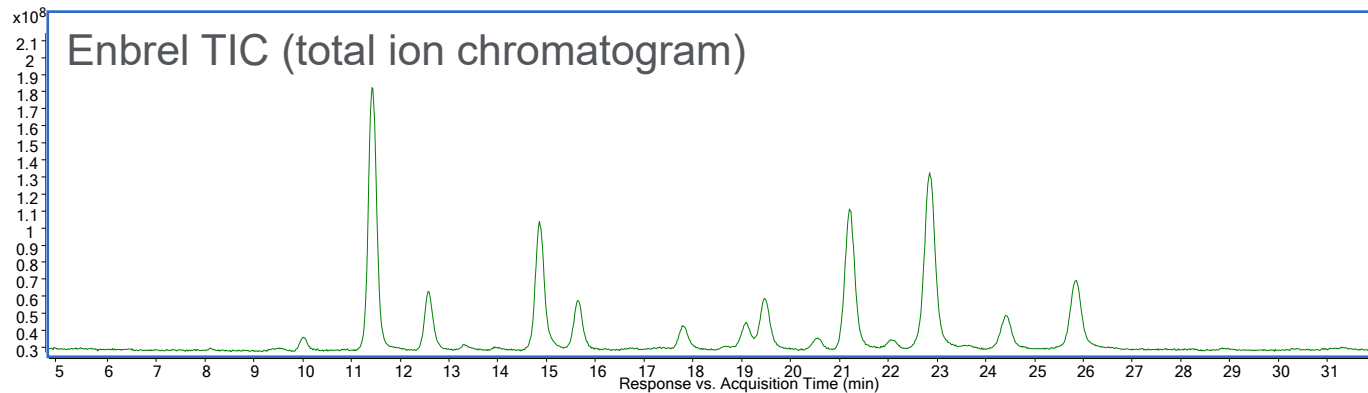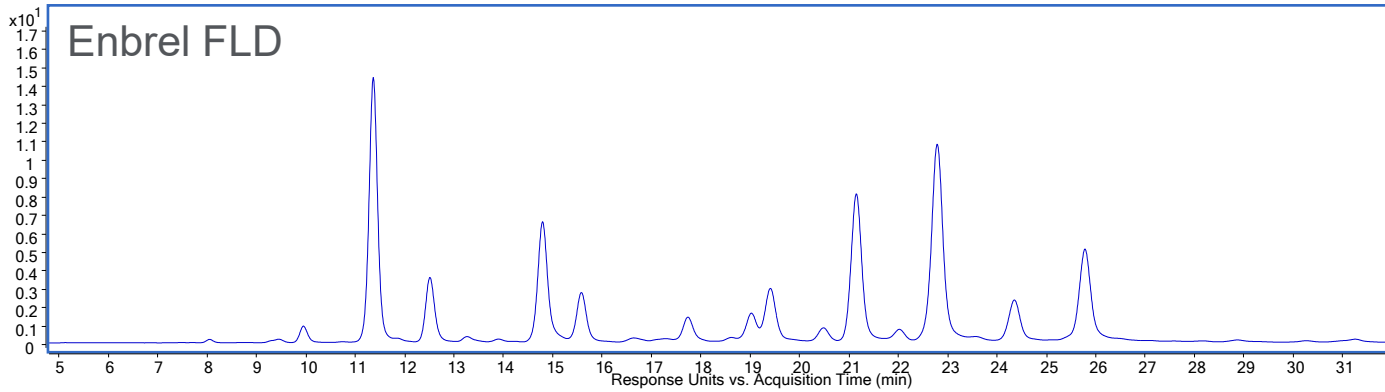

Fluorescence





Q-TOF  
6545XT

(or equivalent MS)

Positive mode MS:

Most biantennary InstantPC N-glycans  
[M+2H]<sup>2+</sup>, larger sialylated will be  
majority [M+3H]<sup>3+</sup>

Man5 mass  
spectrum

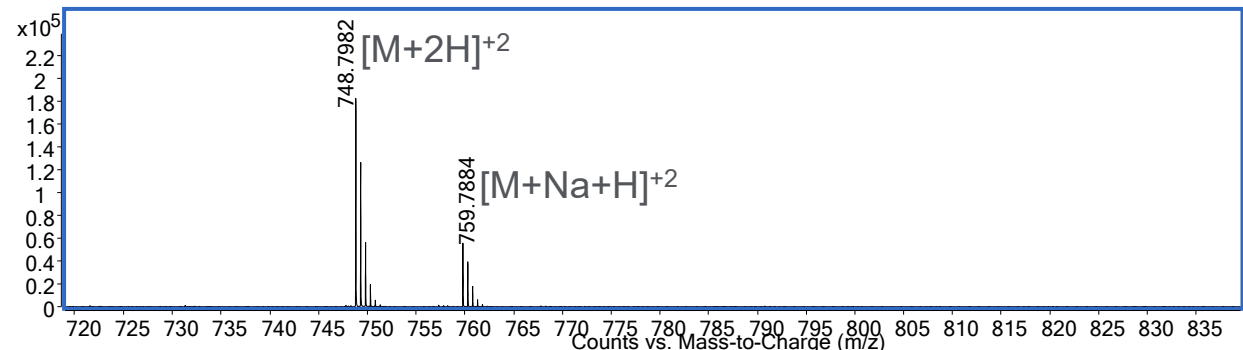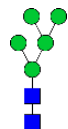

# InstantPC N-Glycan Standards & Libraries

## Individual N-Glycans

Product Code	Description	
GKPC-401	G0-N	
GKPC-301	G0	
GKPC-402	G0F-N	
GKPC-302	G0F	
GKPC-317	G1	
GKPC-316	G1F	
GKPC-304	G2	
GKPC-305	G2F	
GKPC-329	G1S1 (α2,3)	
GKPC-319	G1S1 (α2,6)	
GKPC-330	G1FS1 (α2,3)	
GKPC-320	G1FS1 (α2,6)	
GKPC-321	A1 (α2,3)	

Product Code	Description	
GKPC-311	A1 (α2,6)	
GKPC-325	A1F (α2,3)	
GKPC-315	A1F (α2,6)	
GKPC-322	A2 (α2,3)	
GKPC-312	A2 (α2,6)	
GKPC-323	A2F (α2,3)	
GKPC-313	A2F (α2,6)	
GKPC-103	Man5	
GKPC-104	Man6	
GKPC-105	Man7	
GKPC-106	Man8	
GKPC-107	Man9	

## N-Glycan Libraries

Product Code	Description
GKPC-005	Human IgG N-Linked Glycan Library
GKPC-503	Glucose Homopolymer
GKPC-233	α(2-3) Sialylated Triantennary Library
GKPC-263	α(2-6) Sialylated Triantennary Library
GKPC-234	α(2-3) Sialylated Tetraantennary Library
GKPC-264	α(2-6) Sialylated Tetraantennary Library

**New** CHO mAb glycoprotein & N-glycan library

## Labeled Glycan Standards

2-AB, 2-AA, APTS, InstantPC, InstantAB, InstantQ


Learn more: [www.agilent.com/chem/glycananalysis](http://www.agilent.com/chem/glycananalysis)

**New** Alpha Gal



Glycan Standards Flier [5994-0999EN](#)

# InstantPC Sialylated N-Glycan Standards

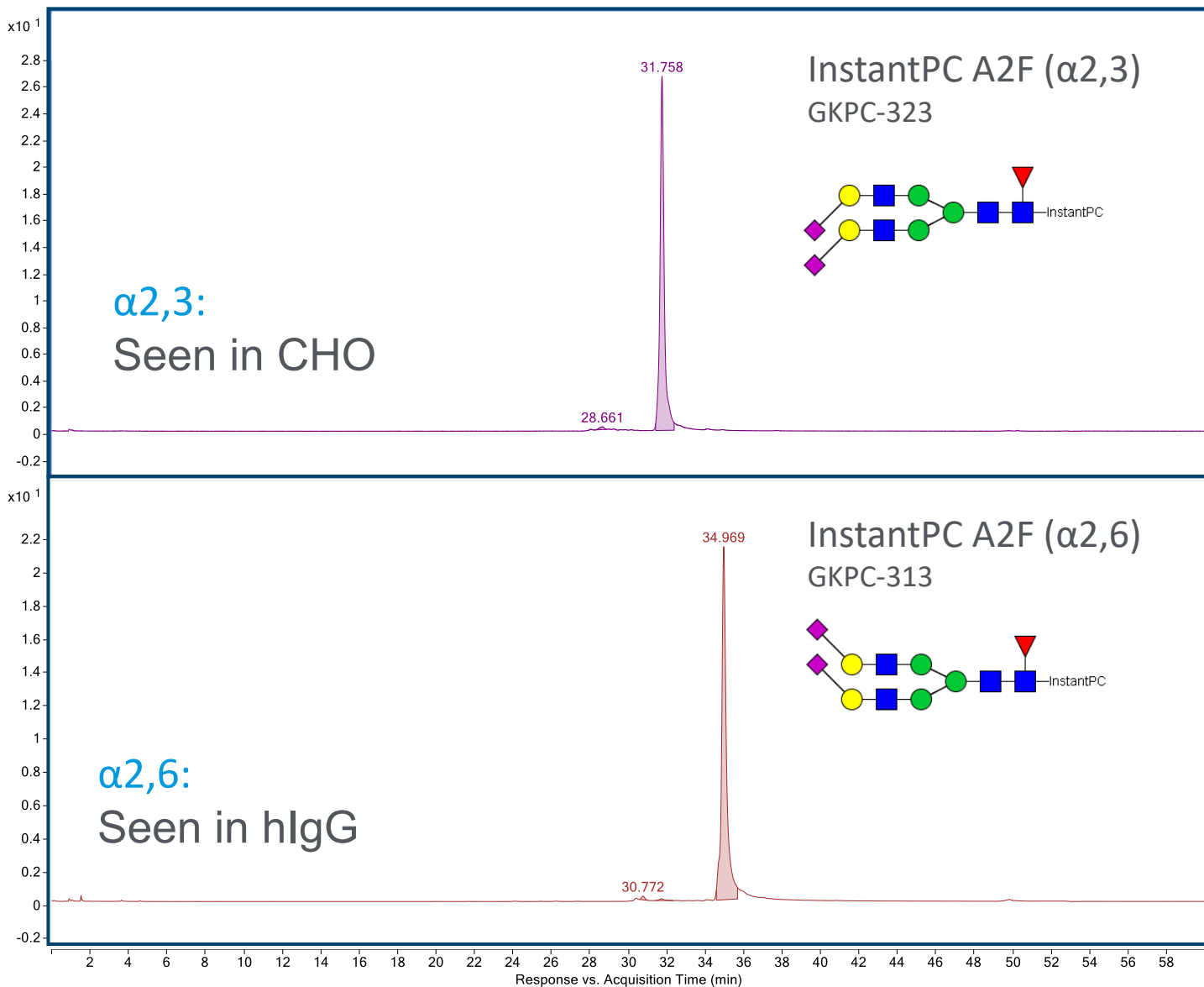


- **CHO glycoproteins:**  $\alpha$ (2,3)-linked sialic acid [1]
- **Human IgG:**  $\alpha$ (2,6)-linked sialic acid [2]
- **$\alpha$ (2,3)-sialylated N-glycans** have a shorter HILIC retention time than isomeric N-glycans with  **$\alpha$ (2,6)** sialic acid linkages [3]

1. Lee EU et al., *J Biol Chem.* 1989;264(23):13848-55.
2. Anthony RM et al. *Science.* 2008;320(5874):373-6.
3. Raymond C et al., *mAbs.* 2015;7(3):571-83.





# InstantPC Sialylated Triantennary N-Glycan Library

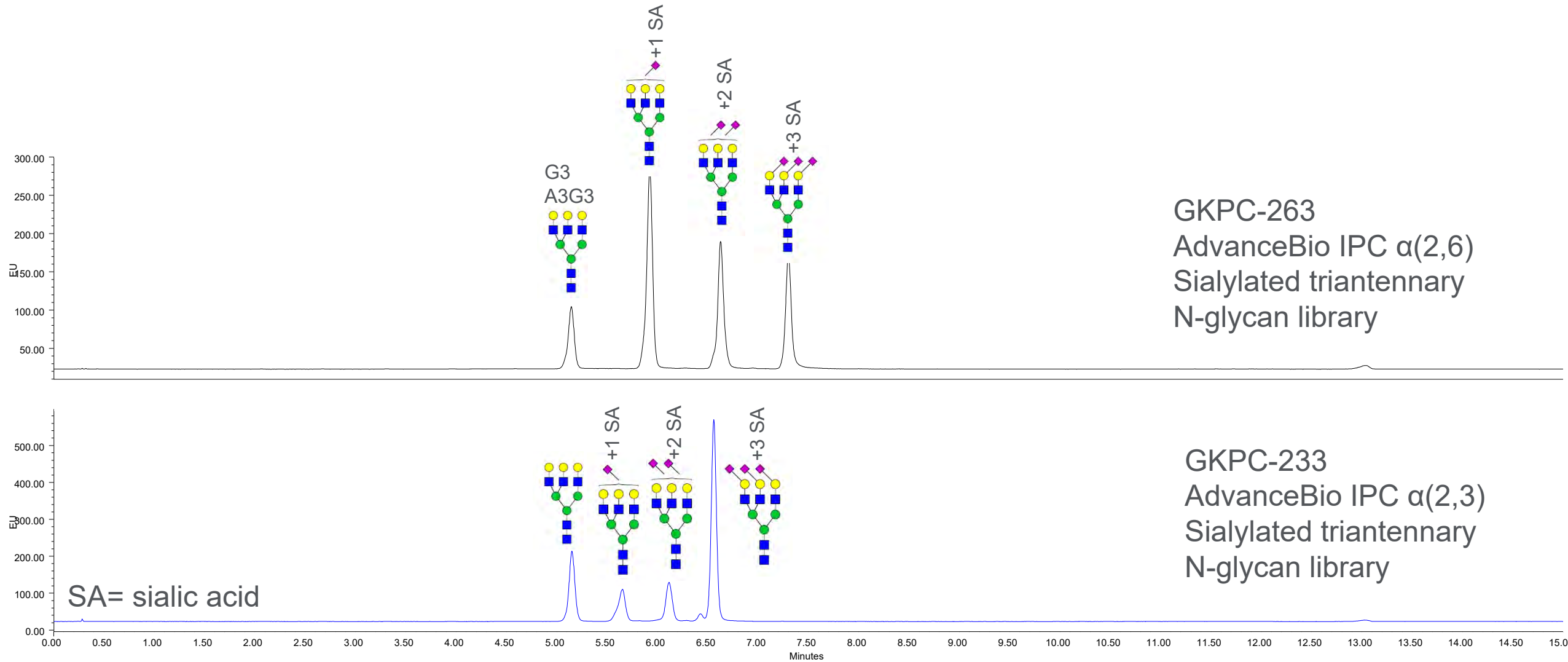

Shorter retention of  $\alpha(2,3)$ -sialylated N-glycans vs  $\alpha(2,6)$  sialic acid linkages

# InstantPC Sialylated Tetraantennary N-Glycan Library

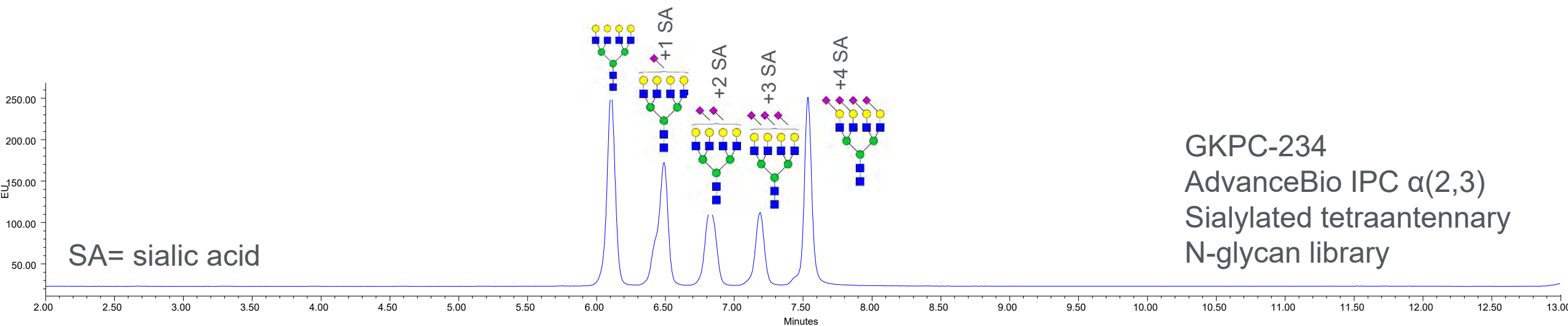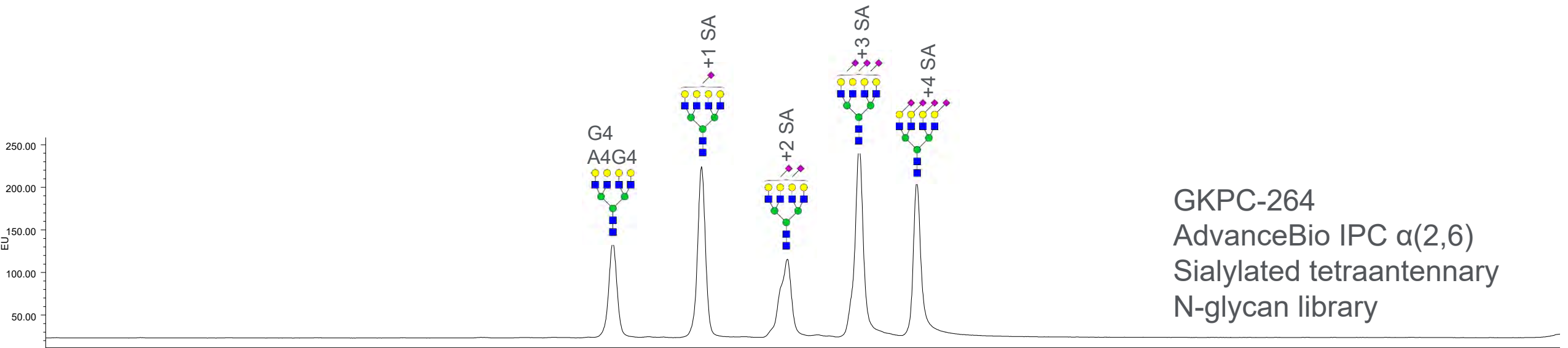

Shorter retention of  $\alpha(2,3)$ -sialylated N-glycans vs  $\alpha(2,6)$  sialic acid linkages

# Additional Resources for InstantPC N-Glycan LC/FLD & MS Analysis

## **Streamlined Workflows for N-Glycan Analysis of Biotherapeutics Using InstantPC with LC/FLD/MS**

*John Yan, Andres Guerrero, Ace G. Galermo, Ted Haxo, Sergey Vlasenko, Justin Hyche, Tom Rice and Aled Jones*

ASMS 2019 poster [ThP697](#)

## **A Comprehensive Approach for Monoclonal Antibody N-linked Glycan Analysis from Sample Preparation to Data Analysis**

*David Wong, Oscar Potter, Jordy Hsaio, Te-Wei Chu*

Technical Note [5991-8550EN](#)

## **Analysis of Monoclonal Antibody N-glycans by Fluorescence Detection and Robust Mass Selective Detection Using the Agilent LC/MSD XT**

*Oscar Potter, Gregory Staples, Jordy Hsaio, Te-Wei Chu*

Technical Note [5991-8071EN](#)

## **Comparison of Relative Quantification of Monoclonal Antibody N-Glycans Using Fluorescence and MS Detection**

*Oscar Potter and Greg Staples, Agilent Technologies*

Technical Note [5991-6958EN](#)

## **Comparison of Common Fluorescent Labels for LC/MS Analysis of Released N-Linked Glycans**

*John Yan, Andres Guerrero, Steven Mast, Ted Haxo, Aled Jones*

Technical Note [5994-0942EN](#)

## **Automation of Gly-X N-Glycan Sample Prep with InstantPC and InstantQ Dyes**

*Loredana Serafini\*, Ted Haxo†, Emily Dale†, Adele Taylor†, Katherine M. Brenda\* (\*Gilead, †ProZyme)*


ProZyme Technical Note [\(copy available upon request\)](#)

## **Development of an Instant Glycan Labeling Dye for High Throughput Analysis by Mass Spectrometry**

*Michael Kimzey, Zoltan Szabo, Vaishali Sharma, Alexander Gyenes, Samnang Tep, Adele Taylor, Aled Jones, Justin Hyche, Ted Haxo, Sergey Vlasenko*

ProZyme Technical Note [\(copy available upon request\)](#)

## 2-AB Or Not 2-AB.....

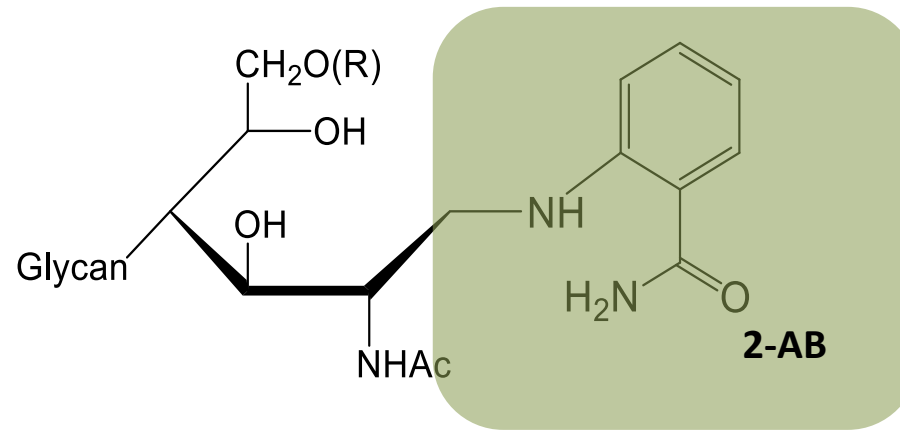

Label makes N-glycan less polar

Traditionally used for HILIC separations of N-glycans with fluorescence detection

Lower FLD and MS signal than InstantPC

2-AB (2-aminobenzamide) dye has been used to generate released N-glycan data for more than 20 years






Well established in many laboratories

May need to match historic data using 2-AB

# Gly-X 2-AB Express Workflow

1-40  $\mu\text{g}$   
0.05-2 mg/ml

Reductive  
Amination

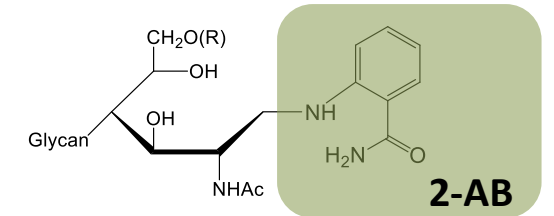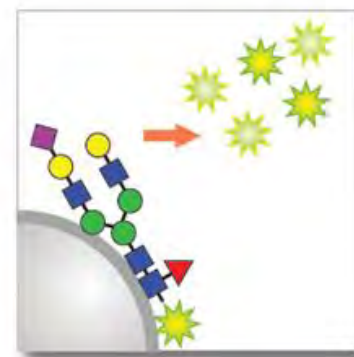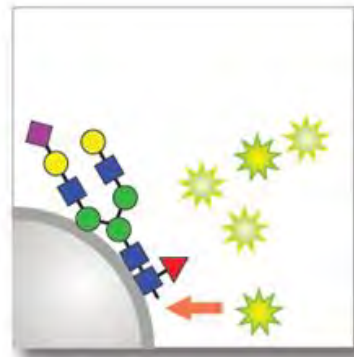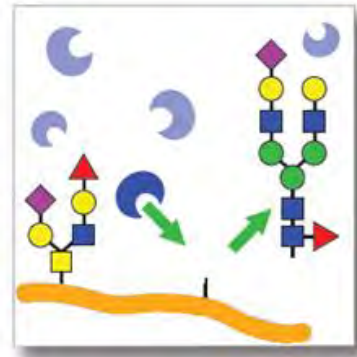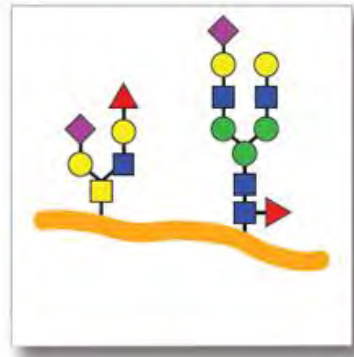


Denature  
3 min at 90°C

Deglycosylate  
5 min at 50°C

Label  
60 min at 65°C  
**No Dry Down**

Clean up  
15-20 min at  
RT

Labeled N-Glycans

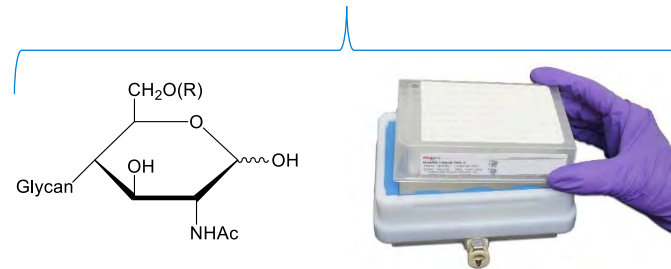


Workflow time: 120 min

# Rituximab Gly-X 2-AB Express N-Glycans

## Analytical Method:

UHPLC-HILIC, 60 minute method, 1 µl injection (equivalent to glycans from 0.4 µg protein), FLD: 260/430 nm

AdvanceBio Glycan Mapping column, 2.1 x 150 mm, 1.8 µm

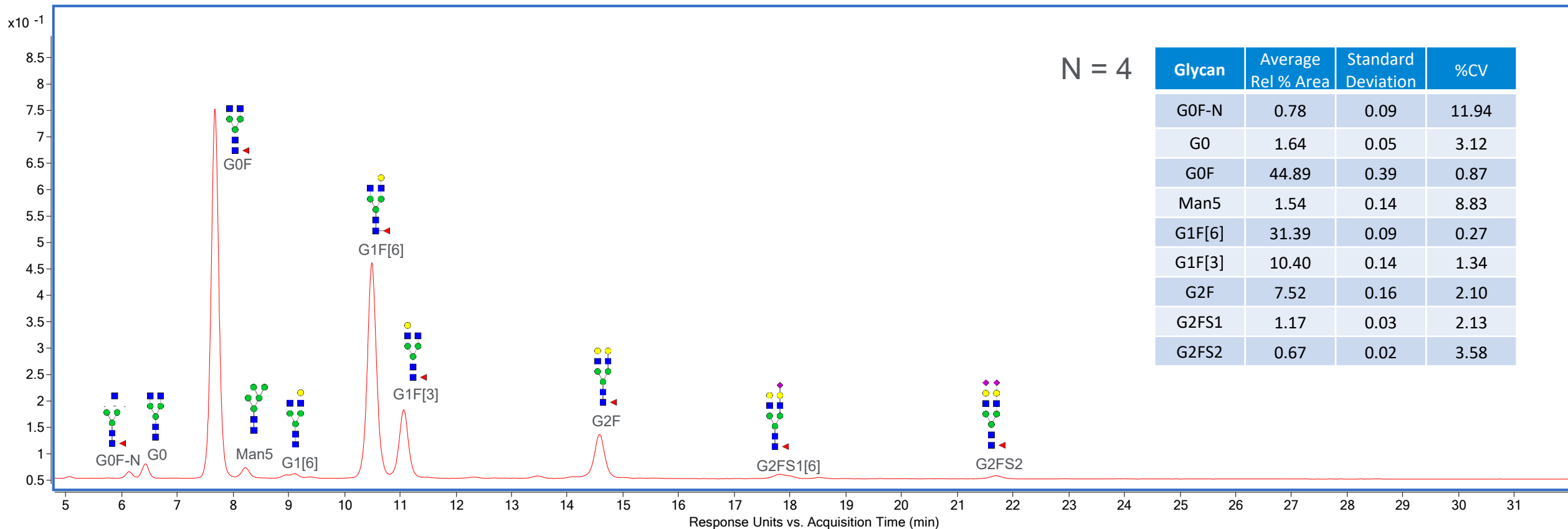

# Gly-X 2-AB vs InstantPC - rituximab



N=4

2-AB

InstantPC


	Average Rel % Area	Standard Deviation	%CV
G0F-N	0.78	0.09	11.94
G0	1.64	0.05	3.12
G0F	44.89	0.39	0.87
Man5	1.54	0.14	8.83
G1F[6]	31.39	0.09	0.27
G1F[3]	10.40	0.14	1.34
G2F	7.52	0.16	2.10
G2FS1	1.17	0.03	2.13
G2FS2	0.67	0.02	3.58

	Average Rel % Area	Standard Deviation	%CV
G0F-N	0.75	0.01	1.55
G0	1.47	0.02	1.18
G0F	46.82	0.07	0.15
Man5	1.21	0.01	0.83
G1[6]	0.75	0.02	2.67
G1F[6]	31.21	0.11	0.35
G1F[3]	9.27	0.05	0.54
G2F	7.04	0.04	0.51
G2FS1[6]	0.67	0.02	2.29
G2FS1[3]	0.37	0.06	15.98
G2FS2	0.45	0.03	6.67

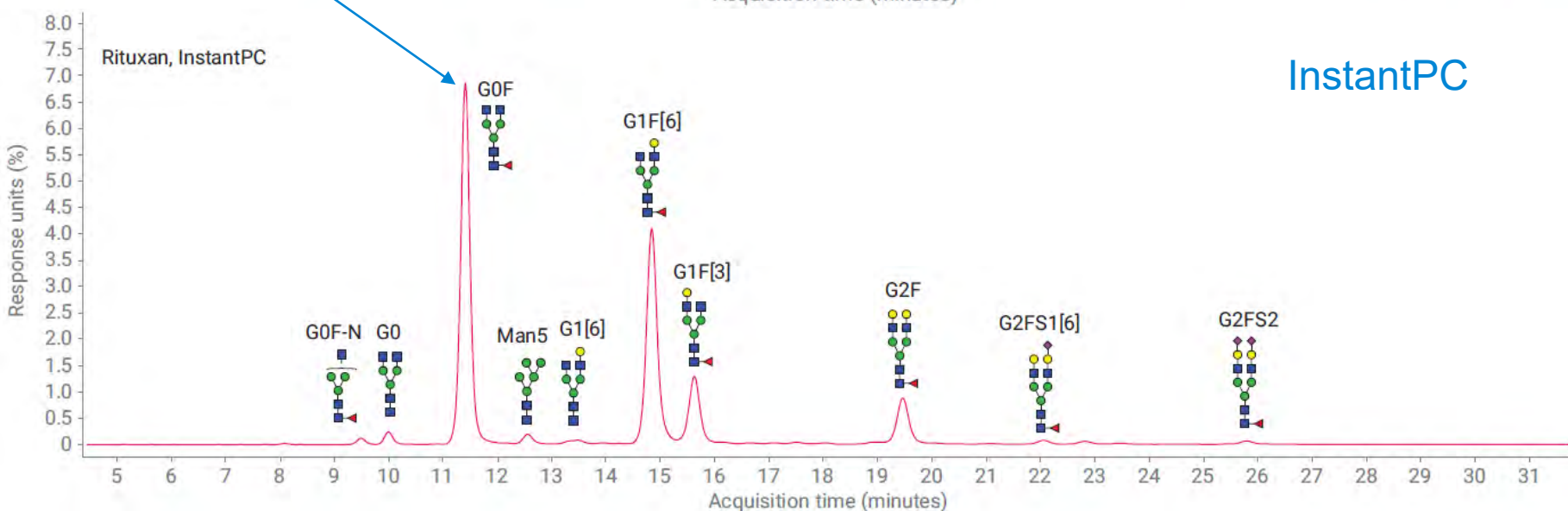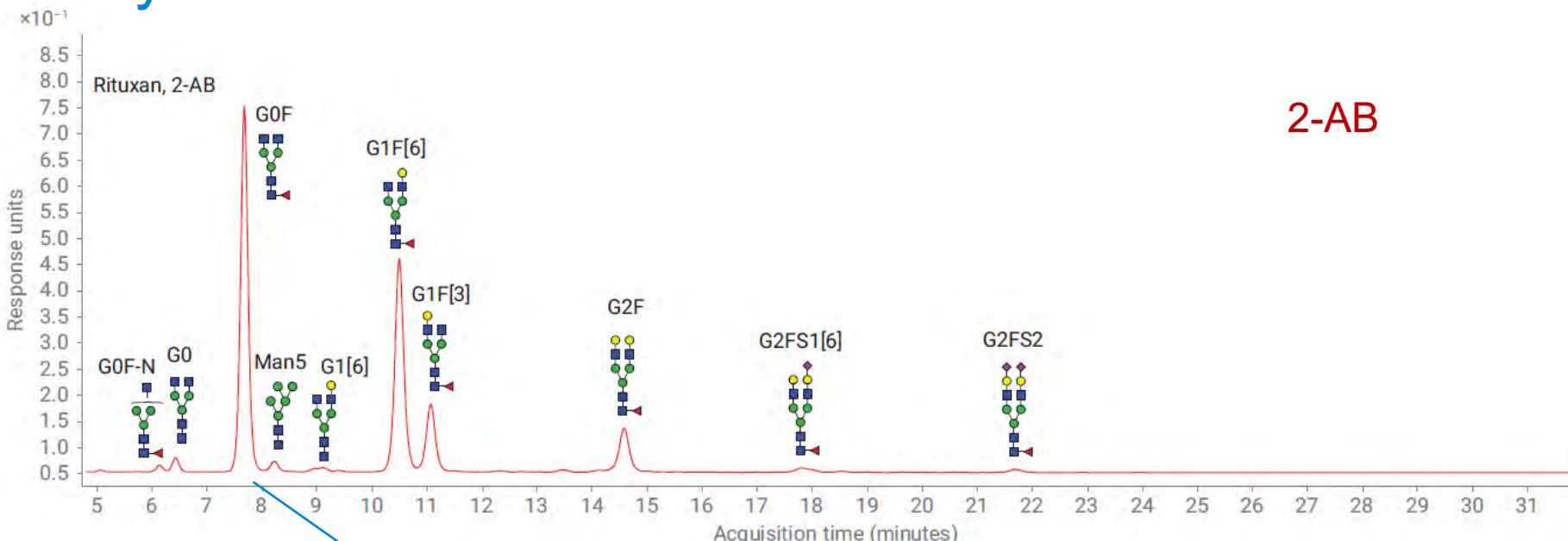


# Gly-X 2-AB vs InstantPC - etanercept

N=4

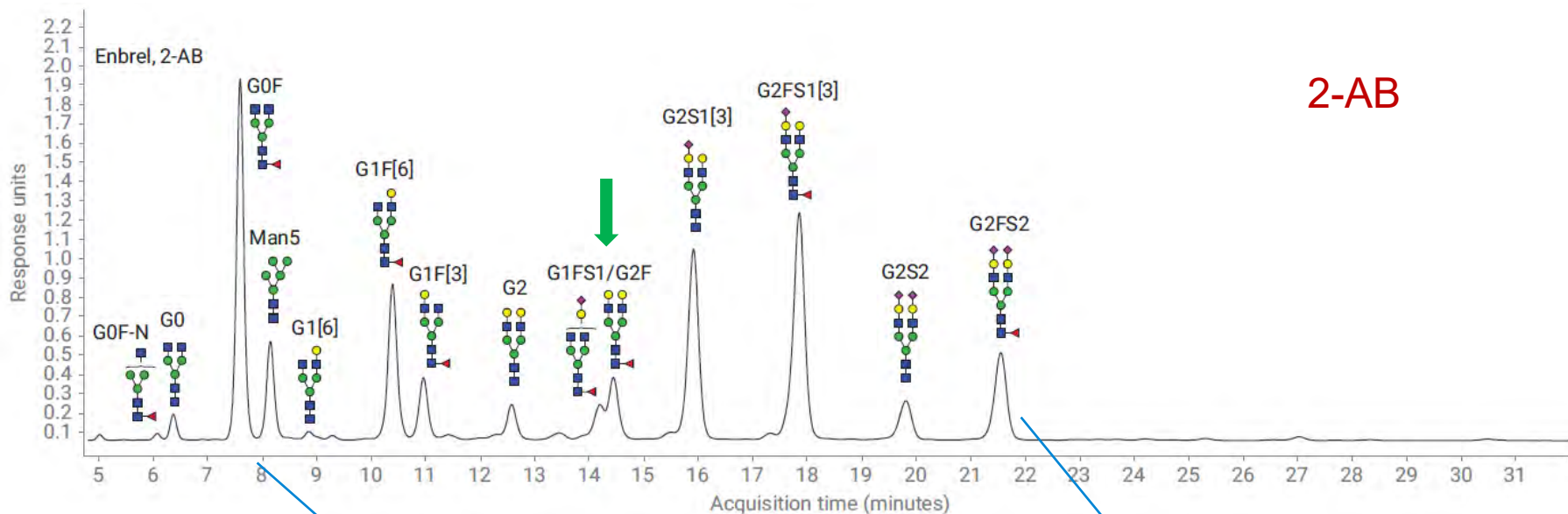

2-AB

	Average Rel % Area	Standard Deviation	%CV
G0F-N	0.32	0.02	7.44
G0	1.27	0.07	5.34
G0F	20.18	0.45	2.22
Man5	5.50	0.34	6.17
G1[6]	0.45	0.02	3.89
G1F[6]	10.35	0.33	3.18
G1F[3]	3.92	0.17	4.39
G2	2.21	0.15	6.78
G2F/G1FS1	7.00	0.25	3.63
G2S1	15.19	0.17	1.09
G2FS1	20.10	0.32	1.59
G2S2	4.19	0.25	5.95
G2SF2	9.35	0.74	7.93




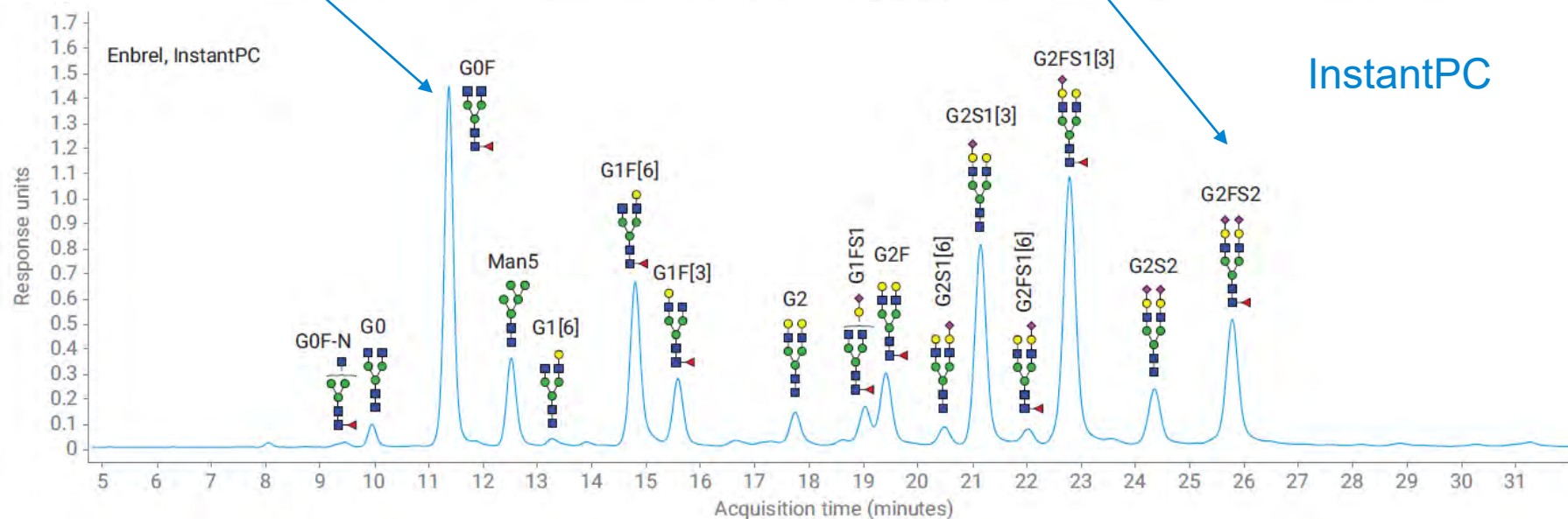


InstantPC

	Average Rel % Area	Standard Deviation	%CV
G0	1.10	0.02	2.09
G0F	19.36	0.16	0.84
Man5	5.08	0.03	0.52
G1[6]	0.48	0.00	0.00
G1F[6]	10.48	0.04	0.39
G1F[3]	3.97	0.01	0.25
G2	2.08	0.01	0.55
G1FS1	1.84	0.05	2.49
G2F	4.26	0.09	1.99
G2S1[6]	1.18	0.01	0.49
G2S1[3]	13.91	0.04	0.31
G2FS1[6]	0.89	0.00	0.00
G2FS1[3]	20.54	0.08	0.37
G2S2	4.26	0.01	0.14
G2FS2	10.54	0.08	0.78





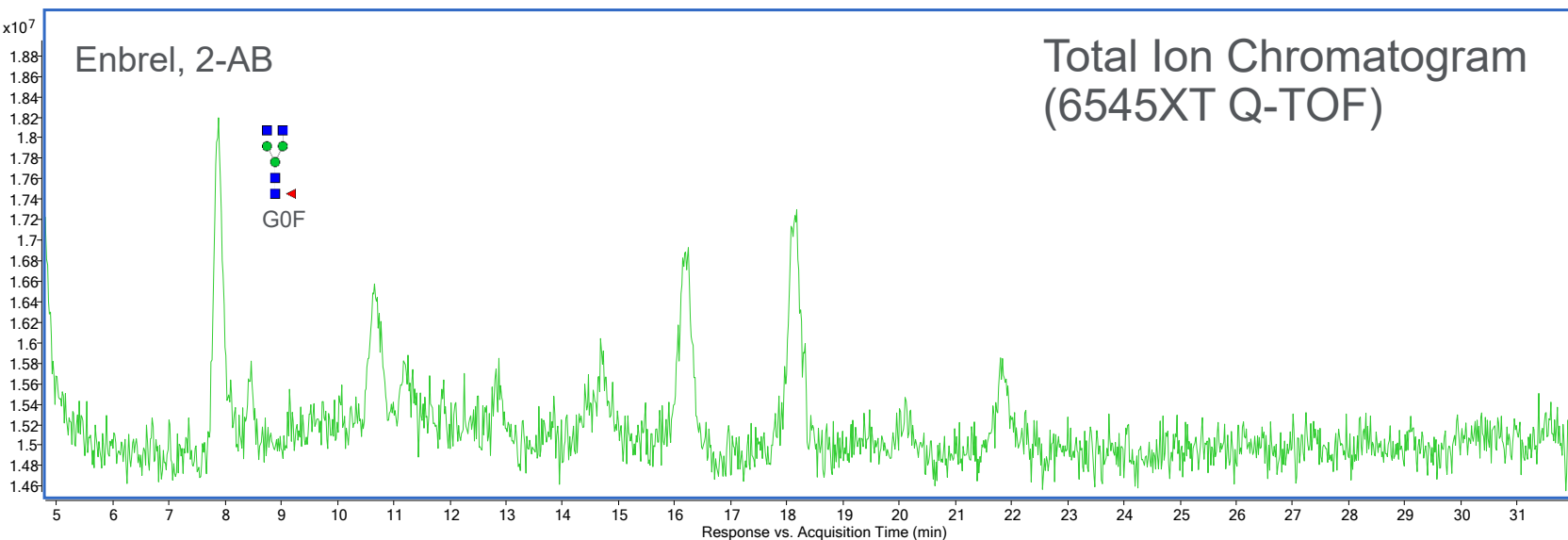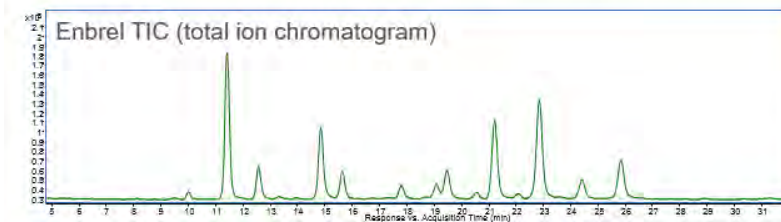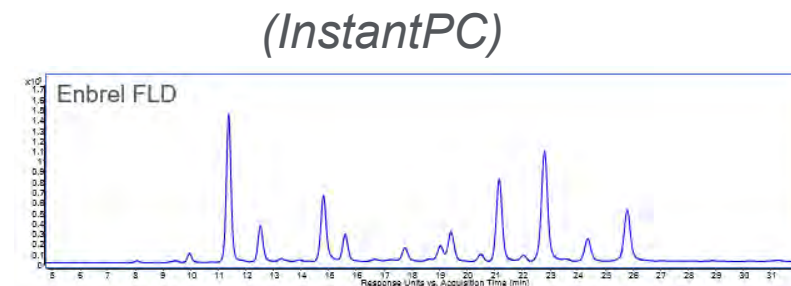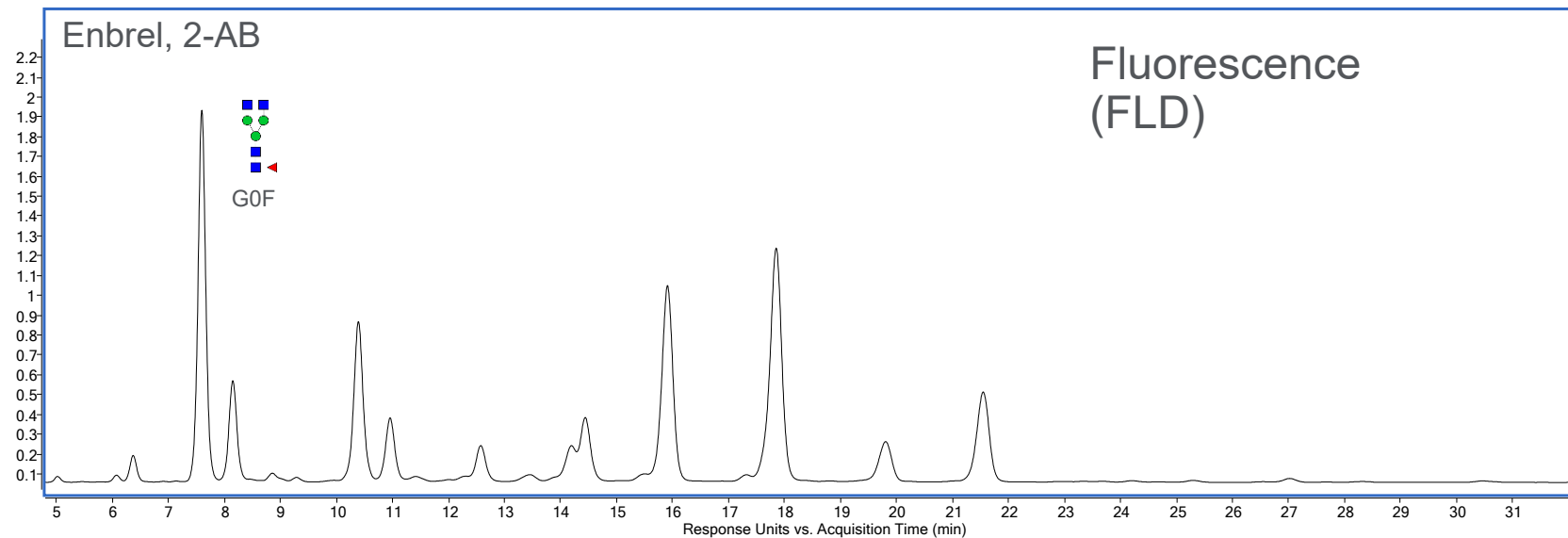





























# 2-AB N-Glycans: MS



# 2-AB N-Glycan Standards & Libraries

## Individual N-Glycans


Product Code	Description
GKSB-401	G0-N 
GKSB-301	G0 
GKSB-402	G0F-N 
GKSB-302	G0F 
GKSB-303	G0FB 
GKSB-317	G1 
GKSB-316	G1F 
GKSB-304	G2 
GKSB-305	G2F 
GKSB-306	G2FB 
GKSB-318	NA2Ga2F 
GKSB-311	A1 
GKSB-315	A1F 
GKSB-312	A2 
GKSB-313	A2F 
GKSB-307	NGA3 

Product Code	Description
GKSB-308	G3 
GKSB-314	A3 
GKSB-309	NGA4 
GKSB-310	G4 
GKSB-111	HYBR 
GKSB-100	NN 
GKSB-101	Man3 
GKSB-102	Man3F 
GKSB-103	Man5 
GKSB-104	Man6 
GKSB-105	Man7 
GKSB-106	Man8 
GKSB-107	Man9 
GKSB-201	GalGalNAc 
GKSB-203	3'-SLN 
GKSB-204	6'-SLN 

## N-Glycan Libraries

Product Code	Description
GKSB-005	Human IgG N-Linked Glycan Library
GKSB-520	Biantennary & High Mannose Partitioned Library
GKSB-001	Human $\alpha$ 1-acid Glycoprotein N-Linked Glycan Library
GKSB-002	Bovine Fetuin N-linked Glycan Library
GKSB-503	Glucose Homopolymer Standard
GKSB-232	$\alpha$ (2-3) Sialylated Biantennary Library
GKSB-262	$\alpha$ (2-6) Sialylated Biantennary Library
GKSB-233	$\alpha$ (2-3) Sialylated Triantennary Library
GKSB-263	$\alpha$ (2-6) Sialylated Triantennary Library
GKSB-234	$\alpha$ (2-3) Sialylated Tetraantennary Library
GKSB-264	$\alpha$ (2-6) Sialylated Tetraantennary Library

# 2-AB Sialylated Triantennary N-Glycan Library

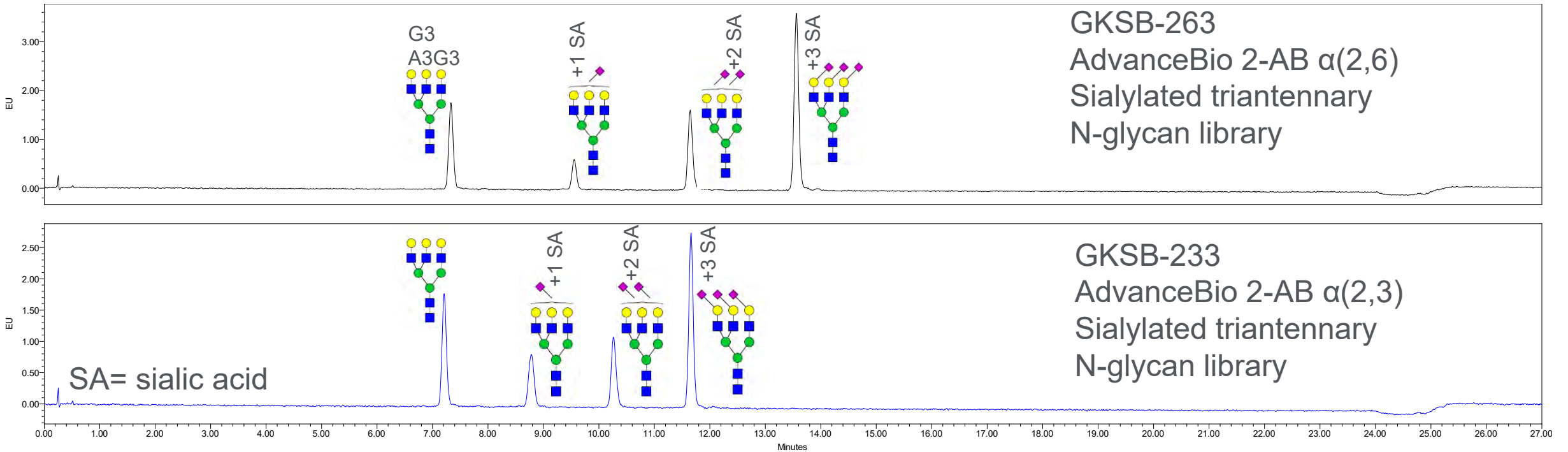


Shorter retention of  $\alpha(2,3)$ -sialylated N-glycans vs  $\alpha(2,6)$  sialic acid linkages

# Importance of Sialic Acid on Biotherapeutics


## Terminal sialic acid

↓ clearance →

↑ anti-inflammatory →

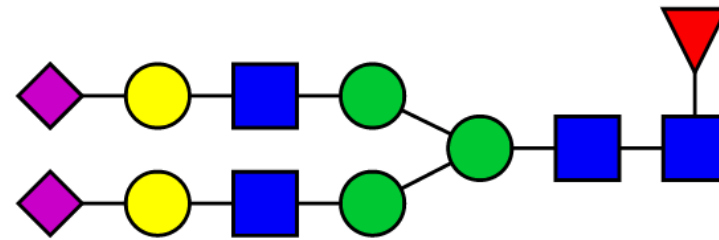

NeuGc  
NGNA →

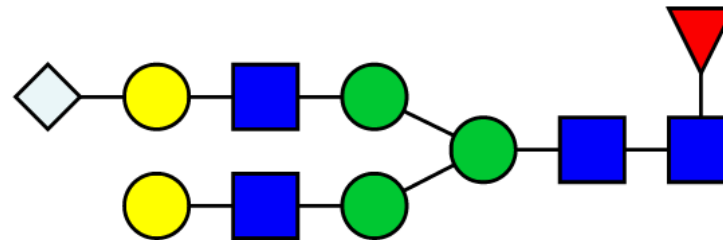


- Galactose
- Mannose
- ▲ Fucose
- N-Acetylglucosamine
- ◆ N-Acetylneuraminic acid
- ◇ N-Glycolylneuraminic acid

N-glycolylneuraminic acid (NeuGc/NGNA): Potentially immunogenic


# Total Sialic Acid Quantitation

- Plate-based assay (GS48-SAQ, GS96-SAQ)
- Rapid quantitation of total sialic acid released from intact proteins by sialidase A
- Fluorescence plate reader (absorbance an option, lower sensitivity)
- Broad range of detection: 40 – 1,000 pmol fluorescence, 500 – 4,000 pmol absorbance


N-Acetylneuraminic aldolase catalyzes the reversible reaction:

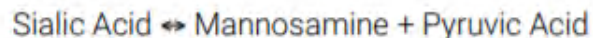

Then pyruvate oxidase catalyzes the reaction:

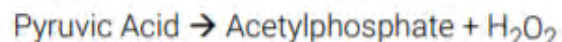

Finally,  $\text{H}_2\text{O}_2$  forms a 1:1 complex with the Dye to form a fluorescent Reporter Dye that may be read by fluorescence or absorbance detection.

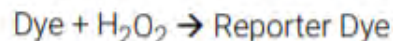

**10  $\mu\text{L}$  Glycoprotein**

30 min Sialidase A Digest



**Released Sialic Acid**

60 min Conversion and Developing



**Read the Plate**

Ex 530 nm / Em 590 nm (FLD)  
530 nm (Absorbance)

# Total Sialic Acid Quantitation: Starting Concentrations and Amounts of Glycoprotein

Glycoprotein	Concentration (mg/mL)	Sample volume (μL)	Sample mass (μg)	MW (kDa)	pmol protein
Fetuin	0.25	10	2.5	48	52
MabThera	10	10*	100	145	690
Enbrel	0.25	10	2.5	150	16.7
Zaltrap	0.5	10	5	115	43
Orencia	0.5	10	5	92	54
EPO alfa	1	10	10	30.4	329

*\*For glycoproteins with lower sialylation such as monoclonal antibodies, up to 30 μL sample may be used with the kit.*

# Intra-Assay Repeatability

- Single operator, 3 replicates per sample
- SA - sialic acid

	Fetuin	MabThera	Enbrel	Zaltrap	Orencia	EPO alfa
Starting concentration (mg/mL)	0.25	10	0.25	0.5	0.5	1
MW (kDa)	48	145	150	115	92	30.4
nmol SA/10 $\mu$ L sample	0.58	0.08	0.55	0.45	0.59	0.16
nmol SA/mg protein	232.53	0.79	220.93	90.27	117.27	15.50
mol SA/mol protein	11.16	0.11	33.14	10.38	10.79	0.47
%CV (n=3)	0.55%	2.65%	1.81%	0.46%	0.60%	3.59%
SD (mol SA/mol protein)	0.06	0.00	0.60	0.05	0.06	0.02



# Operator to Operator Repeatability

- 6 operators
- Samples: Fetuin, Enbrel, Orencia, Zaltrap, MabThera, EPO alfa
- 3 replicates per sample


	Fetuin	MabThera	Enbrel	Zaltrap	Orencia	EPO alfa
Operator 1	11.35	0.18	33.24	11.41	10.56	0.45
Operator 2	11.16	0.11	33.14	10.38	10.79	0.47
Operator 3	12.07	0.18	35.58	12.65	11.52	0.49
Operator 4	10.34	0.17	33.76	10.92	11.17	0.49
Operator 5	12.08	0.16	33.20	9.91	9.50	0.46
Operator 6	10.44	0.17	31.66	9.74	10.32	0.45
Average mol SA/mol protein	11.24	0.16	33.43	10.83	10.64	0.47
%CV (n=6)	6.72%	15.28%	3.79%	10.03%	6.63%	3.86%
SD	0.76	0.02	0.02	0.71	1.09	1.27




# 100 $\mu$ M Sialic Acid Standard Lot Comparison (target 500 pmol)





- WS0182: older NANA standard included with GS300/GF57
- WS0377: new NANA standard for GS48/96-SAQ, uses USP material

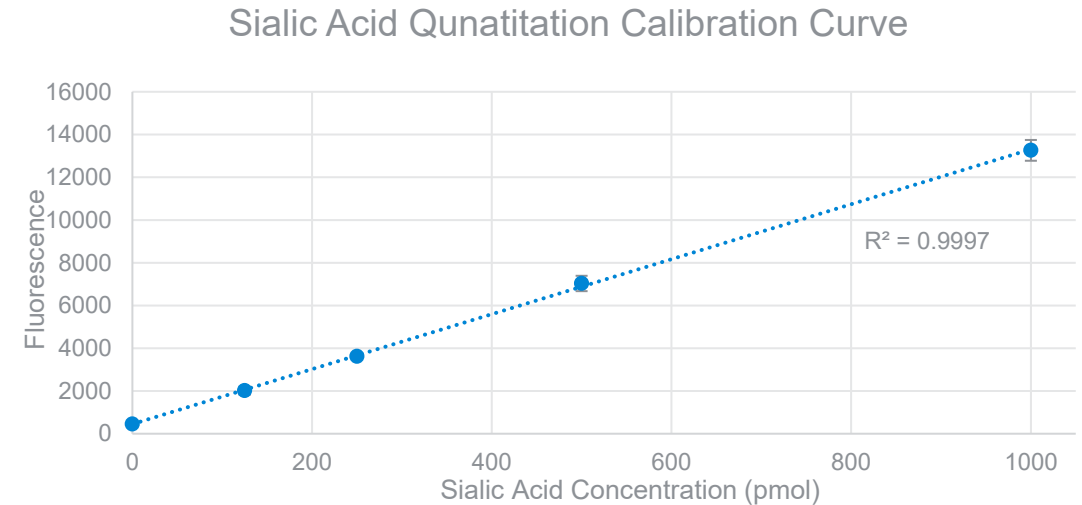



# Neu5Ac and Neu5Gc

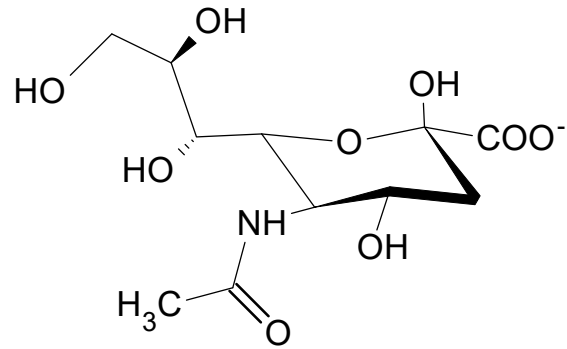

Neu5Ac  
NANA

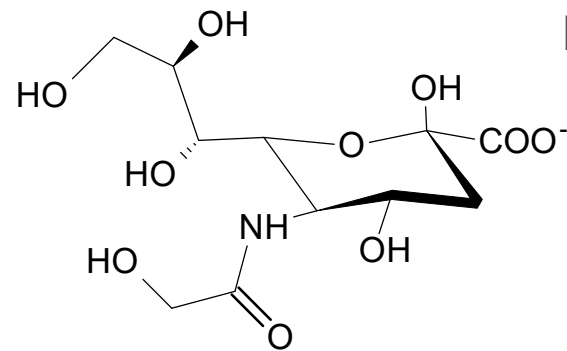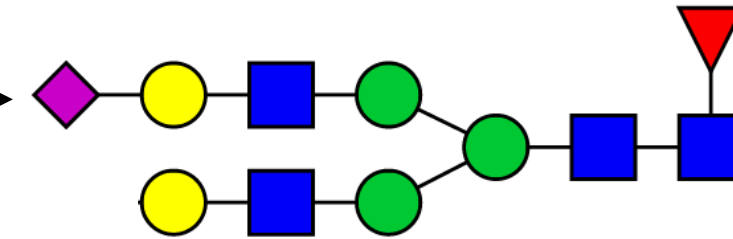

Neu5Gc  
NGNA

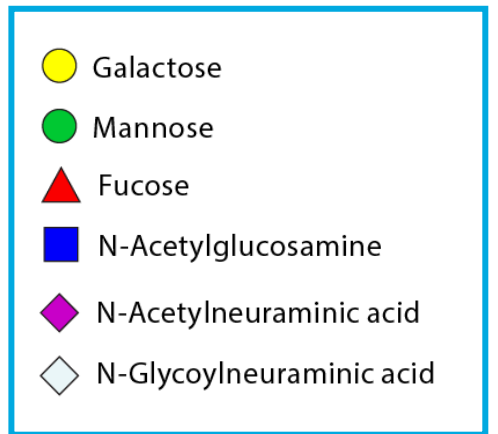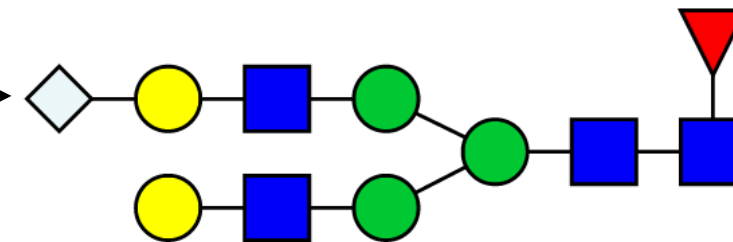


Potentially immunogenic


Ghaderi et al., *Biotechnol Genet Eng Rev.* 2012; 28:147-75

Jones, *BioPharm Intl.* 2017; 30:20-25


# DMB Labeling for Profiling and Quantitation of Sialic Acid

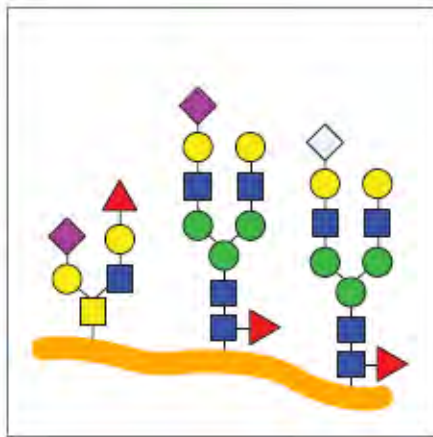


Sialylated  
glycoprotein

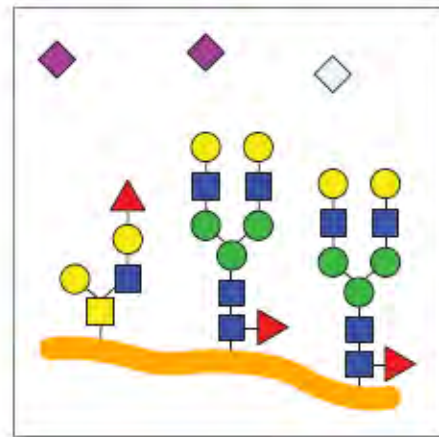

Sialic acid  
release

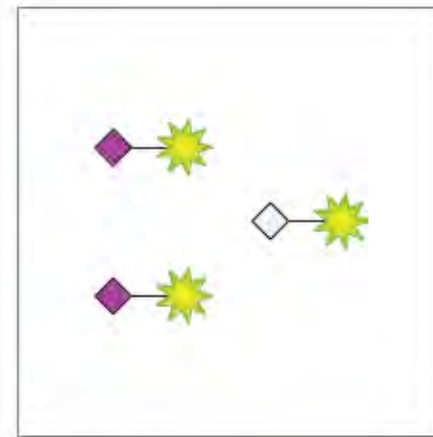


DMB labeling

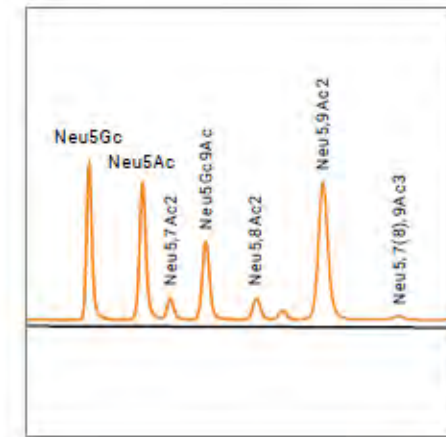

RPLC separation  
FLD or MS detection  
Quantitation

# DMB Labeled Sialic Acid Reference Panel (SARP)

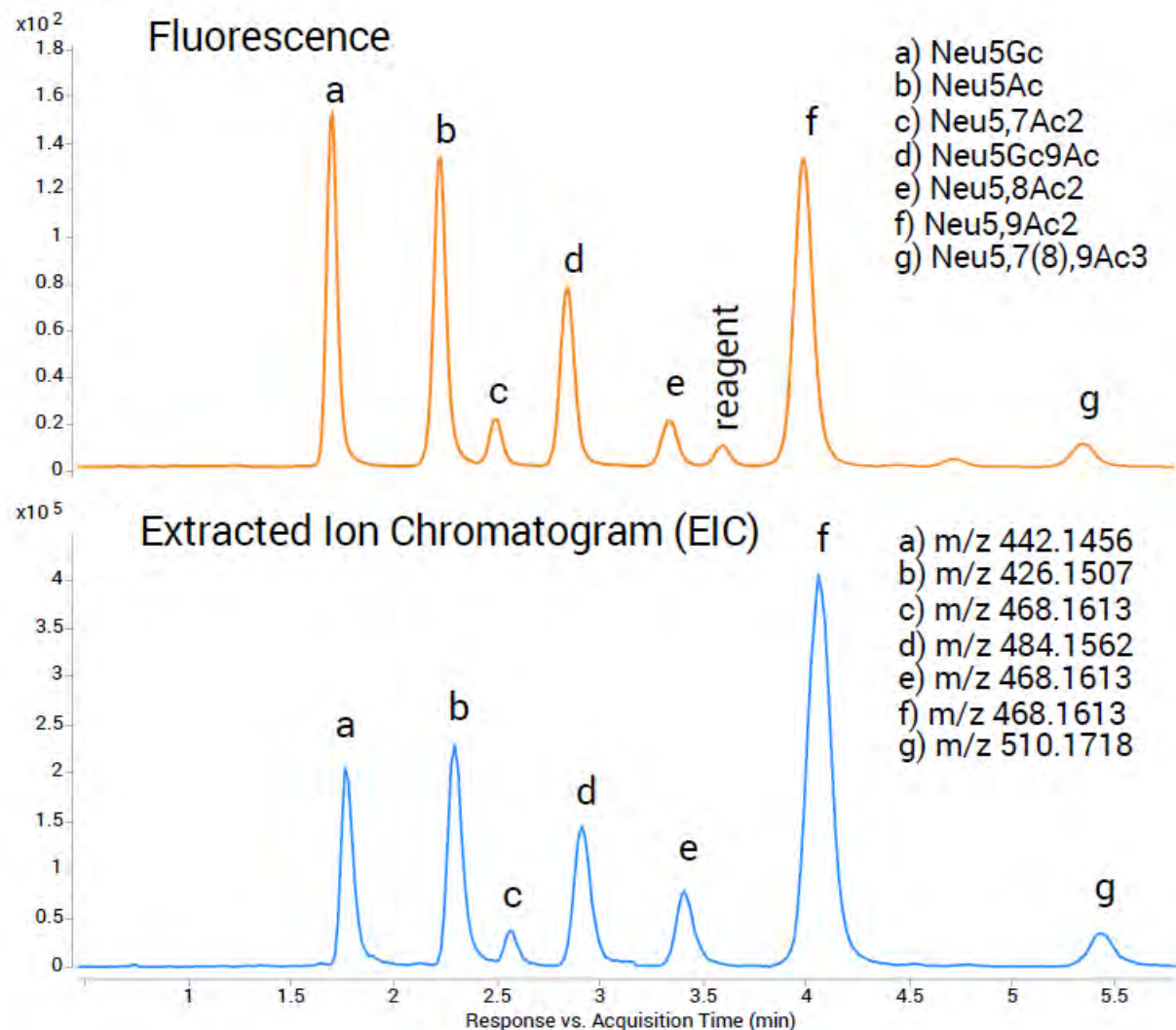


Parameter	Value	6545XT Q-TOF																																				
Instrument	Agilent 1290 Infinity II LC System	Source	Dual AJS ESI																																			
Column	Agilent InfinityLab PoroShell 120 EC-C18, 2.1 x 75 mm, 2.7 μm (p/n 697775-902)	Gas Temperature	350 °C																																			
Column Temp	30 °C	Drying Gas Flow	11 L/min																																			
Mobile Phase	A) methanol: acetonitrile: water (4:8:88) B) acetonitrile	Nebulizer	15 psi																																			
Gradient Program	<table border="1"> <thead> <tr> <th>Time (min)</th> <th>%A</th> <th>%B</th> <th>Flow rate (mL/min)</th> <th></th> </tr> </thead> <tbody> <tr> <td>0.00</td> <td>100</td> <td>0</td> <td>0.4</td> <td>Isocratic</td> </tr> <tr> <td>6.00</td> <td>100</td> <td>0</td> <td>0.4</td> <td>Elution</td> </tr> <tr> <td>6.25</td> <td>20</td> <td>80</td> <td>0.4</td> <td></td> </tr> <tr> <td>7.30</td> <td>20</td> <td>80</td> <td>0.4</td> <td>Wash</td> </tr> <tr> <td>7.50</td> <td>100</td> <td>0</td> <td>0.4</td> <td>Re-equilibration</td> </tr> <tr> <td>10.00</td> <td>100</td> <td>0</td> <td>0.4</td> <td></td> </tr> </tbody> </table>	Time (min)	%A	%B	Flow rate (mL/min)		0.00	100	0	0.4	Isocratic	6.00	100	0	0.4	Elution	6.25	20	80	0.4		7.30	20	80	0.4	Wash	7.50	100	0	0.4	Re-equilibration	10.00	100	0	0.4		Sheath Gas Temperature	400 °C
Time (min)	%A	%B	Flow rate (mL/min)																																			
0.00	100	0	0.4	Isocratic																																		
6.00	100	0	0.4	Elution																																		
6.25	20	80	0.4																																			
7.30	20	80	0.4	Wash																																		
7.50	100	0	0.4	Re-equilibration																																		
10.00	100	0	0.4																																			
Injection Volume	10 μL	Sheath Gas Flow	12 L/min																																			
Detection	Agilent 1260 Infinity II FLD λ <sub>Ex</sub> 373 nm, λ <sub>Em</sub> 448 nm	Vcap	1400 V																																			
		Nozzle Voltage	2800 V																																			
		Fragmentor	120 V																																			
		Skimmer	65 V																																			
		Mass Range	m/z 400-1000																																			
		Scan Rate	1 spectra/sec																																			
		Acquisition Mode	High resolution (4 GHz)																																			

Pressure > 400 bar

Extracted ion chromatogram of DMB labeled sialic acid species, [M+H]<sup>1+</sup>

# DMB Labeled Sialic Acids of Biotherapeutics & NISTmAb

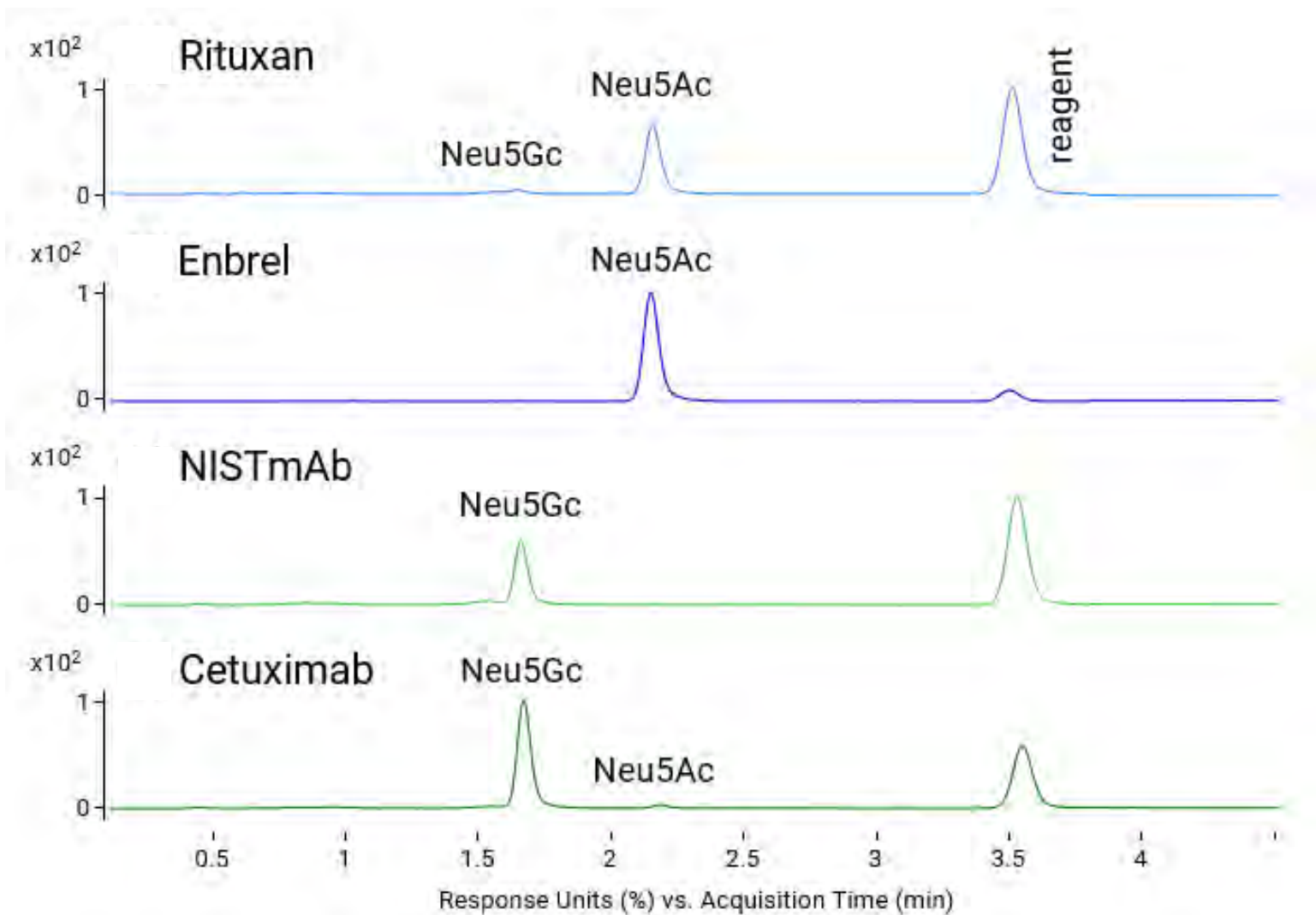

NISTmAb – available from Agilent


- 25  $\mu$ L (5191-5744)
- 4 x 25  $\mu$ L (5191-5745)

RP UHPLC (Poroshell EC-C18)


Fluorescence detection

# Mass spectra of DMB NeuAc and NeuGc

Enbrel

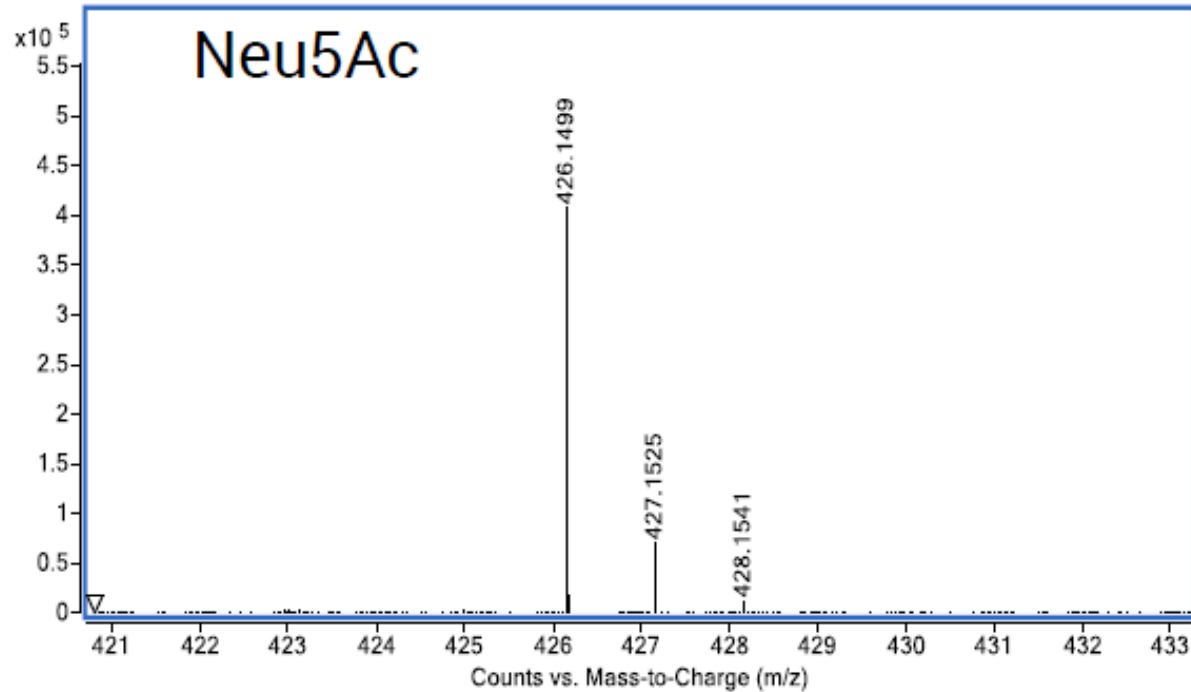

Cetuximab

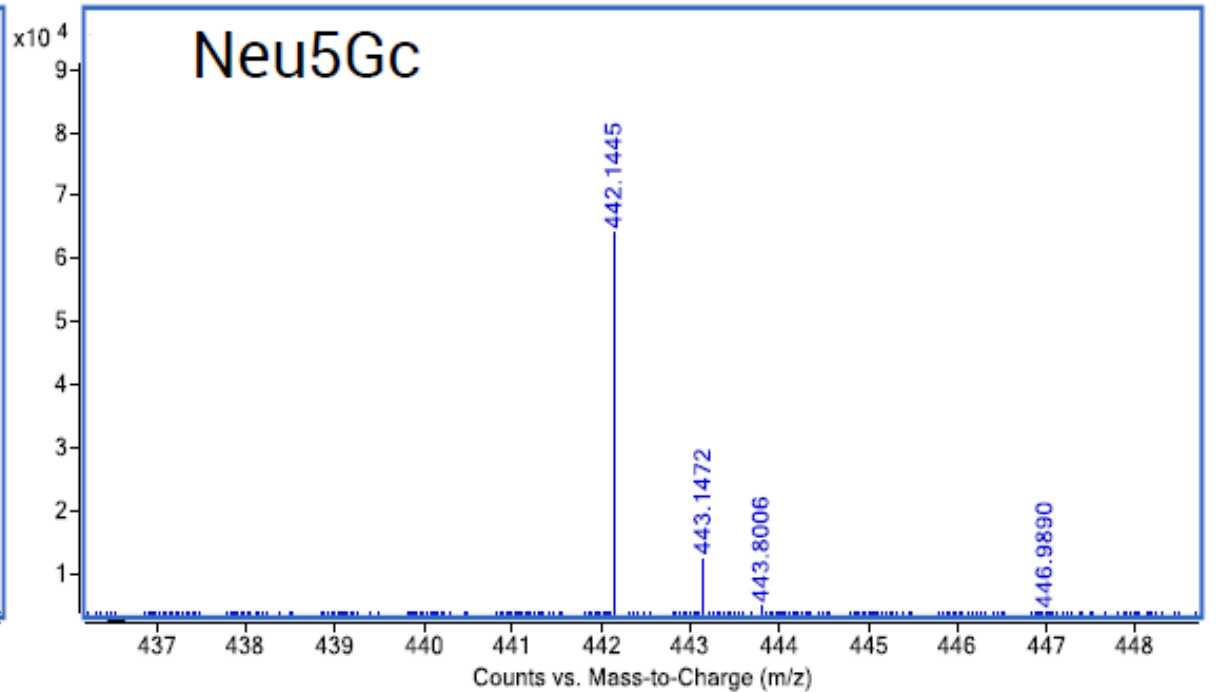

# Summary

Gly-X: rapid, simple, automatable sample preparation for N-glycan analysis

Choice of dyes:

- InstantPC (LC/FLD/MS): >1 hour
- 2-AB Express (LC/FLD/MS): 2 hours

Fluorescence detection for relative % area quantitation

MS detection for the assignment of glycan structures to peaks

Sialylated glycan standards

Sialic acid:

- Total sialic acid quantitation (SAQ), plate based
- Sialic acid profiling (DMB), LC/FLD/MS

# Collaboration

Talk to us about:

- N-Glycan sample prep
- Glycan standards
- Endoglycosidases & Exoglycosidases
- Sialic acid quantitation
- Streptavidin and phycobiliproteins
- Analytical services

Product demos

Beta testing

New technologies



# Acknowledgements

## Agilent Technologies

John Yan

John Barich

Hamutal Bonen

Archana Datt

Anna Fong

Ace G. Galermo

Ted Haxo

Justin Hyché

Aled Jones

Tom Mannion

Steve Mast

Beth Morgan

Shiva Pourkaveh

Marco Rhotert

Jim Torrence

Sergey Vlasenko

Anne Blackwell

Linda Lloyd

Oscar Potter

Suma Ramagiri

Greg Staples

