


Assembly Instruction

Table Stand
For Stereo Microscope

Model A605

MicroscopeNet.com

1. Parts identification

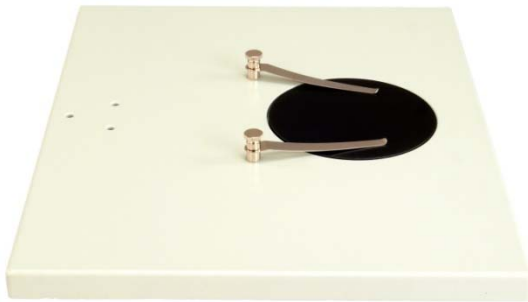

Stand Base



Pillar




Focus Block



Small Parts

2. Assembly instructions

- 1) Turn the base upside down and insert the 3 screws through the holes in the base with the washers on.

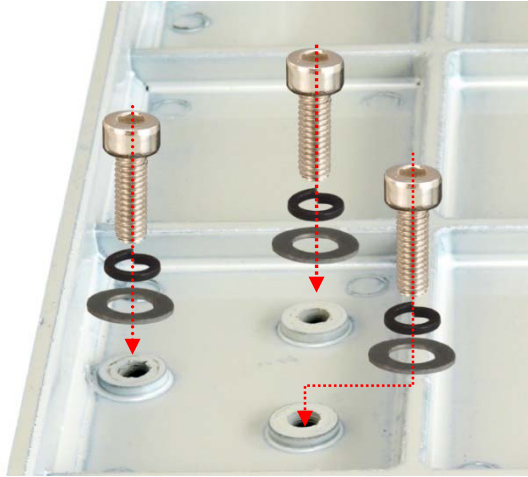

- 2) Align the 3 holes in the bottom of the pillar with the 3 holes in the base and tighten the 3 screws with the Allen key.



3) Slip the focus block through pillar and let it down against the collar.



4) Install the knob onto the focus block.

