

User Manual

Digital Compound Binocular LED Microscope

Model MD82EE10



MicroscopeNet.com

Table of Contents

i. Caution	1
ii. Care and Maintenance.....	2
1. Components Illustration	3
2. Installation.....	4
3. Operation.....	4
4. Specifications.....	7
5. Troubleshooting Guide.....	8


i. Caution

1. Open the carton carefully with a knife or paper cutter. Find the “UP” sign and place the Styrofoam container on the side that makes the arrow upward. If the “UP” sign is missing, please open the Styrofoam container gently to prevent any accessory, i.e. objectives or eyepieces, from dropping and being damaged.
2. Do not discard the molded Styrofoam container; the container should be retained should the microscope ever requires reshipment.
3. Keep the instrument out of direct sunlight, high temperature or humidity, and dusty environments. Ensure the microscope is located on a smooth, level and firm surface.
4. If any specimen solutions or other liquids splash onto the stage, objective or any other component, disconnect the power cord immediately and wipe up the spillage. Otherwise, the instrument may be damaged.
5. All electrical connectors (power cord) should be inserted into an electrical surge suppressor to prevent damage due to voltage fluctuations.
6. Confirm that the input voltage indicated on your microscope corresponds to your line voltage. The use of a different input voltage other than that as indicated will cause severe damage to the microscope.
7. **Note:** please read the instruction of the operation of camera in manual **3.8** below and the CD in the package before you start to use it.

ii. Care and Maintenance

1. Do not attempt to disassemble any component including eyepieces, objectives or focusing assembly.
2. Keep the instrument clean; remove dirt and debris regularly. Accumulated dirt on metal surfaces should be cleaned with a damp cloth. More persistent dirt should be removed using a mild soap solution. **Do not use organic solvents for cleansing.**
3. The outer surface of the optics should be inspected and cleaned periodically using an air stream from an air bulb. If dirt remains on the optical surface, use a soft cloth or cotton swab dampened with a lens cleaning solution (available at camera stores). All optical lenses should be swabbed using a circular motion. A small amount of absorbent cotton wound on the end of a tapered stick makes a useful tool for cleaning recessed optical surfaces. Avoid using an excessive amount of solvents as this may cause problems with optical coatings or cemented optics or the flowing solvent may pick up grease making cleaning more difficult. Oil immersion objectives should be cleaned immediately after use by removing the oil with lens tissue or a clean, soft cloth.
4. Store the instrument in a cool, dry environment. Cover the microscope with the dust cover when not in use.

1 Components Illustration

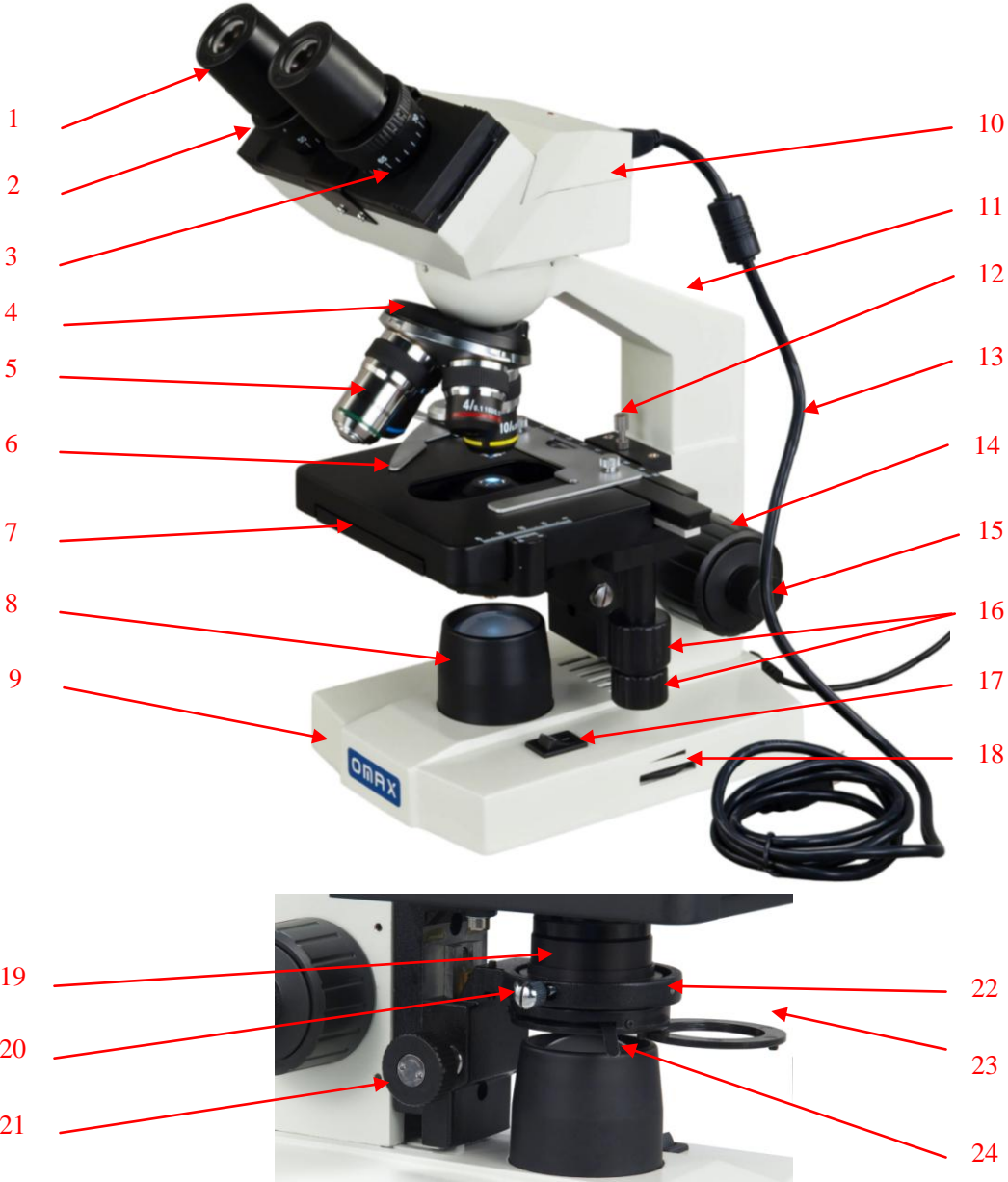

- | | |
|----|----|
| 1 | 13 |
| 2 | 14 |
| 3 | 15 |
| 4 | 16 |
| 5 | 17 |
| 6 | 18 |
| 7 | 19 |
| 8 | 20 |
| 9 | 21 |
| 10 | 22 |
| 11 | 23 |
| 12 | 24 |
-
- | | | |
|----|----|----------------------------|
| 1 | 21 | Condenser Control Knob |
| 2 | 22 | Condenser Holder |
| 3 | 23 | Color Filter Holder |
| 4 | 24 | Iris Diaphragm Lever |
| 5 | | |
| 6 | | |
| 7 | | |
| 8 | | |
| 9 | | |
| 10 | | |
| 11 | | |
| 12 | | |
| 13 | | USB Cable |
| 14 | | Coarse Focus Knob |
| 15 | | Fine Focus Knob |
| 16 | | X-Y Stage Moving Knobs |
| 17 | | Power Switch |
| 18 | | Brightness Intensity Dial |
| 19 | | Abbe Condenser |
| 20 | | Condenser Lock Thumb Screw |
| 21 | | Condenser Control Knob |
| 22 | | Condenser Holder |
| 23 | | Color Filter Holder |
| 24 | | Iris Diaphragm Lever |

2 Installation

2.1 Installation of the eyepieces

- 1) Remove the protective caps from the eyepiece tubes.
- 2) Insert the eyepieces into the eyepiece tubes.

2.2 Installation of the glass filter

- 1) Swing out the color filter holder under the condenser.
- 2) Place the filter into the holder as shown in **Fig. 1**, swing the holder in.

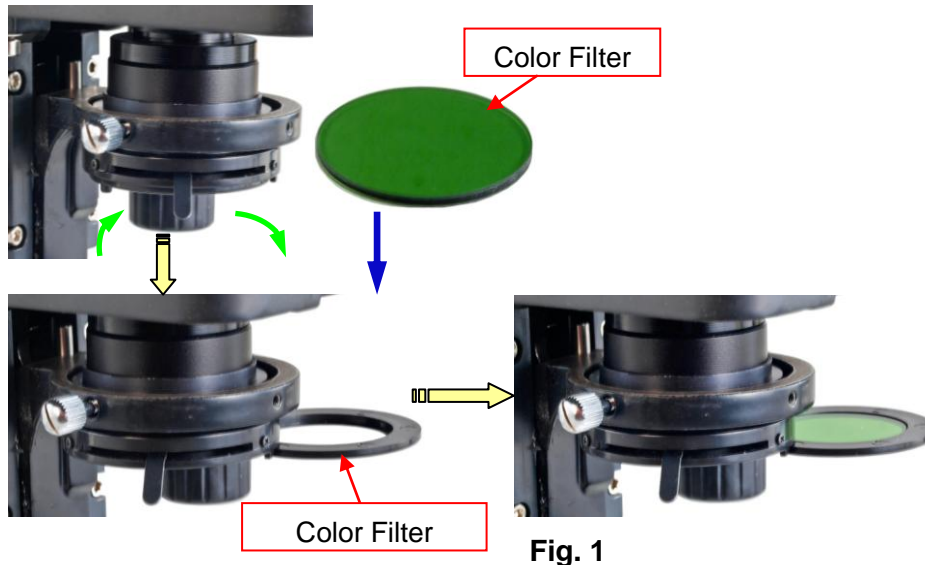


Fig. 1

- ### 2.3 Installation of the dry darkfield condenser (optional, may not included in your package)
- See section 6 (*Dry Darkfield Condenser Installation & Operation Instruction*)

3 Operation

3.1 Adjusting illumination

- 1) Plug the power cord into the power socket on the microscope and connect it to the power outlet.
- 2) Turn on the power switch.
- 3) Rotate the brightness intensity dial to increase or decrease the brightness.

Caution:

A diffusion filter is attached beneath the condenser to get uniform light and protect your eyes from strong LED light when a low power objective applies. The diffusion filter can be swung out to make the view field brighter when observing with a high power objective, such as 100X objective.

3.2 Placing specimen

- 1) Place the slide on the mechanical stage.
- 2) Use the slide holder to gently secure the slide.
- 3) Turn the X and Y stage moving knobs to position the specimen in the center of viewing field.

Caution:

Be sure not to allow an objective to touch a specimen slide when changing objectives.

3.3 Adjusting interpupillary distance

While observing with both eyes, hold the left and right eyepiece tubes and slide the tubes in and out. The interpupillary distance is correct when the left and right fields of view converge completely into one image.

3.4 Adjusting eyepiece diopter

- 1) Rotate the 10X objective into position.
- 2) Rotate the diopter ring on the right eyepiece tube until its numerical value is the same as your interpupillary distance, for example, **70** in the right figure (See **Fig. 2**).
- 3) Close your left eye and bring the specimen into focus following the focusing procedures in 3.5.
- 4) Close your right eye and bring the same specimen into clear sharp focus by adjusting the diopter ring on left eyepiece tube only. Do not use focus knobs at this step.
- 5) Since both sides are adjustable, you may also do the above in the opposite way, in other words, left eye first and right eye second.

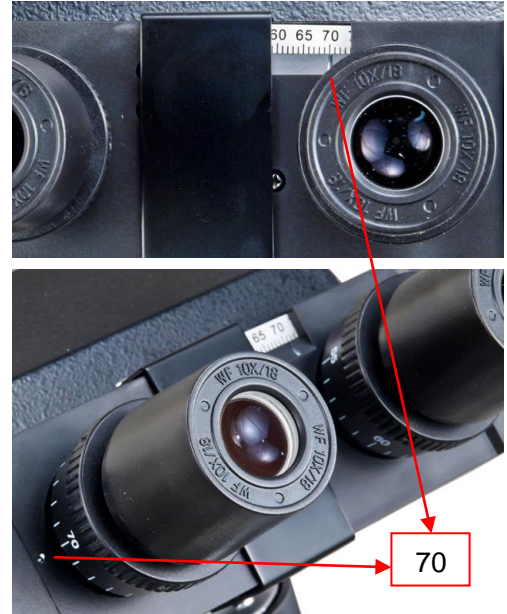


Fig. 2

3.5 Focusing

- 1) With the 10X objective in position, raise the mechanical stage using the coarse focus knob until the specimen is close to the objective
- 2) Turn the coarse focus knob until the specimen is in focus. Use the fine focus knob to obtain a sharp image. You may now switch to another magnification objective.
- 3) To get a good focused image, you may need to combine the focus knob adjustment and interpupillary distance adjustment, along with eyepiece diopter adjustment stated in **3.3** and **3.4**.

Tips:

To prevent your specimen slide from making contact with an objective, turn the 100X objective in position and adjust the focusing rack stop screw (**Fig. 3**) so that the 40X objective will not contact the specimen while the stage is adjusted to its highest position. Give the stage a tiny extra moving space to ensure the objective can be focused every time.



Fig. 3

3.6 Adjusting condenser

- 1) Turn the condenser focus knob to raise or lower the condenser.
- 2) Raise the condenser when using high power objectives and lower it when using low power objectives.

Note:

- The centering of the condenser and the light axis of the objective are factory adjusted. Do not attempt re-adjustment.
- The highest position of the condenser has been factory adjusted. Do not attempt re-adjustment.

3.7 Adjusting iris aperture diaphragm

Swing the iris diaphragm lever (**Fig. 4**) left or right to adjust the aperture size.

Note:

The iris diaphragm is designed to adjust the aperture size, not to adjust the brightness although the brightness will be changed when it's adjusted. When aperture is adjusted to smaller size, the contrast will be increased and the depth of field will be increased as well. Turn up the intensity of the light if the image is too dim.

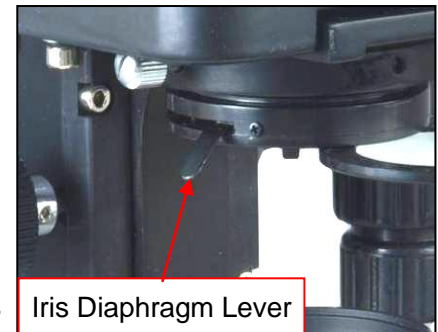


Fig.4

3.8 Photo/video observing, capturing and recording

- 1) Bring the microscope into focus by following the procedures in 3.5.
- 2) Insert the USB cable into the USB port (**Fig. 5**) on the back of viewing head, and the other end to the computer.
- 3) Turn on the computer; install the camera following the manual in the mini CD.
- 4) Open image observing software to examine (more details see camera's manual).
- 5) Capture images with manual white balance for image size of 1280X1024 or 1024X768 resolution when using auto white balance.
- 6) You also can record live videos through the software.

Note:

- Please refer to the manual in the camera's CD for the details of installation and operation of the camera.
- Do not capturing images of 1280X1024 resolution in auto white balance mode. The captured images may have problems with color rendering at this mode.



Fig. 5

4 Specifications

General	
Model	MD82EE10
Total Magnification	40X, 100X, 200X, 400X
Viewing Head	Binocular, inclined 45°, 360° swiveling with Built-in camera
Interpupillary Distance	Sliding adjustment, 2-3/16" ~ 2-15/16" (55mm ~ 75mm)
Diopter Adjustment	On both eyepiece tubes
Eyepieces	1 pair of wild field WF10X/18
Nosepiece	Revolving quadruple
Objectives	Achromatic DIN 4X/0.10 160/0.17 10X/0.25 160/0.17 20X/0.40 160/0.17(spring) 40X/0.65 160/0.17(spring)
Condenser	Abbe, NA=1.25, w/ iris diaphragm and diffusion filter Rack and pinion adjustment
Focus Mechanism	Coaxial coarse and fine focusing knobs on both sides with rack stop
Mechanical Stage	Double layer, dimension: 4-1/2" x 4-15/16"(115 mm x 125 mm) X-Y translational range: 2-13/16" x 1-3/16" (70 mm x 30 mm)
Illumination	Transmitted, LED, variable intensity
Camera	Built-in USB2.0 1280 x 1024 pixel (1.3MP) Driver and Software included in the CD Compatible with Windows 2000, XP, Vista and Windows 7 32/64-bit
Power Supply	AC/DC adapter, 100V-240V
Dimension	9-1/16" x 7-1/8" x 13" (23cm x 18cm x 33cm)
Net weight	7 lbs 5 oz (3.32 kg)

5 Troubleshooting Guide

Problem	Cause	Solution
Lamp does not light when switched on	No electrical power	Check power cord connection
	LED or power unit dead	Contact seller for service
Darkness at the periphery	Revolving nosepiece not in click stop position	Revolve the nosepiece to click-stop position by swinging the objective correctly into the optical path
Uneven brightness in the field of view	The diffusion filter is not in the light path	Check the diffusion filter and make sure it's in the light path
Dirt or dust on the view	Dirt or dust on the lens eyepiece, condenser, objective, collector lens or specimen	Clean the lens with a lens cleaning paper
	Dirt or dust on the specimen	Clean the specimen
Poor image quality or not able to get focused image	No slide cover attached to the slide	Attach a 0.17mm slide cover
	Slide cover is too thick or thin	Use a slide cover of the appropriate thickness (0.17mm)
	Slide may be upside down (specimen at the bottom)	Turn slide over so the cover-glass faces up
	Diopter adjustment is not set properly	Readjust the diopter settings
	Specimen rises from stage surface	Secure the specimen in the slide holder
	Condenser aperture is closed or open too much	Open or close properly
	Condenser is positioned too low or too high	Move condenser upper or lower
	Lamp intensity is too high or low	Adjust the light intensity by rotating the intensity control dial
Image is black screen through	No enough lighting or driver not installed properly	Check the camera manual
Image is totally white through	The light is too bright	Manually adjust the exposure in the software. See manual for details.