

VFC 400 Vaccine Monitoring Data Logger kit

Step 1: Install the probe in the fridge/freezer

Presented by
Control Solutions
(503) 410-5996

Install the probe in the fridge/freezer

- Place acrylic stand and probe vial and near the middle of the fridge/freezer
- Route the cable underneath the wire rack and secure it with a zip tie
- Route the cable toward the wall of the hinge side and secure with a zip tie
- Continue to route the cable towards front of the fridge/freezer on the hinge side and secure
- (see the following slides for details)
- Note: If you have a cable port then run it through there and secure to rack with zip ties



Install the probe in the fridge/freezer



Zip Tie



Install the probe in the fridge/freezer



Zip Tie



Install the probe in the fridge/freezer

- Place the logger into the wall mounting bracket, plug in the sensor wire to the logger (left side)
- Adhere the mounting bracket for the VFC 400-1 to either the side or front of the door in a spot that is easy to view & operate.
- Approx. 6 inches underneath the logger, Adhere cable tie bracket and secure the cable with a Zip tie.
- Leave enough slack in the cable so you can plug and unplug the VFC400-1 easily

