

# VFC 400 Vaccine Monitoring Data Logger kit

## Step 2 VTMC Software Download

Presented by  
Control Solutions  
(503) 410-5996

# VFC 400 software download

- From the internet, go to our website [www.vfcdataloggers.com](http://www.vfcdataloggers.com) to download the latest software revision.
- Click on **Software Downloads** and then click on “VFC400 software download”

www.vfcdataloggers.com

CSI Manufactures Personal Google

**VFC Dataloggers**  
888-311-0636

HOME MY ACCOUNT SIGN IN OR CREATE AN ACCOUNT VIEW CART SEARCH FOR A PRODUCT

VFC COMPLIANT LOGGERS ALL PRODUCTS SUPPORT CALIBRATION **SOFTWARE DOWNLOADS**

Welcome to our newly updated website!

**MEET THE NEW STANDARD**

Introducing the VFC 400

[Learn more](#)

**VFC 400**  
Vaccine Temperature Data Logger  
Control Solutions, Inc.  
www.vfcdataloggers.com  
(888) 311-0636

TIME 13:29  
TEMP 12.9°

REVIEW MARK  
START CLEAR STOP




# VFC 400 software download

- Scroll down and click on “Click to Download”




When the box shows up as shown below, click on “Save File”.

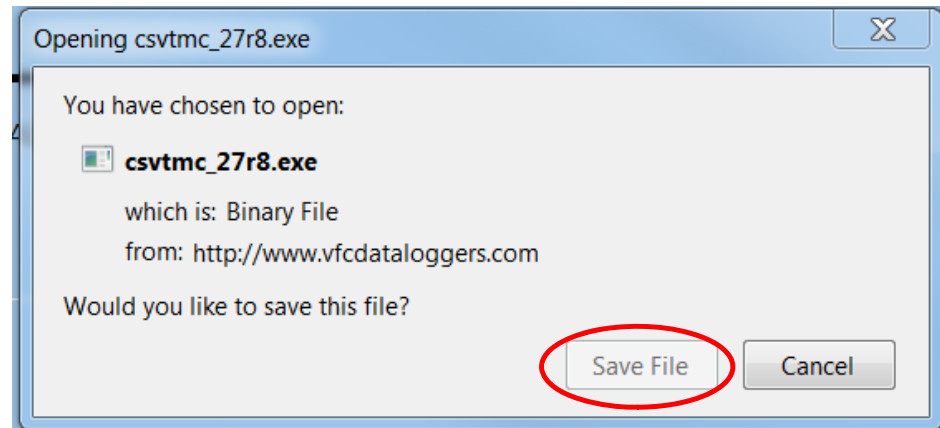


# VFC 400 software download

- Go to where you saved the file to and click on the file to start the download. You will need administration rights to do this and may need IT support.
- The file is saved as a zip file. Double click on the zip file and then double click on file “csvtmc\_27r8.exe”

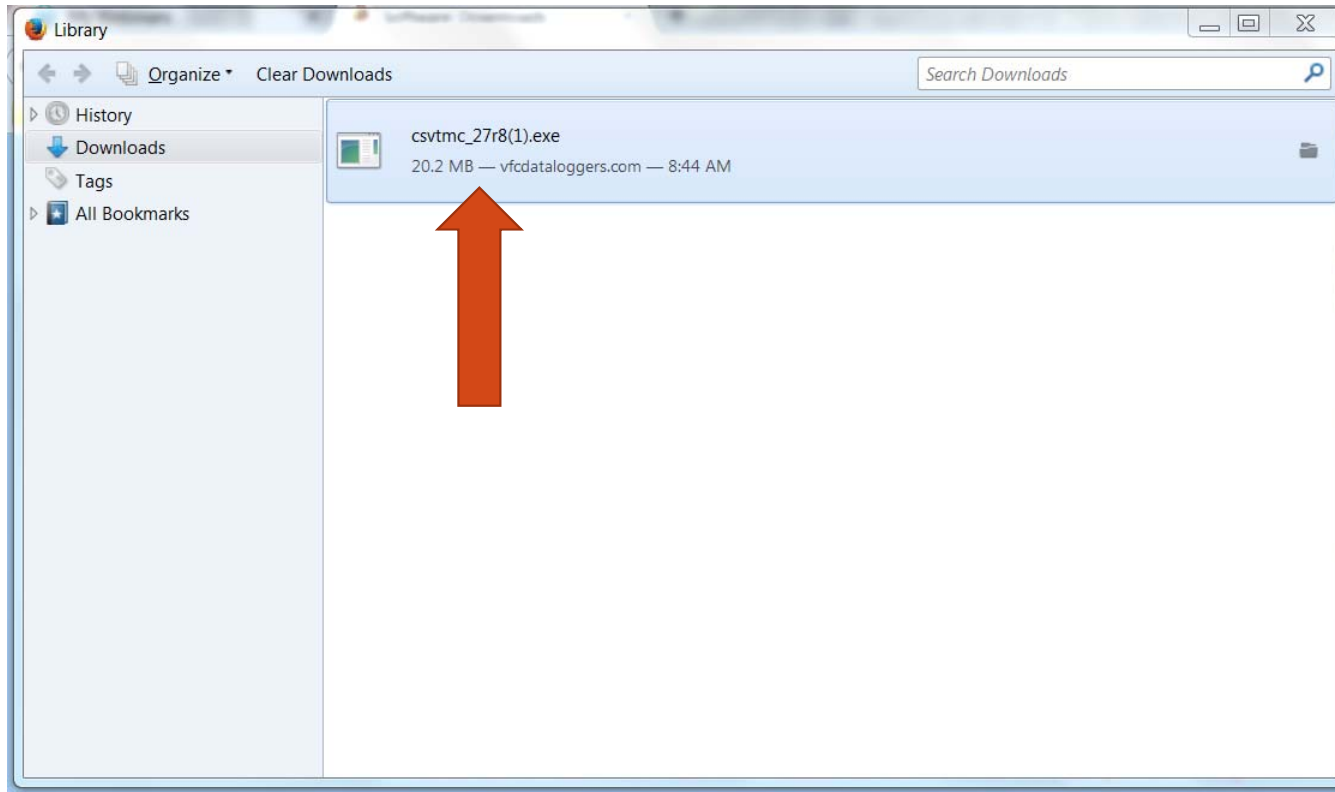


# VFC 400 software download



- A pop-up will appear that asks “Do you want to allow the following program to make changes to this computer?”. Click yes.
- The next series of pop-ups click as shown by the arrows

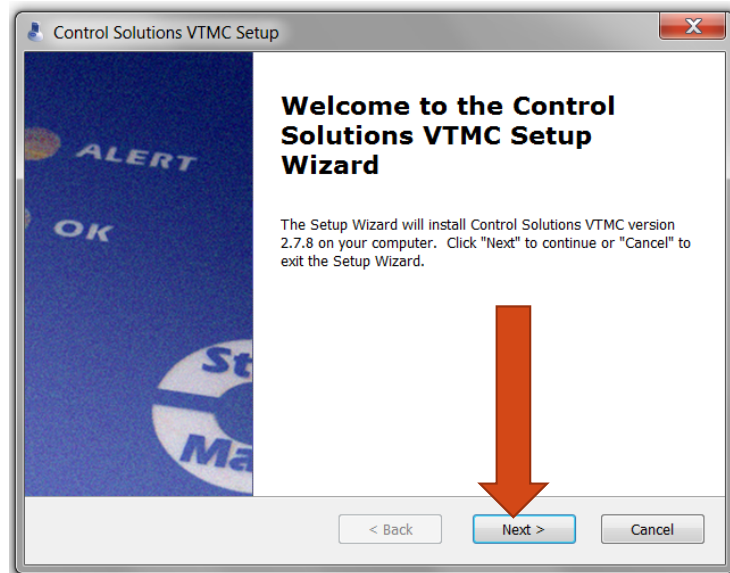


# VFC 400 software download

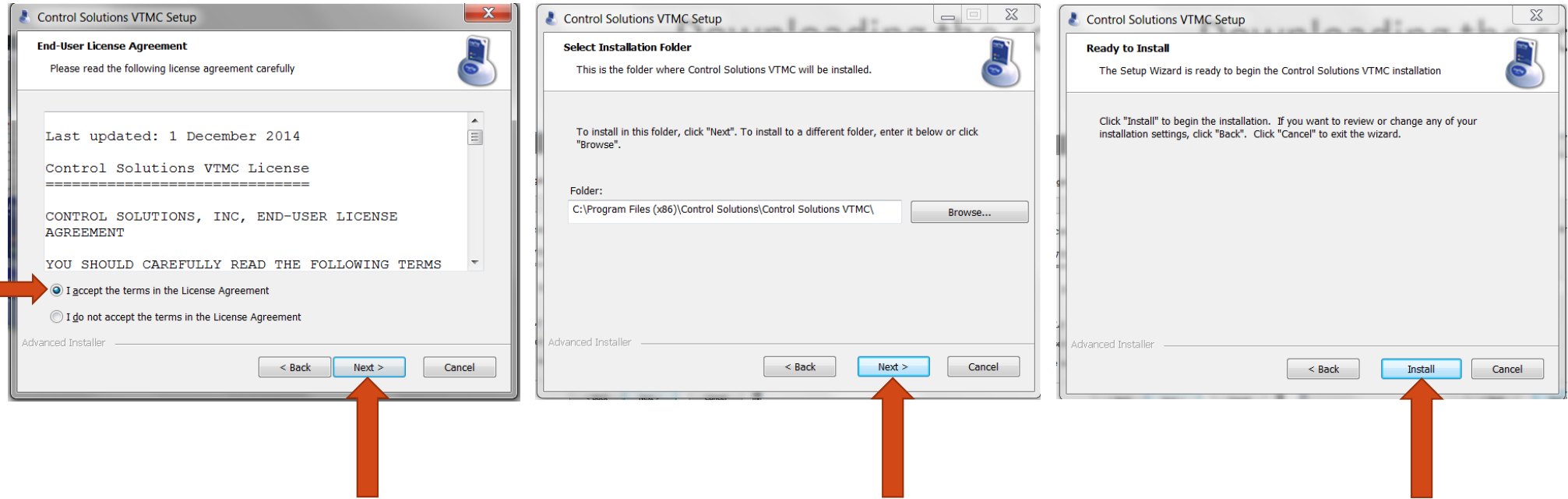

# VFC 400 software download

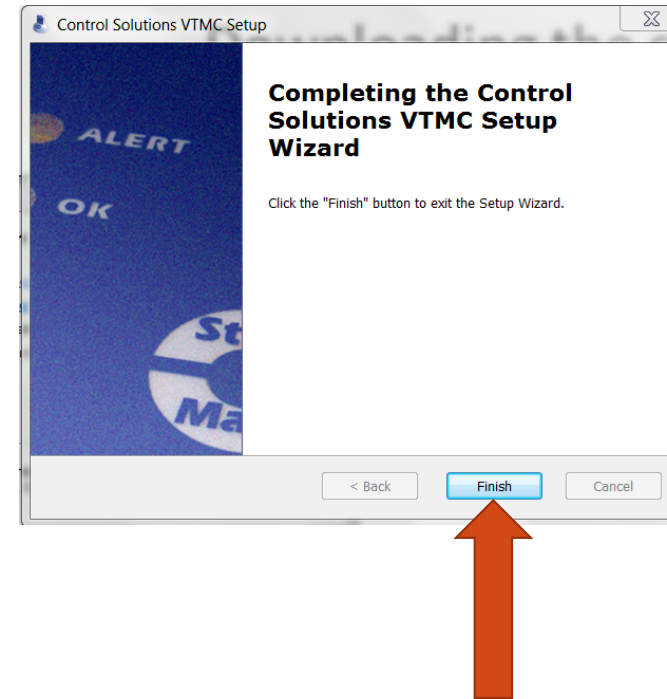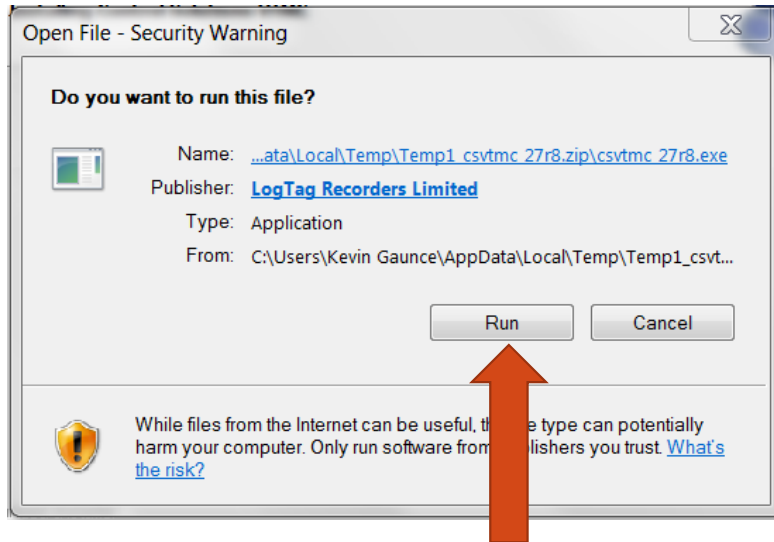

# VFC 400 software download

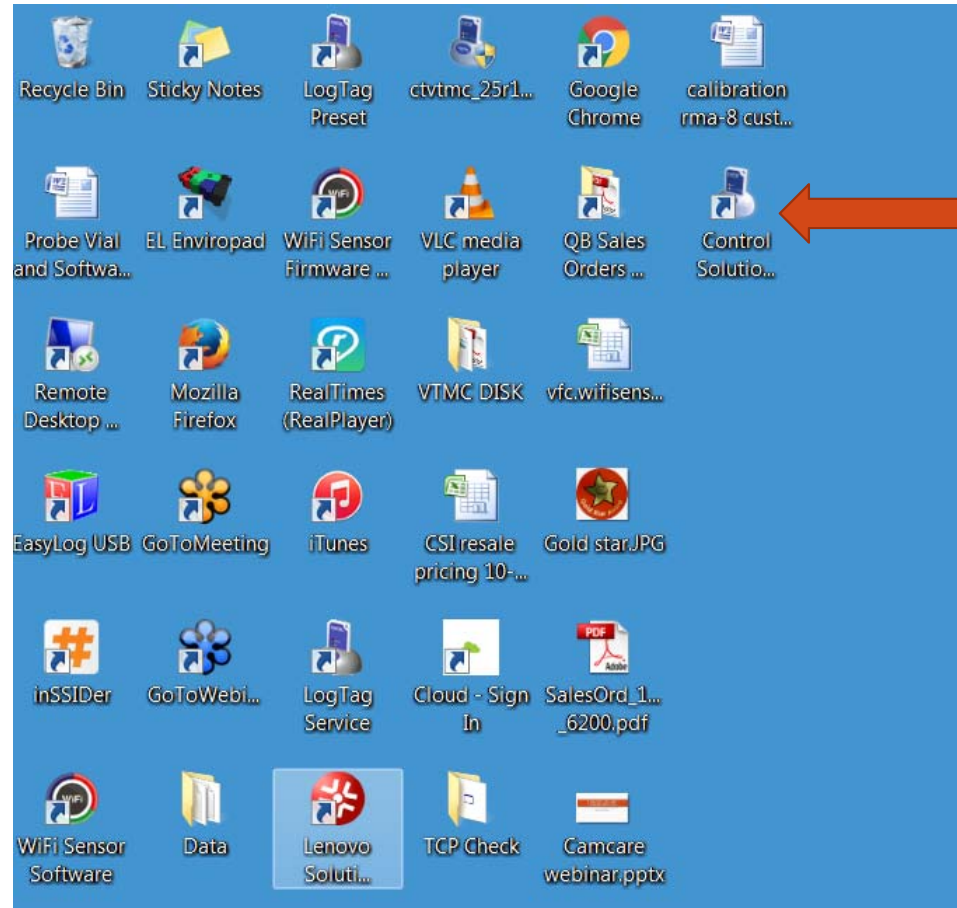

You should now have an icon on your desktop that looks like this.

