VFC400 Installation



Your kit includes:

- VFC400 Data Logger
- Stainless Steel Probe encased in glycol
- Acrylic stand for probe and mounting equipment for logger
- Adhesive backed zip tie mounts and zip ties for securing cable
- Spare battery
- 2 year NIST traceable Certificate of Calibration compliant to ISO 17025: 2005






1

Place acrylic stand and probe vial near the middle of the fridge/freezer Route the cable under the wire rack and secure with zip tie

Route the cable toward the wall of the hinge side and secure with zip tie

Route the cable toward the front of the fridge/freezer on the hinge side and secure

**Place the glycol bottle in the fridge/freezer for <u>at least 1.5 hours</u> prior to starting your logger to allow the solution to reach the appropriate temperature.

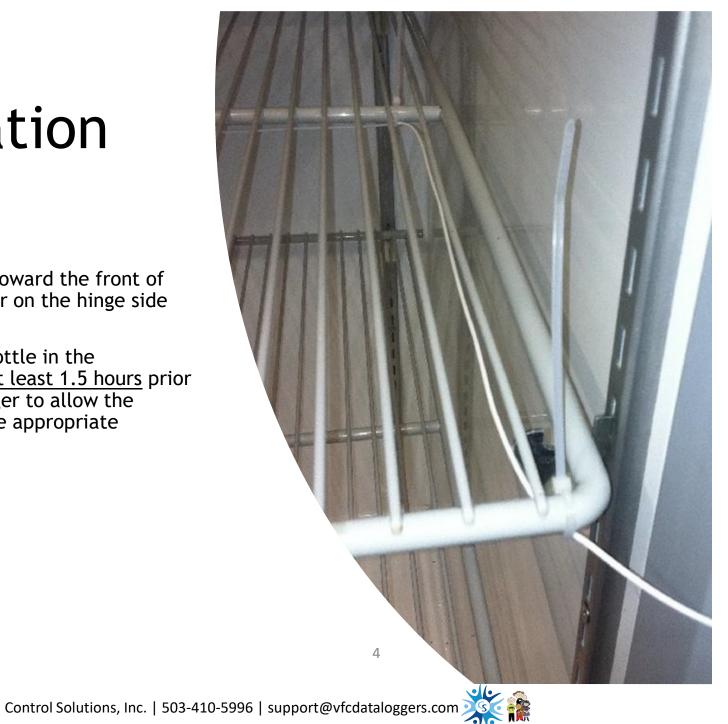



- Adhere the mounting bracket to the side or front of your fridge/freezer
- Place the logger into the mounting bracket and plug the sensor wire into the logger (left side)
- Approx. 6 inches underneath the logger, adhere cable tie bracket and secure the cable with a Zip tie. Leave enough slack in the cable so you can plug and unplug the VFC400 easily

Questions about installation?

Visit our support page

Send us an email support@vfcdataloggers.com

Jump to:

Data Logger Installation
Software Download and Setup
Data Logger Configuration
Data Logger Operation
FAQs and Troubleshooting



VFC400 Software Download and Setup



Software Download and Set-Up

Visit our website vfcdataloggers.com And choose Software Downloads VFC COMPLIANT LOGGERS ALL PRODUCTS SOFTWARE DOWNLOADS CALIBRATION SUPPORT CONTACT US

Home > Software Downloads

SOFTWARE DOWNLOADS

Control Solutions

VFC 400 SOFTWARE (VTMC)

Download the most current software

NEW Control Solutions VTMC Software. Version 2.9 revision 5 (Click to download) **NEW**


LOGTAG SOFTWARE (VTMC)


After clicking the software link your download will begin immediately


NEW Control Solutions VTMC Software. Version 2.9 revision 5 (Click to download) **NEW**


Click on the download in progress to prompt the install wizard to open when download has completed


The following steps will appear; advance through the installation process.

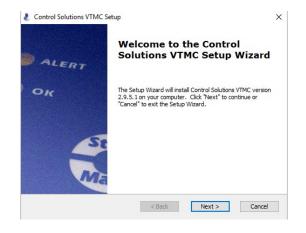

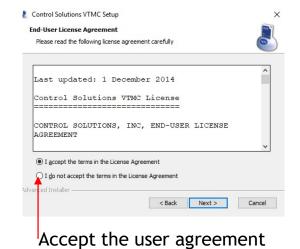

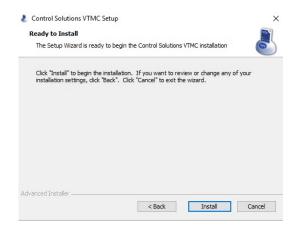

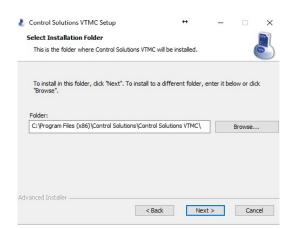

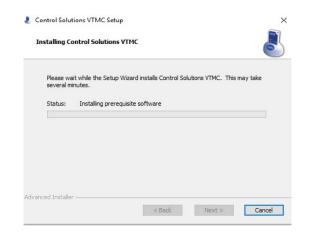




If the install wizard does not appear automatically after download navigate to your downloads folder and choose the newly downloaded file by double clicking



Desktop Icon

After successful software download and installation you will have a Control Solutions icon on your desktop

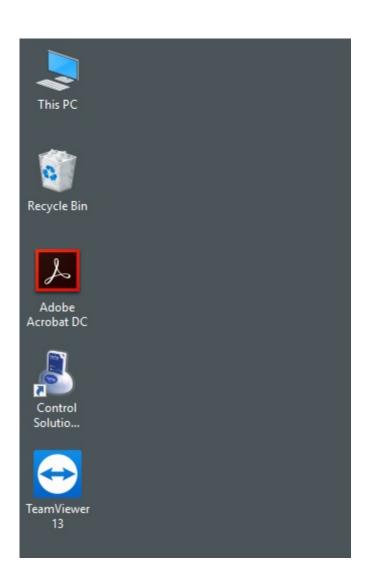

Questions about software download?

Visit our support page

Send us an email support@vfcdataloggers.com

Jump to:

VFC400 Installation

VFC400 Software Download

VFC400 Software Set Up

VFC400 Logger Configuration

Data Logger Operation

