



VFC400 Software Setup

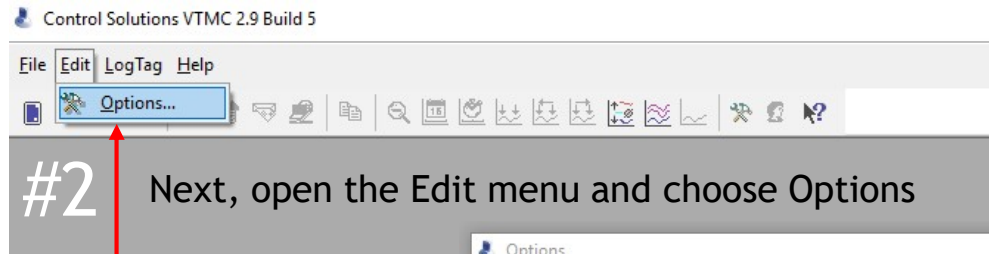



Open Control Solutions VTMC

by clicking on the icon on your desktop

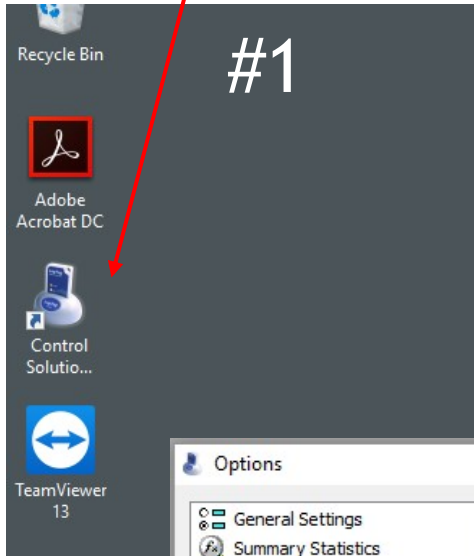

Next, open the Edit menu and choose Options

#3 From the Options menu choose General Settings and select your preferred temperature scale: F° or C°*

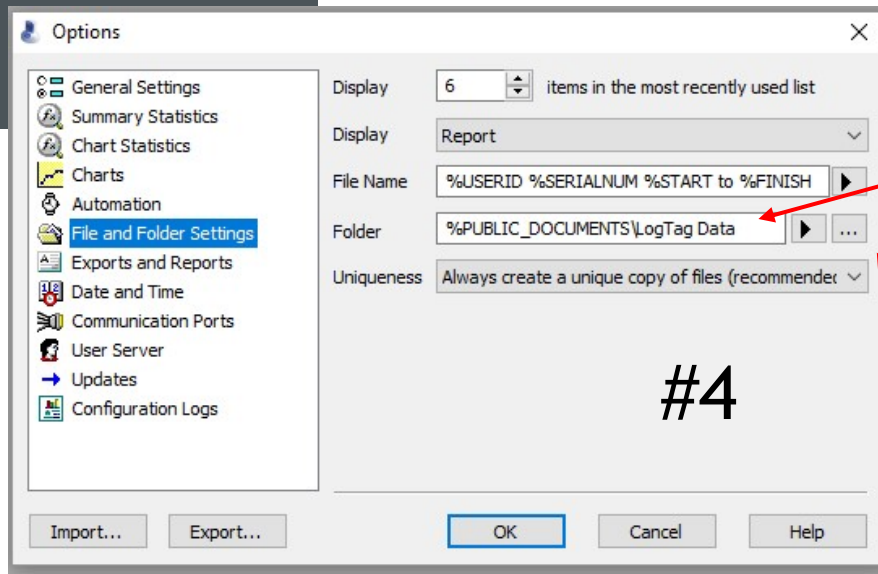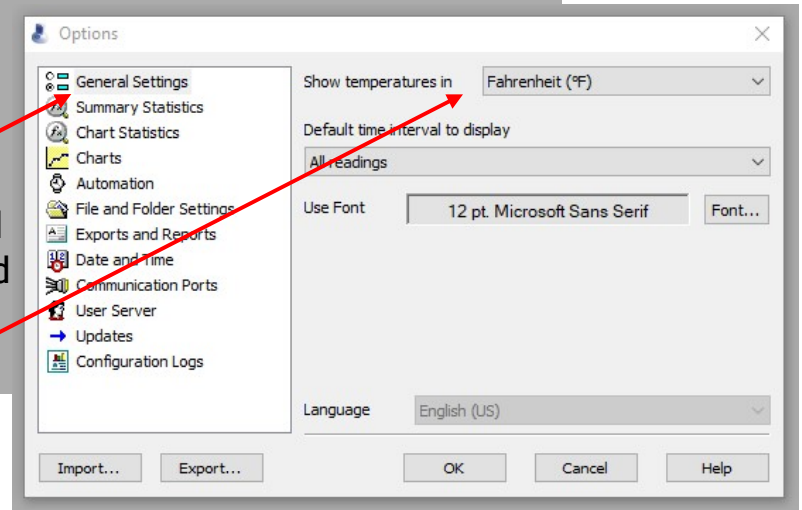

Data will automatically be saved to the newly created folder Public>Documents>LogTag Data every time your logger is placed in the docking station. Anyone logged into the PC can access the data.

If you prefer to save the data to a network or shared drive, click the button with 3 dots(...). A box will appear saying "Locate Folder". Simply go to the folder you want to save the data to and click "OK".




Choose OK

to complete software configurations

the remaining options do not require changes

Other options are configurable but not necessary

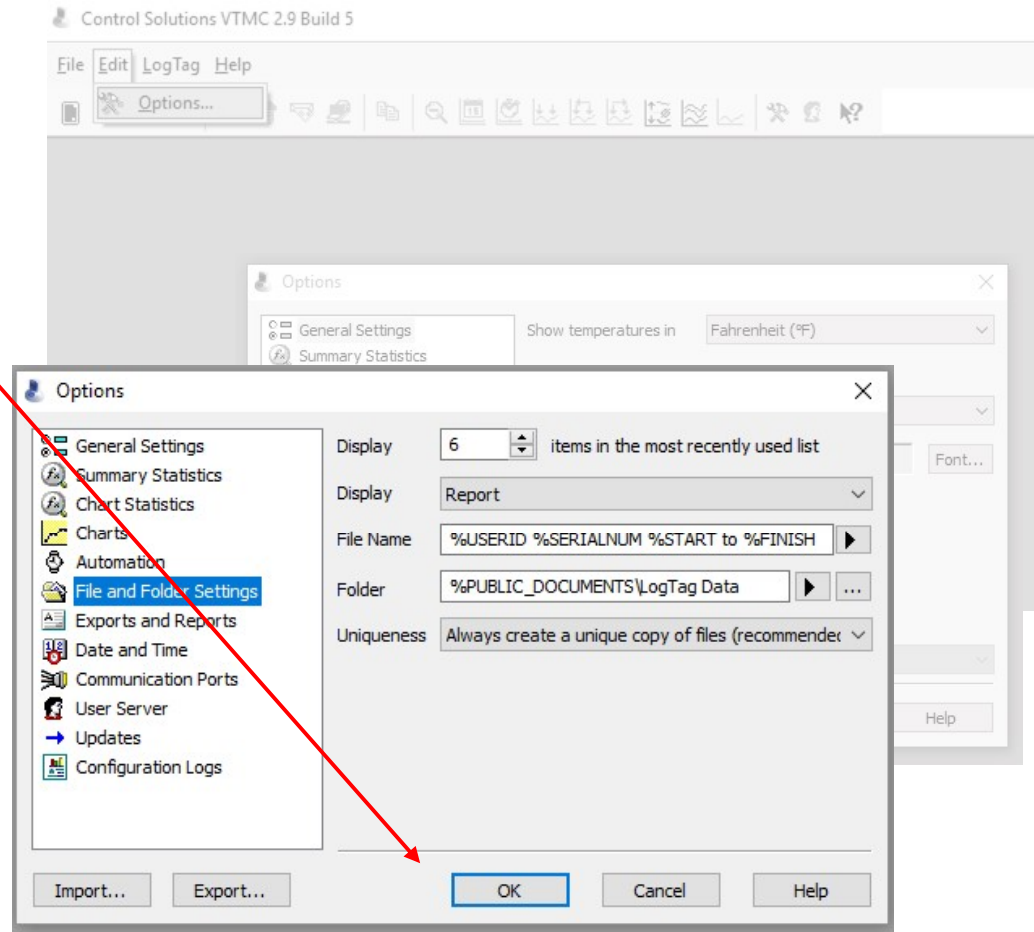

Questions about software configuration?

Visit our [support page](#)

Send us an email
support@vfcdataloggers.com

Jump to:

- Data Logger Installation
- Software Download and Setup
- Data Logger Configuration
- Data Logger Operation
- FAQs and Troubleshooting

